


Unison UK-C18

100 x 4.6 mm


Application

### Chlorophyll and Carotenoid


壬生菜アセトン抽出物中のクロロフィルとカロテノイド

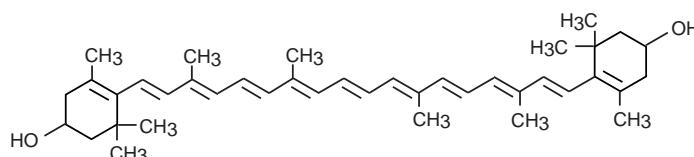


1 lutein (xanthophyll)

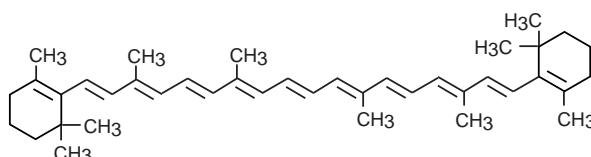


4 β-carotene

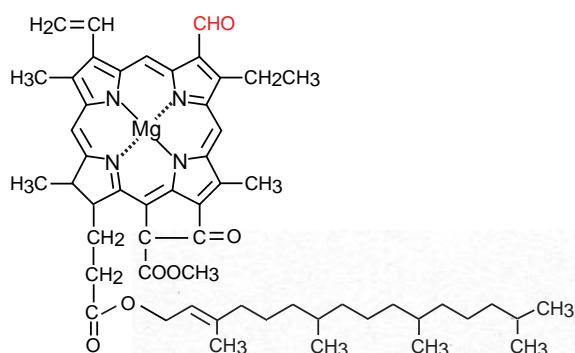

2 chlorophyll b

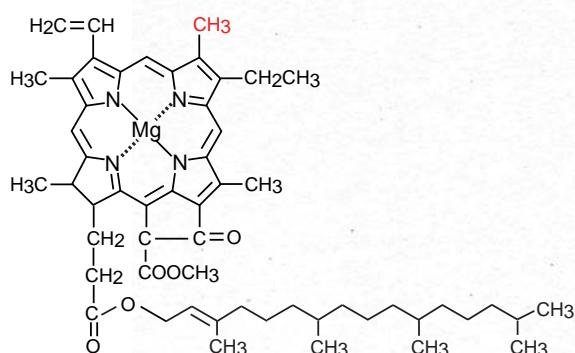

3 chlorophyll a

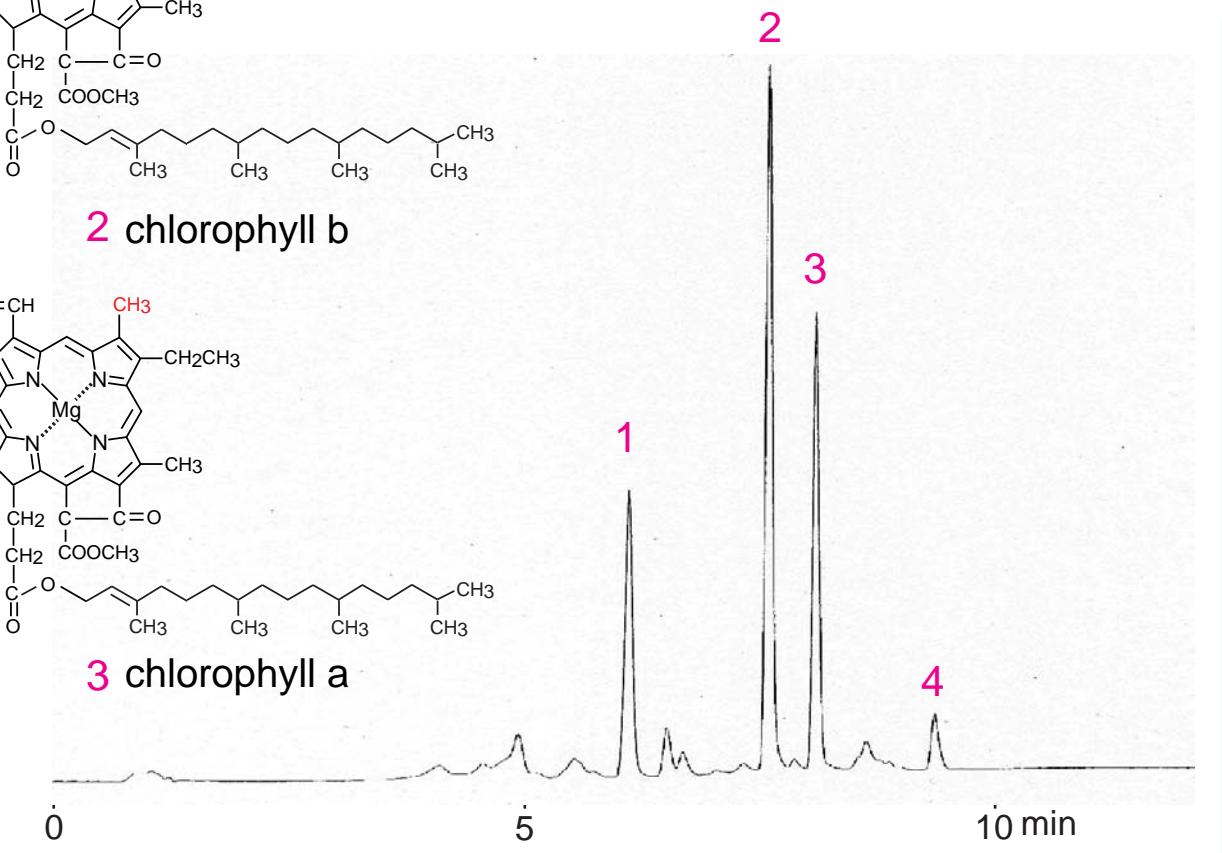

Unison UK-C18, 100 x 4.6 mm

A: methanol / water = 90 / 10, B: ethyl acetate

0-100% B (0-7min), 100% B (7min - ), 1 mL/min, ambient, VIS 450 nm  
acetone extracts from *Brassica campestris*

Courtesy of K. Nakanishi, Tokushima Prefectural Industrial Technology Center, Japan