





**Scherzo SM-C18**  
Unison UK-C18

150 x 3 mm


Application

**Cold medicine compounds**

風邪薬成分

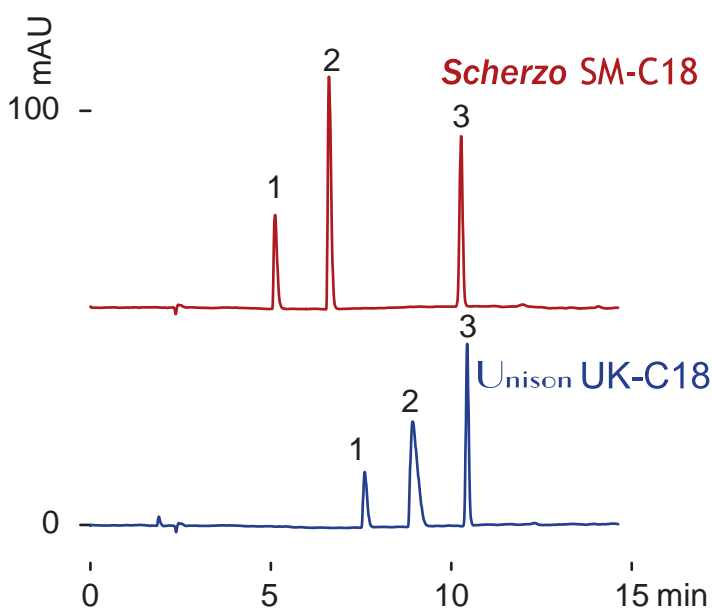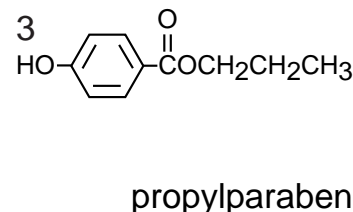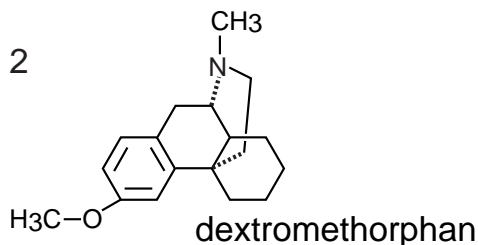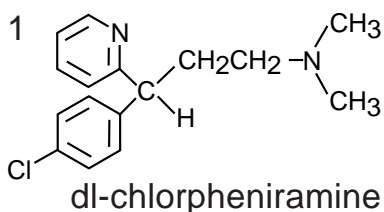

A: water / formic acid = 100 / 0.1  
B: acetonitrile / formic acid = 100 / 0.5  
15-80 %B (0-10min)  
80 %B (10-13min)  
15 %B (13-30min)  
mixer: 600uL  
0.4 mL/min (8MPa)  
37 deg.C, 280 nm  
1 uL (0.2-0.9ug)

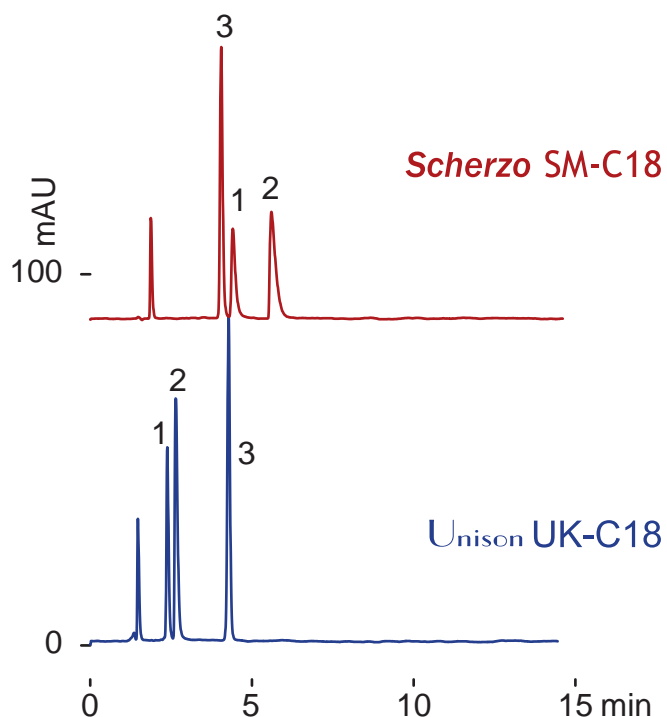

100mM ammonium formate / acetonitrile = 45 / 55  
0.4 mL/min (8MPa)  
37 deg.C, 235nm,  
1 uL (0.2-0.9ug)