


**Intrada** Amino Acid

10 x 2 mm

Application

LC-MS: 1min High Throughput Analysis amino acids

アミノ酸の超高速LC-MS分析

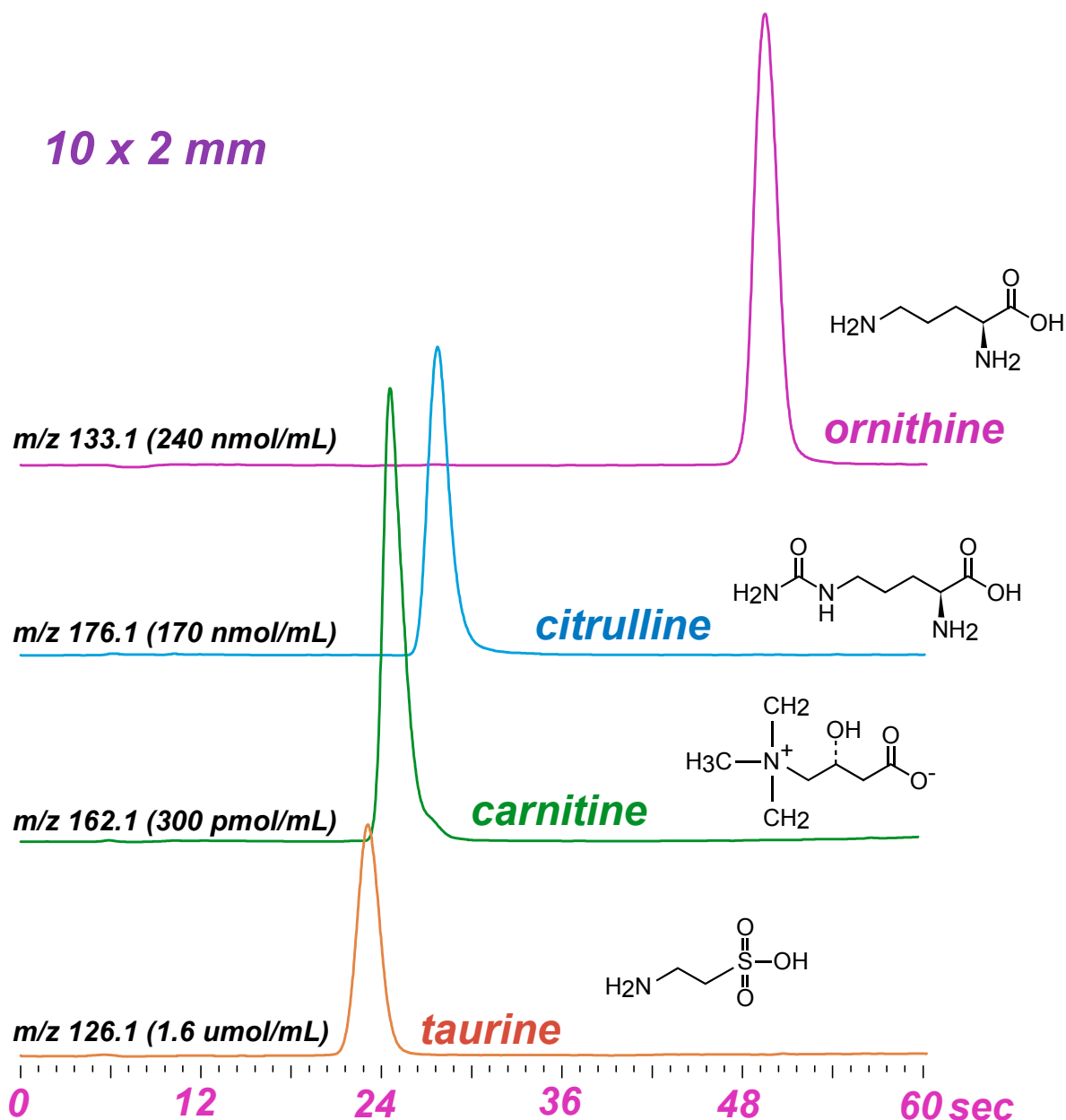

Intrada Amino Acid, 10 x 2 mm

A: acetonitrile / formic acid = 100 / 0.1

B: 100mM ammonium formate

15-100%B (0-0.8min), 100 %B (0.8-1.0min)

0.4 mL/min (1.6MPa), 35°C, 1 uL (0.1N HCl)

Single Quad. MS (ESI, positive)