


Scherzo SM-C18

50 x 3 mm

Application

LC-MS: biogenic polyamines

生体ポリアミンのLC-MS分析

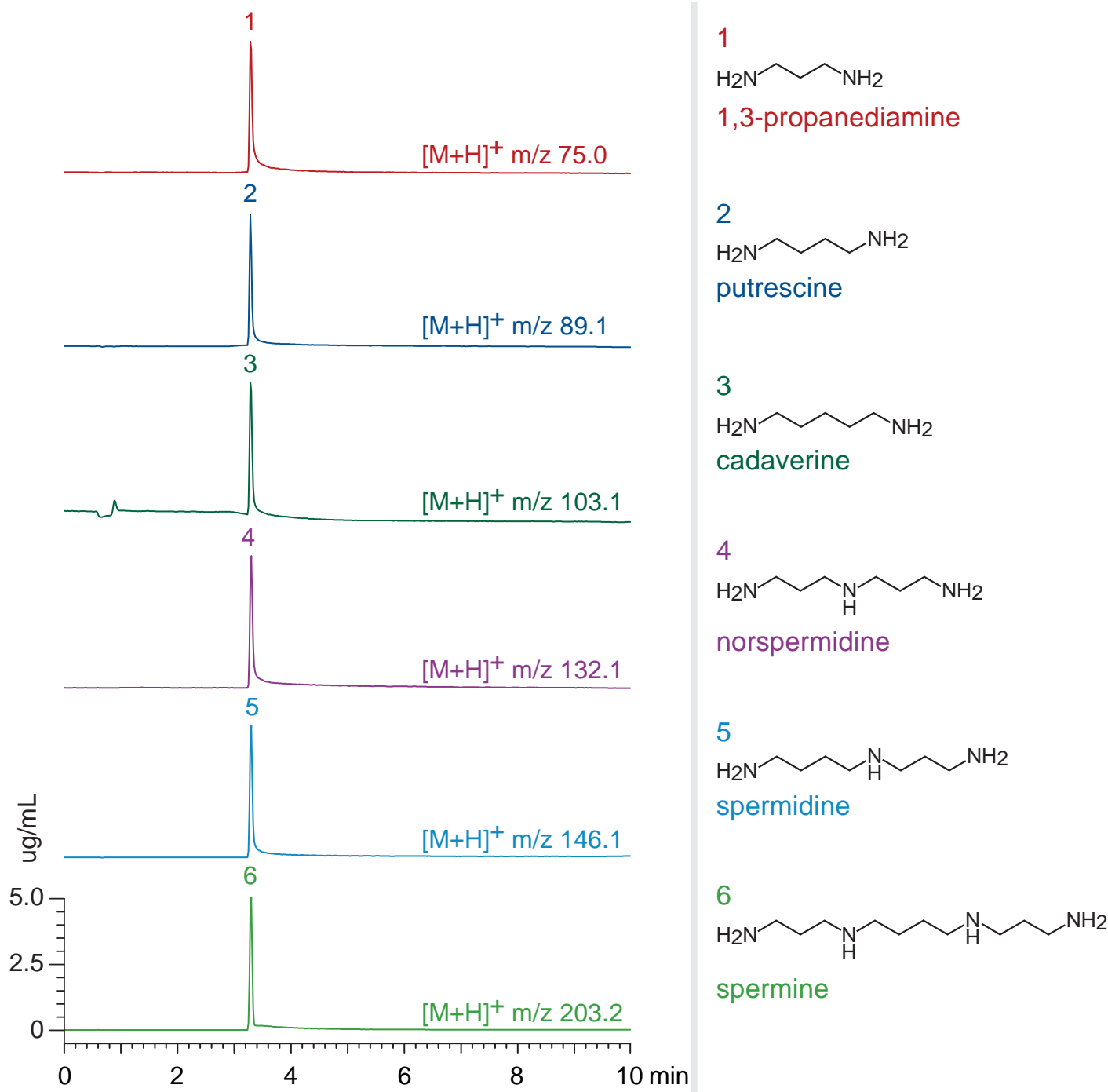

Scherzo SM-C18, 50 x 3 mm
 A: 5 mM ammonium formate
 B: acetonitrile / water / formic acid = 20 / 80 / 0.8
 0 %B (0-1min), 0-70 %B (1-8min), 100 %B (8-10min)
 0.4 mL/min (4 MPa), 37°C, 10uL (5ug/mL, 50mM AcONH₄)
 Single Quad. MS (ESI, positive)