


Unison UK-C18  
Cadenza CD-C18

50 x 2 mm


Application

LC-MS/MS: Insecticide (Fipronil)

フェニルピラゾール系殺虫薬 フィプロニルのLC-MS/MS分析

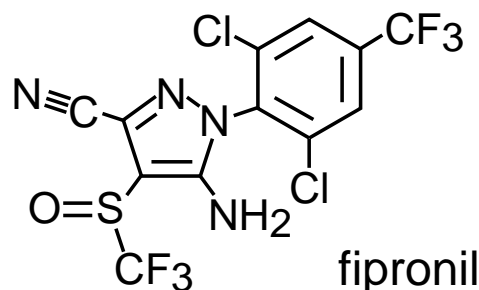

MRM 436.8 > 367.8 m/z

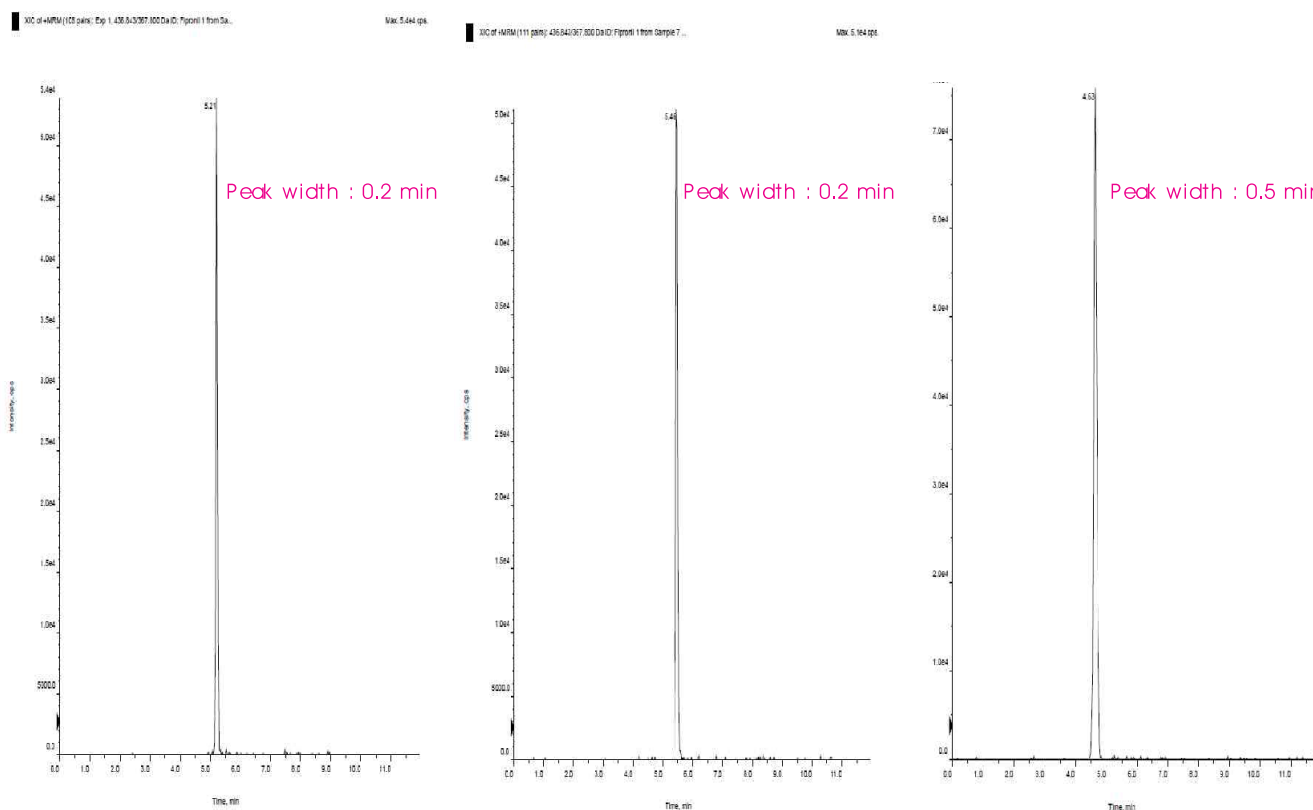

Unison UK-C18

Cadenza CD-C18

BRAND-PX

50 x 2 mm

A: 0.1% formic acid in 5mM ammonium acetate

B: 0.1% formic acid in methanol

5-90%B (0-3min), 90%B (3-7min), 90-5%B (7-7.1min), 5%B (7.1-12min)

0.2 mL/min

2 µL (0.1 µg/mL)

Triple Quad. MS (m/z 436.8 / 367.8)

Courtesy of PEAKMAN SP, KOREA