






Scherzo SM-C18 MF

100 x 2 mm

Application

LC-MS: Organophosphorus pesticides/herbicides

有機リン系農薬のLC-MS分析

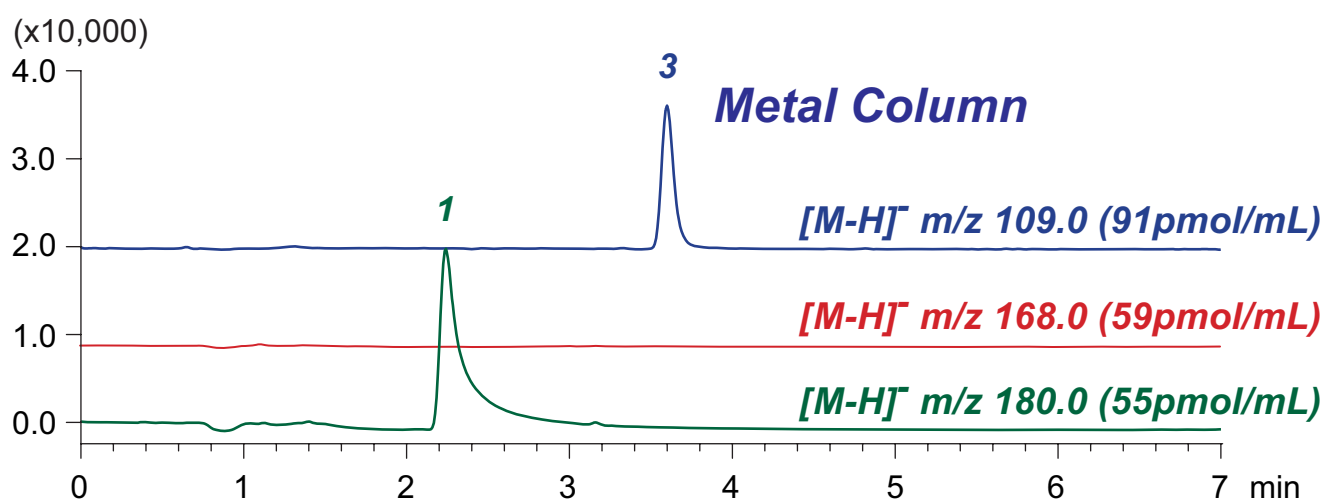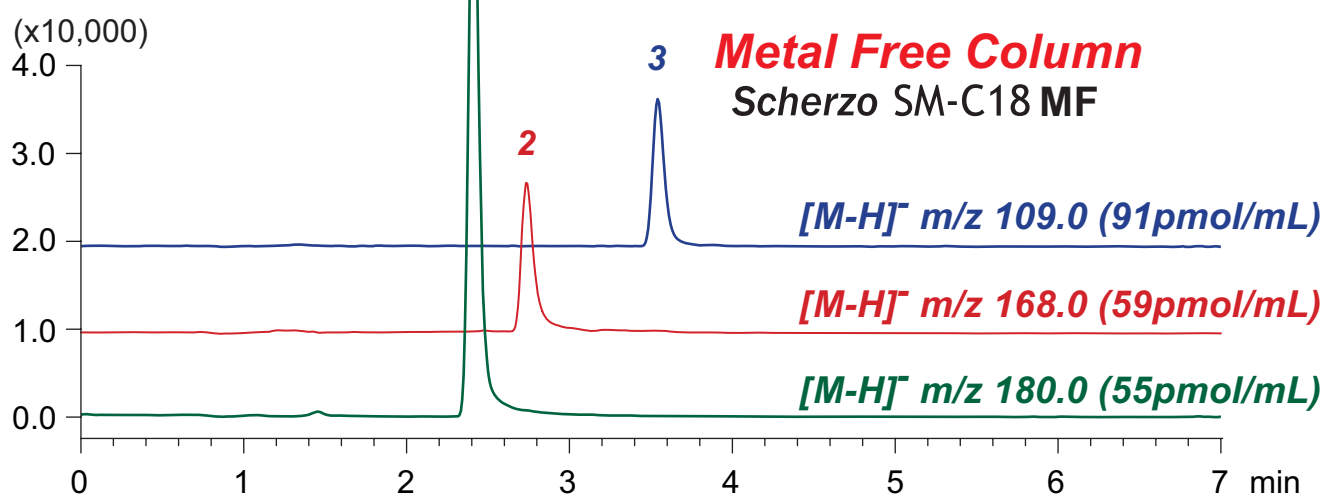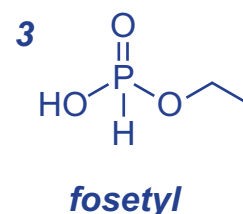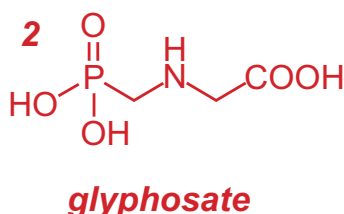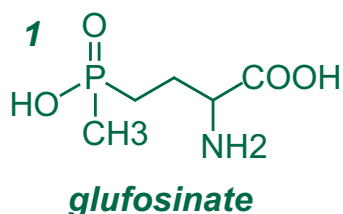

Scherzo SM-C18 MF, 100 x 2 mm
 A: methanol / water / formic acid = 20 / 80 / 0.1
 B: methanol / 50mM ammonium formate = 20 / 80
 0-100%B (0-5min), 100%B (5-7min)
 0.2 mL/min (10MPa), 37°C, 20 μL (55-91pmol/mL, water)
 Single Quad. MS (ESI)