


**Intrada Organic Acid**

150 x 2 mm

Application

**LC-MS: 44 Organic Acids Analysis (IEX mode)**

44種類有機酸のLC-MS分析(イオン交換モード)

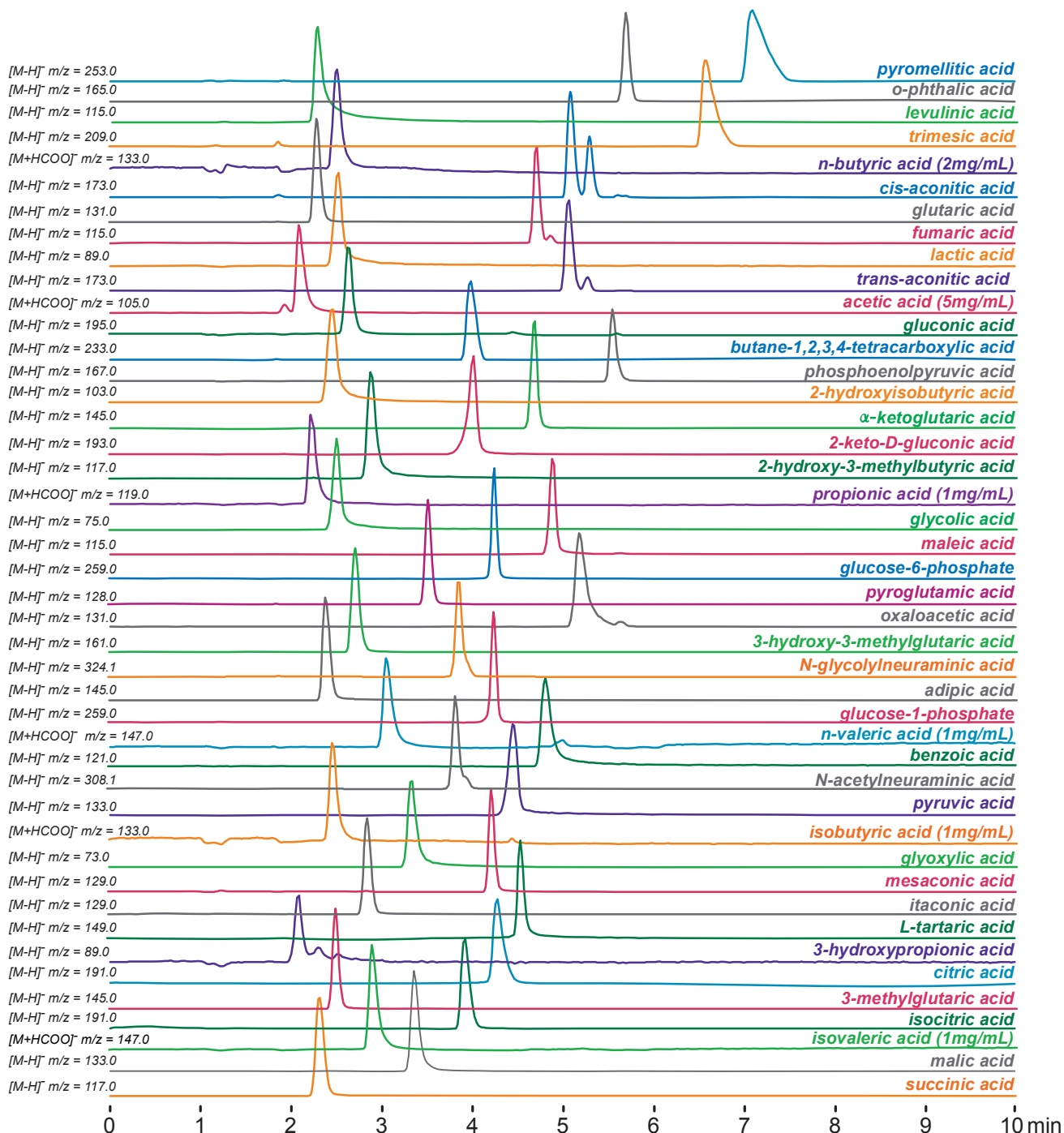

**Intrada Organic Acid, 150 x 2mm**

A: acetonitrile / water / formic acid = 10 / 90 / 0.1

B: acetonitrile / 100mM ammonium formate = 10 / 90

0%B (0-1min), 0-100%B (1-7min), 100%B (7-10min)

0.2mL/min, 37°C, MS (ESI, SIM, Negative)

5µL (100µg/mL with 0.1% HCOOH)