


Intrada Organic Acid

150 x 2 mm

Application

LC-MS: Food preservatives

食品保存料のLC-MS分析

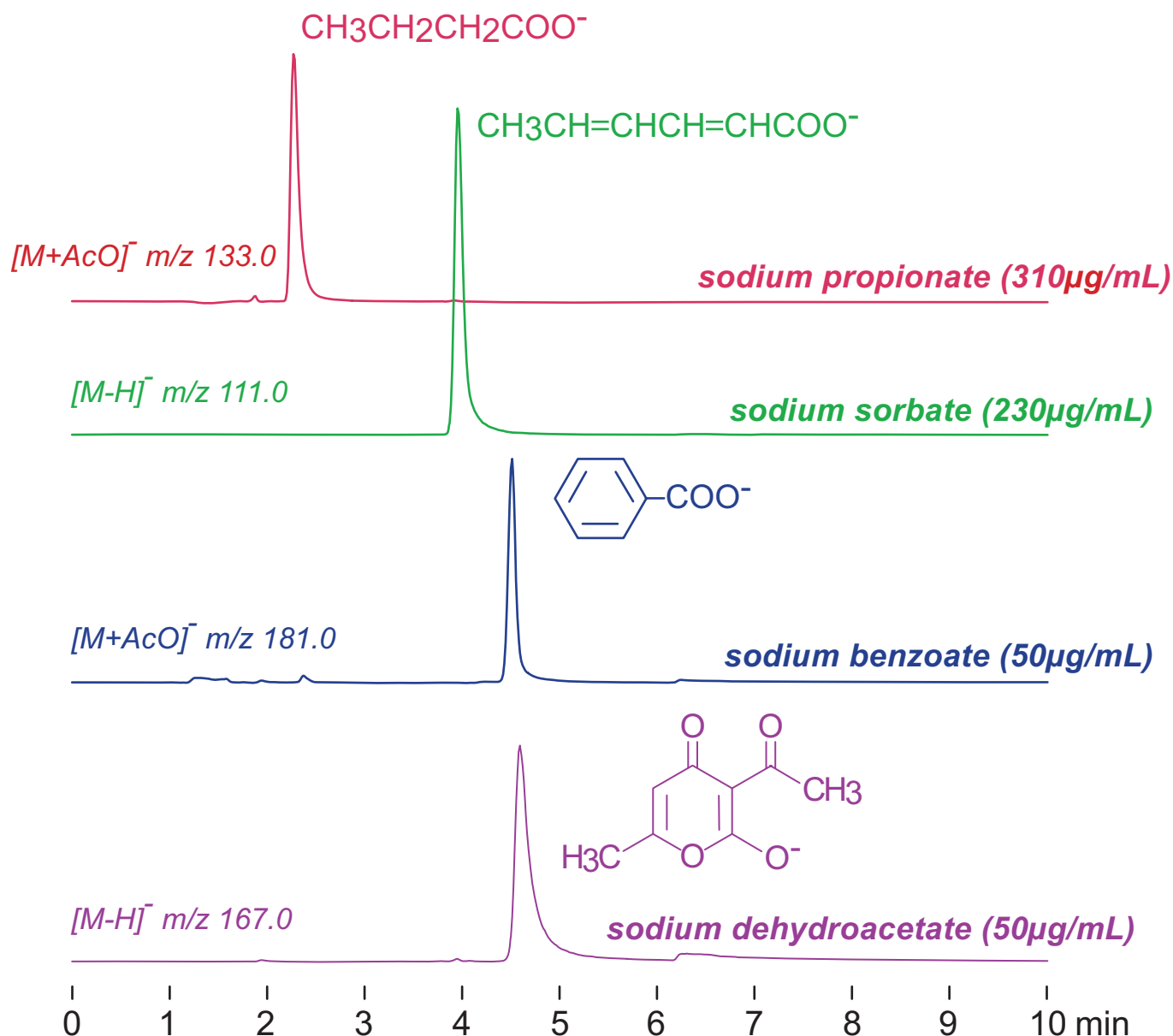

Intrada Organic Acid, 150 x 2 mm

A: acetonitrile / water / acetic acid = 10 / 90 / 0.1

B: acetonitrile / 100mM ammonium acetate = 10 / 90

0%B (0-1min), 0-30%B (1-7min), 100%B (7-10min)

0.2mL/min (12MPa), 37°C, MS (ESI, SIM, Negative)

5µL (0.2N HClO4 in 15% ACN)