






Scherzo SS-C18
 Scherzo SM-C18
 Scherzo SW-C18

100 x 3 mm

Application

LC-MS: Methylimidazole isomers

メチルイミダゾール異性体のLC-MS分析

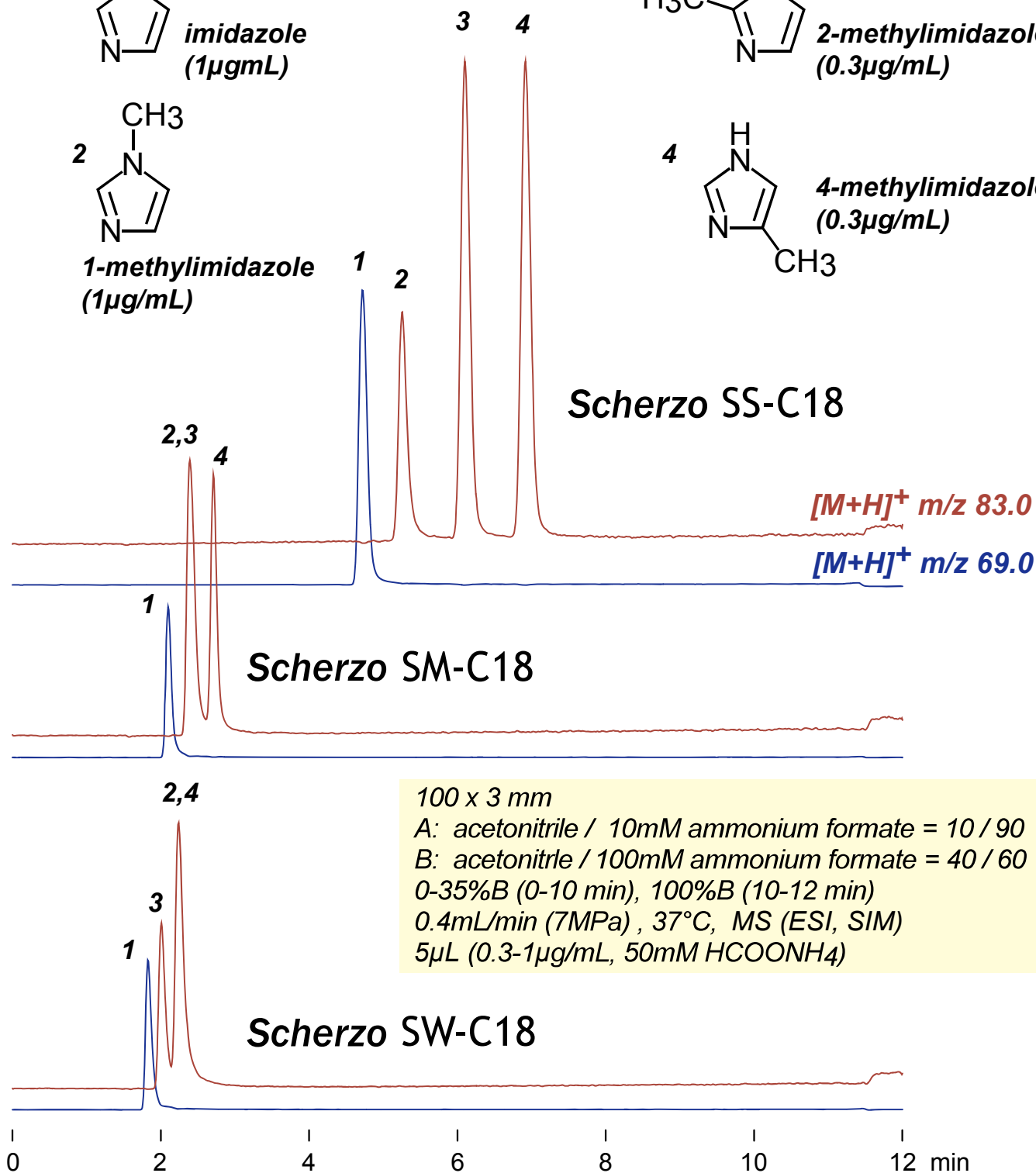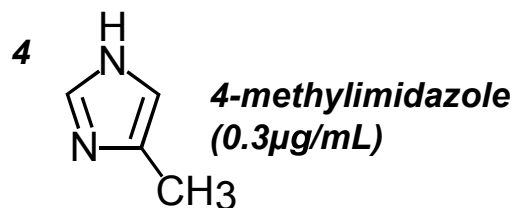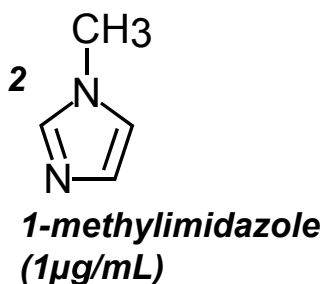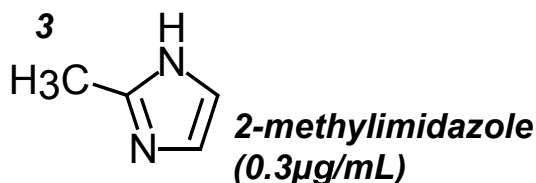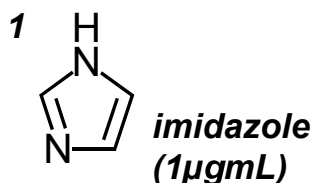

100 x 3 mm
 A: acetonitrile / 10mM ammonium formate = 10 / 90
 B: acetonitrile / 100mM ammonium formate = 40 / 60
 0-35%B (0-10 min), 100%B (10-12 min)
 0.4mL/min (7MPa), 37°C, MS (ESI, SIM)
 5µL (0.3-1µg/mL, 50mM HCOONH₄)