
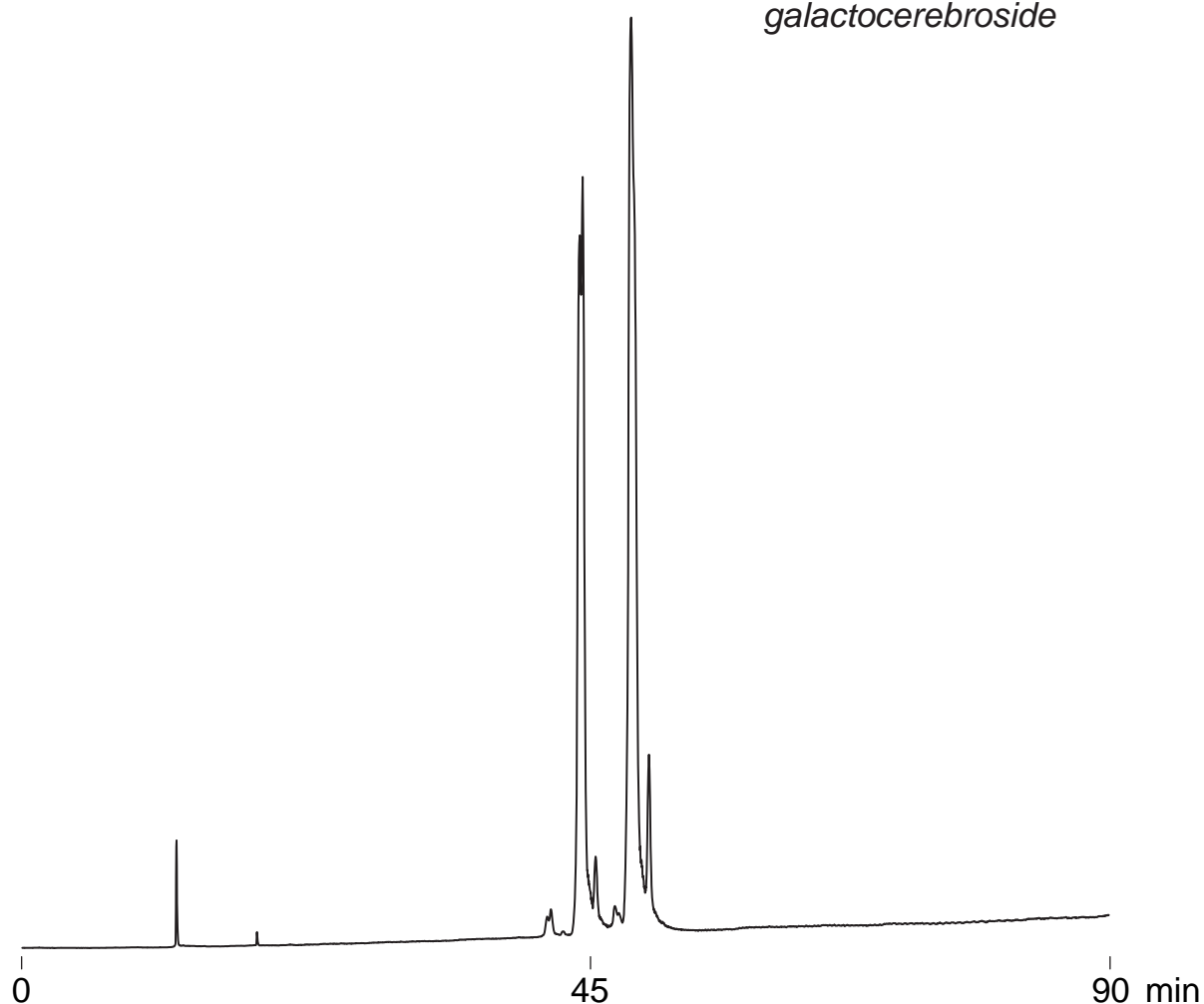


Unison UK-Amino

250 x 4.6 mm

Application

**Galactosylceramide from bovine brain**  
 スフィンゴ糖脂質 (ガラクトセレブロシド, 牛脳由来)

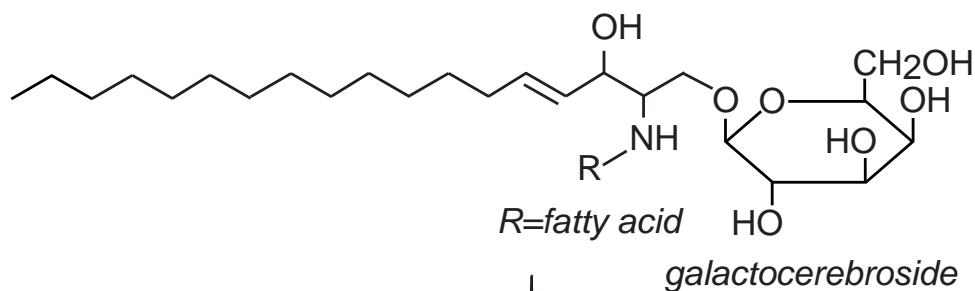

Unison UK-Amino, 250 x 4.6 mm

A: hexane / water / formic acid = 100 / 0.1 / 0.1

B: tetrahydrofuran / methanol / formic acid = 80 / 20 / 0.1

0-100% (0-90min)

1mL/min (5MPa), 37°C, ELSD (spray chamber 50°C, drift chamber 100°C)

5µL (4mg/mL, chloroform / methanol = 2 / 1)