



Cadenza CX-C18

150 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Triazine herbicides

トリアジン系除草剤の LC-MS 分析

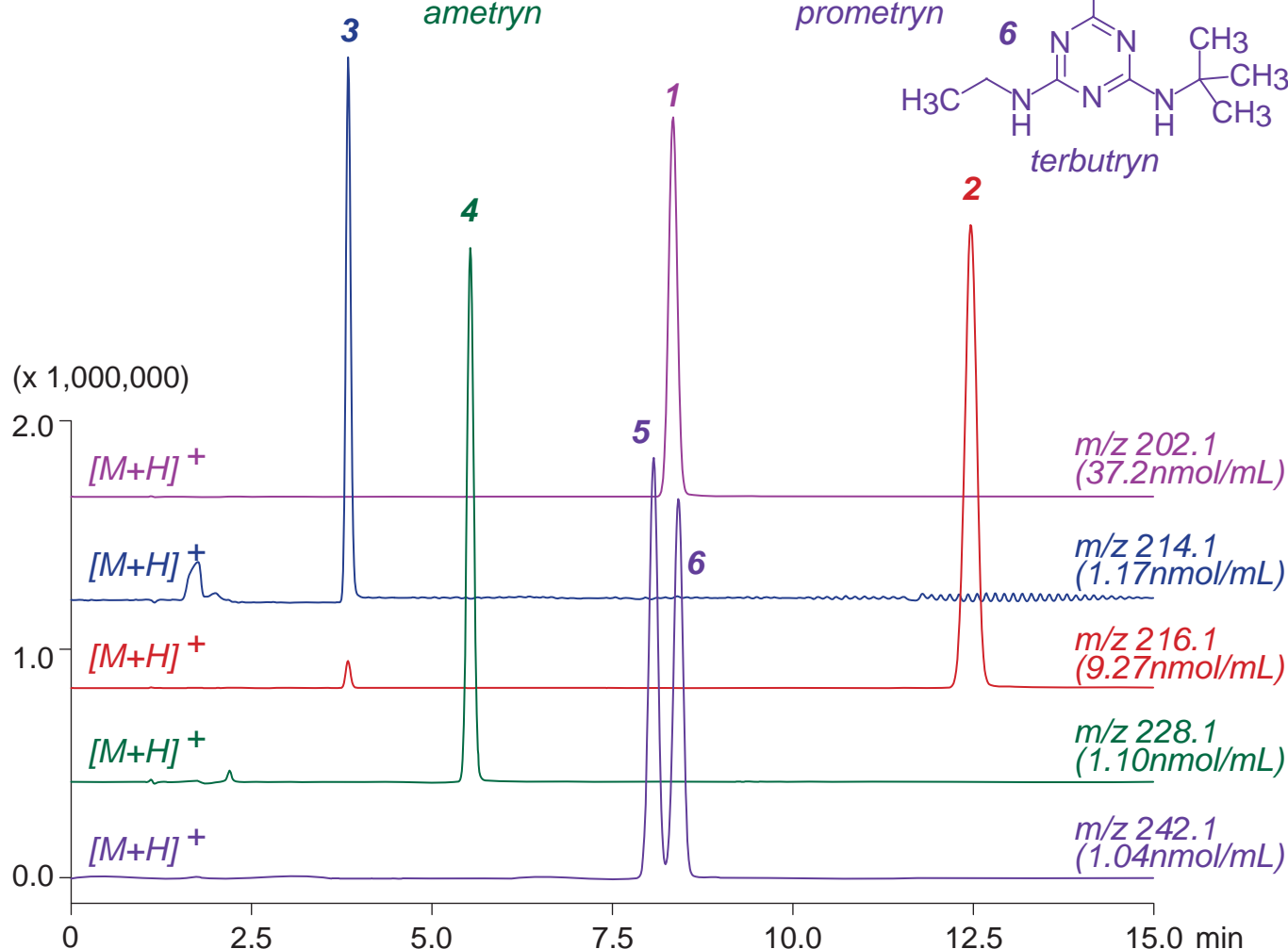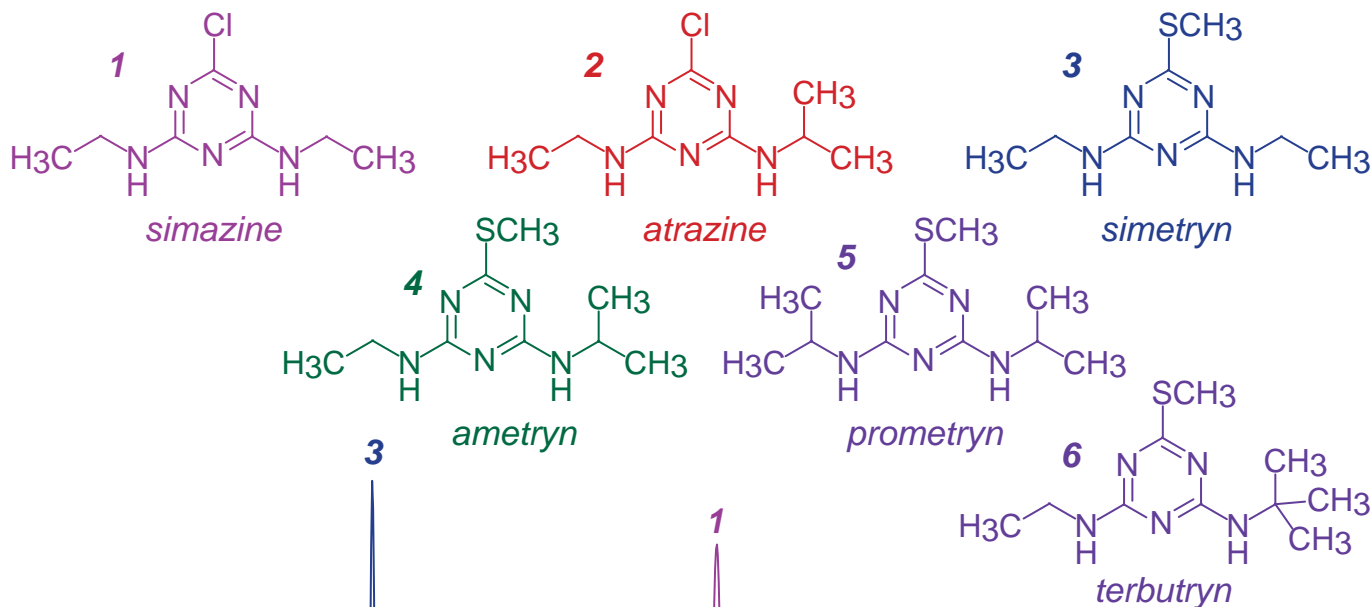

Cadenza CX-C18, 150 x 3 mm
 A: water / formic acid = 100 / 0.1
 B: acetonitrile / formic acid = 100 / 0.1
 30-40%B (0-15min)
 0.4mL/min (12MPa), 37°C, Single Quad. MS (ESI)
 5µL (1.04-37.2nmol/mL, MeOH/ water/ HCOOH=50/ 50 /0.1)