


Presto FF-C18

150 x 4.6 mm

HPLC COLUMN APPLICATION

Polyoxethylene lauryl ether (Nonionic surfactant)

ポリオキシエチレンラウリルエーテル (非イオン性界面活性剤)

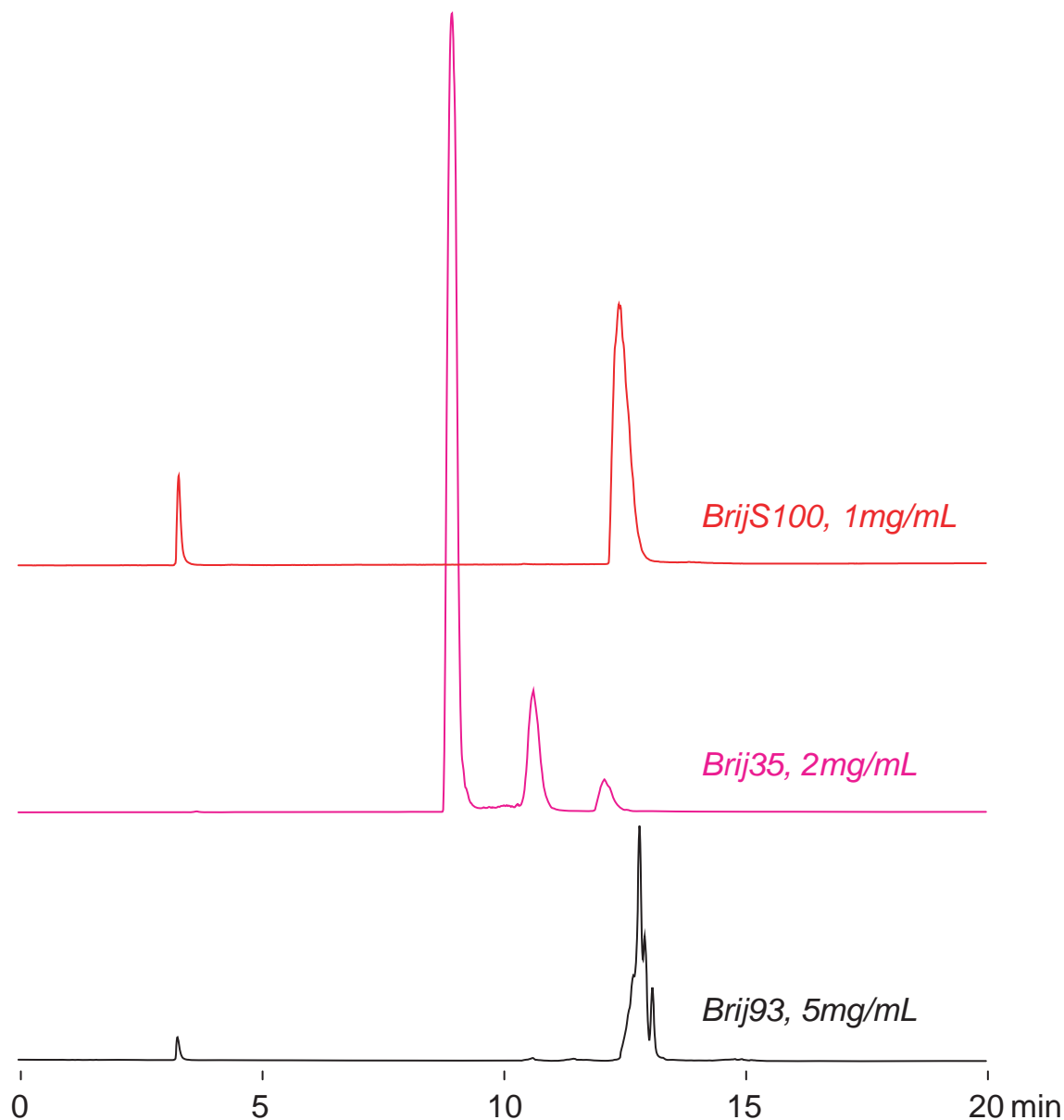

Presto FF-C18, 150 x 4.6 mm

A: water / acetic acid = 100 / 0.5

B: acetonitrile / acetic acid = 100 / 0.5

40-100%B (0-15min), 100%B (15-20min)

0.3mL/min (15MPa), 37°C

ELSD (spray chamber 30°C, drift chamber 60°C)

10µL (1-5mg/mL, acetonitrile / 100mM ammonium formate= 50 / 50)