## Compact mass spectrometer (CMS) analysis of uniformly <sup>13</sup>C-labeled amino acids to identify authentic therapeutic drugs and protect against counterfeits


# 

#### **Overview:**

Cost effective, multidimensional GRAS (generally recognized <u>as safe</u>) counterfeit protection of drugs based on uniformly <sup>13</sup>C labeled amino acids with compact mass spectrometry detection.


#### Introduction:


- Counterfeit drugs are a significant health risk to consumers and an economic burden to the pharmaceutical industry.
- strategy for the unique A novel serialization of genuine drugs can be achieved by addition of uniformly <sup>13</sup>C labelled amino acids (AA) with highly patterned isotopic m/z signals [1,2].
- Isotopic ratio outlier analysis (IROA) utilizing compact mass spectrometry (CMS) is a simple and cost effective analysis approach (Figure 1).

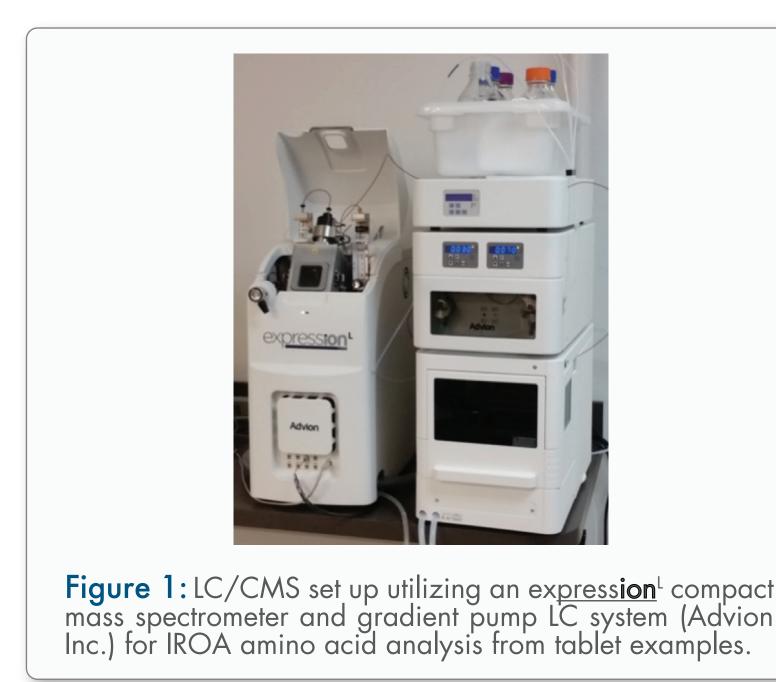

### Methods and Materials:

Uniformly <sup>13</sup>C labeled glutamine and cysteine were selected as GRAS compliant amino acids. They were dissolved in ethanol at one mg/mL and Tic Tacs<sup>®</sup> (Ferrero Inc.) were placed shortly into the mixture, air dried and stored in an Eppendorf tube for further processing (Figure 2). Samples were generated by dissolving the Tic Tac<sup>®</sup> in 0.1 N HCl at 50 mg/mL. 2-200 µL sample was injected onto an Intrada Amino Acid 100 mm x 3 mm column (Imtakt USA, OR) with the following LC gradient settings: 0-3 min at 5 %B; 3-20 min from 5 to 95 %B; 20-22 min from 95 to 5 %B; 22-25 min at 5 %B, with Solvent A (acetonitrile with 0.1 vol% formic acid) and Solvent B (100 mM ammoniumformate) at a total flow rate of 500 µL/min.

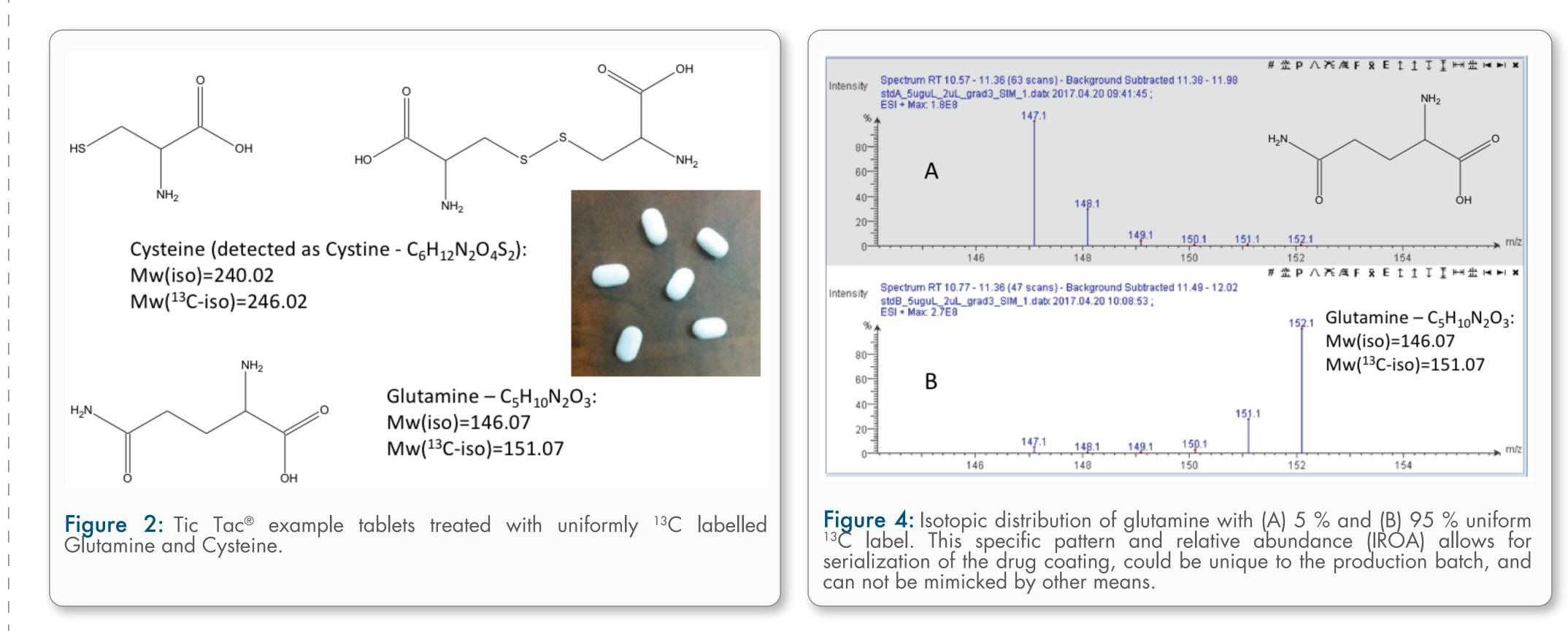


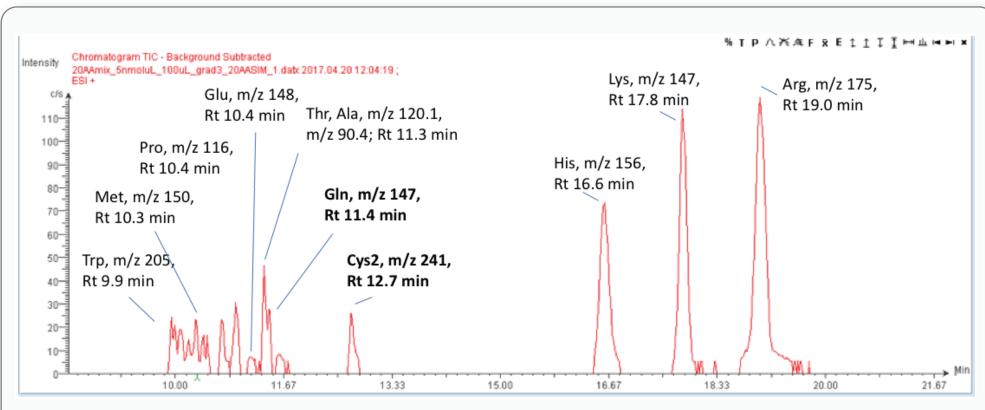

Under these conditions, the amino acid Glutamine eluted at Rt of 11.4 min and Cysteine eluted and was detected as the oxidation product Cystine at Rt 12.7 min (Figure 3) with characteristic <sup>13</sup>C enriched isotopic patterns (example for Glutamine shown in Figure 4). The express**ion<sup>L</sup> CMS** was operated in scan mode from m/z 120 to 250 with a scan time of 300 ms or in SIM mode with m/z 147.1, 148.1, **Figure 3:** Example of a 20 AA mixture (500 nmol each on column) separated chromatogram shows Glutamine and Cysteine (as Cys2 - Cystine) elute at Rt 11.4 and 12.7 min, respectively, (compare [3] for method development and 149.1, 150.1, 151.1, 152.1 for Glutamine and 241.1, [4] for discovery ion chromatogram (DIC) analysis). 242.1, 243.1, 244.1, 245.1, 246.1 and 247.1 for Cysteine (as oxidized Cystine) at dwell times of 50 ms and span of m/z 0.3 each.

 Relative abundance ratios are defined by the uniformity of the label and the mixture ratio of the two batches of e.g. Glutamine and can be used to serialize tablets by various combinations.


• SIM mode shows higher sensitivity of the AA detection and simpler visualization of the isotopic pattern.

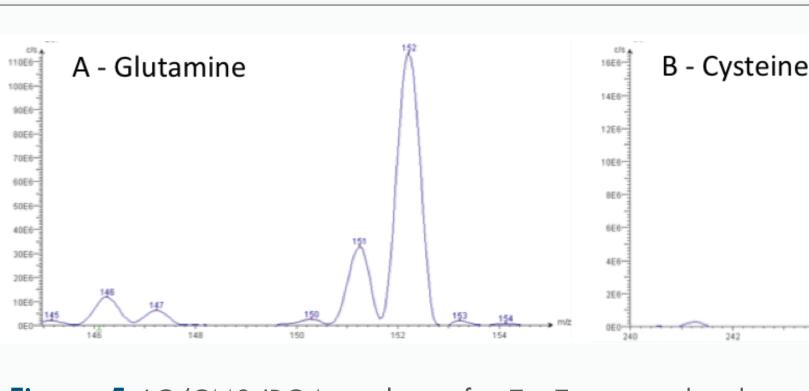



#### **Results:**




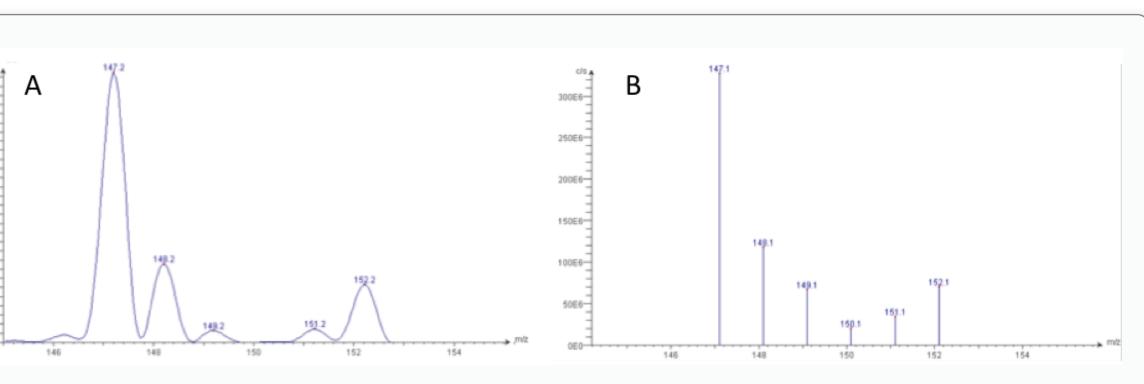

**Figure 6:**LC/CMS IROA analysis of a Tic Tac<sup>®</sup> coated with a mixture of 5% U-<sup>13</sup>C Glutamine and 95% U-<sup>13</sup>C Glutamine in (A) full scan and (B) SIM scan.

% A 90- 70- 50-	A/A+1 ratio: 7 % (theo: 6.4 %)	0066-11 9066-11 9066-11 9066-11 7066-11	A/A+1 ratio: 30 %
40- 30- 20- 10- 146	149.1 149.1 150.1 151.1 152.1 149 150 152 154	40E6- 30E6- 20E6- 10E6- 10E6-	149.1 149.1 149.1 149.1 149.1 152.1 148 150 152 154
<b>igure 7</b> 5% unifo erializati nimicked	: Isotopic distribution pattern of ( orm <sup>13</sup> C label. This specific p on of the drug coating, could by by other means.	Glutamine wit attern and r be unique to	th (A) natural abundance of <sup>13</sup> C and (B) elative abundance (IROA) allows for the production batch, and can not be


- analysis.
- different levels of enrichment.

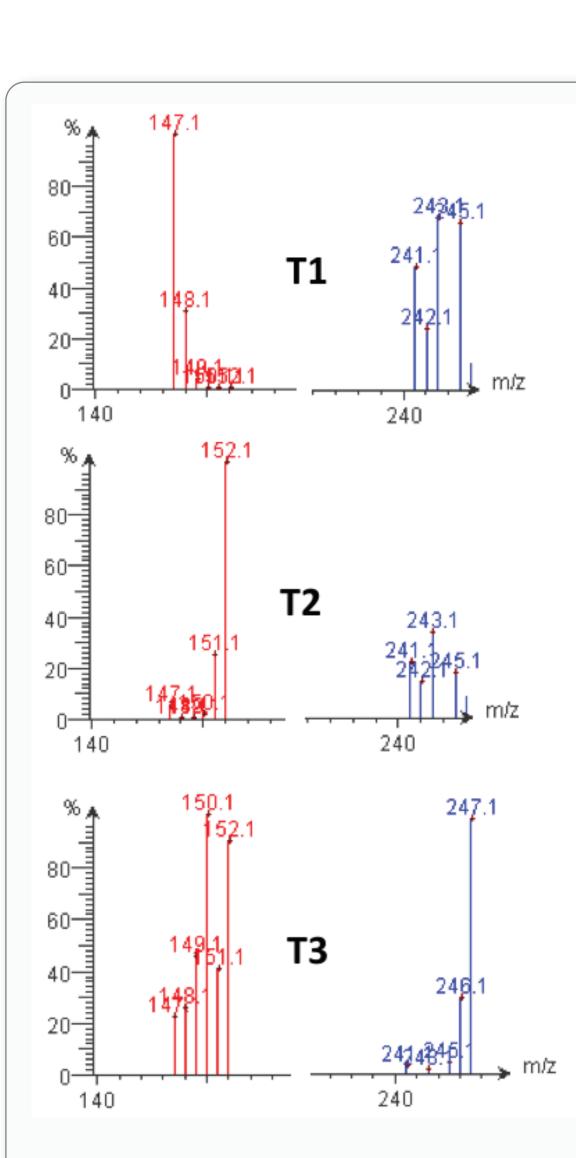

**Figure 8:** Six Tic Tac® Tablets (T1 to T6) analyzed by LC/CMS for 13C-labeled Glutamine (red, left side) and <sup>13</sup>C- labeled Cysteine (blue, right side).

 LC/CMS shows the unique isotopic pattern and enriched A+1 signal, suggesting a 5% level enrichment can reliably be distinguished by IROA

• Example Tic Tac<sup>®</sup> tablets coated with uniformly <sup>13</sup>C labeled Glutamine and/or Cysteine can easily be distinguished from each other by using only two

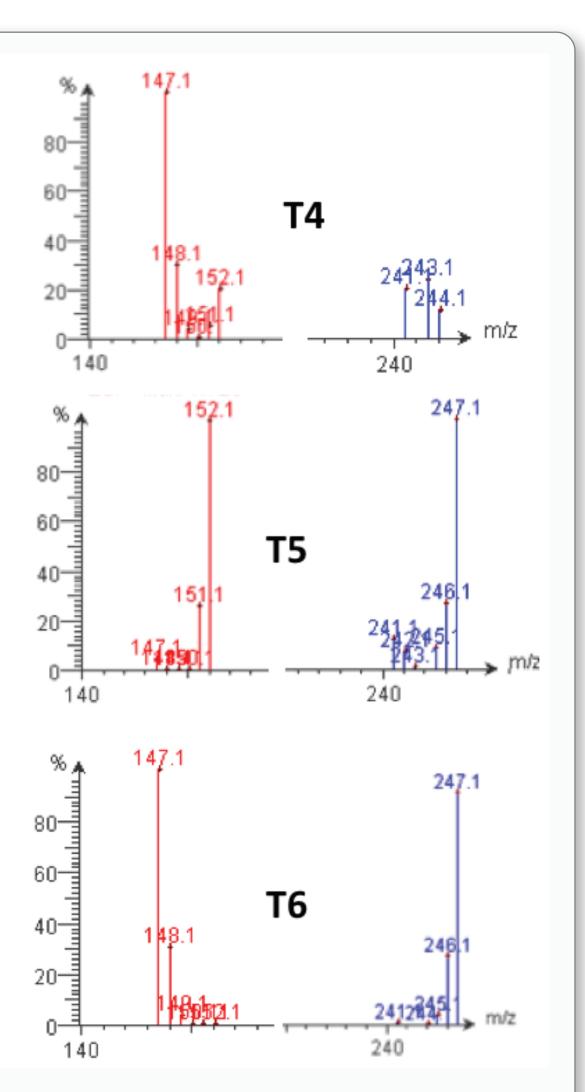

#### Conclusions:

- LC/CMS is a cost-effective IROA analysis approach for uniformly <sup>13</sup>C labeled amino acids.
- Tic Tac<sup>®</sup> example tablets can easily be distinguished using just two different enrichment levels of the same AA or using two different uniformly <sup>13</sup>C AA, resulting labeled in unique relative patterns isotopic and abundances.
- Initial studies of different enrichment levels suggest that LC/CMS can distinguish 5% different enrichment levels which would translate into millions of serialization steps based on 20 amino acids and 20 enrichment levels

#### Literature and Acknowledgements:

<sup>[1]</sup> US Patent 8,969,251 Generation and use of isotopic patterns in mass spectral phenotypic comparison of

<sup>[2]</sup> US Patent 8,536,520 Method for generation and use of isotopic patterns in mass spectral data of simple

<sup>[3]</sup> Eikel D, Prosser SJ and Henion JD: Determination of amino acids using a novel liquid chromatography compact mass spectrometry system (LC/CMS) with mixed-mode separation media. 63rd Meeting of the American Society for Mass Spectrometry (ASMS) 2015, St. Louis, MO, USA

<sup>[4]</sup> Klecha L and Eikel D: Discovery Ion Current – A novel approach to plot total ion current real-time by enhancing signal from newly detected ions. 65th Meeting of the American Society for Mass Spectrometry (ASMS) 2017, Indianapolis, IN, USA

We would like to thank Imtakt USA Corp. for the generous donation of an Intrada Amino Acid LC column.

This poster was presented at the 65th annual conference of the American Society of Mass Spectrometry (ASMS) 2017 in Indianapolis, IN, USA