
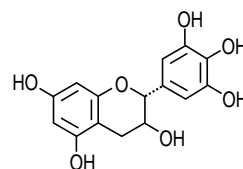


Cadenza CD-C18


75 x 4.6 mm

Application


Catechins

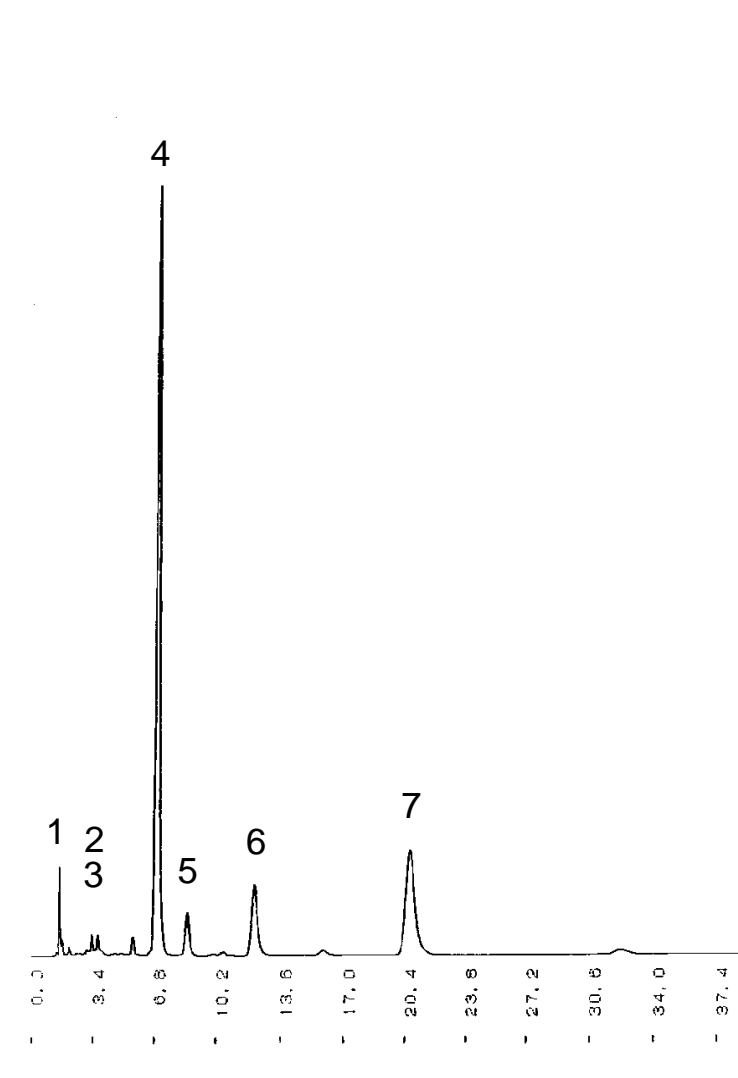


1
galocatechin
GC(+)

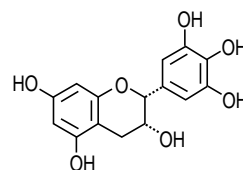


2
epigallocatechin
EGC(-)

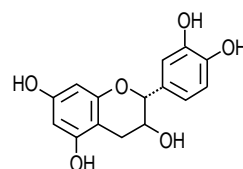


3
catechin
C(+)

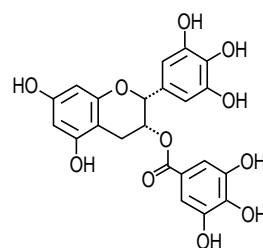


4
epigallocatechin gallate
EGCg(-)

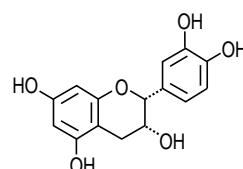

5
epicatechin
EC(-)

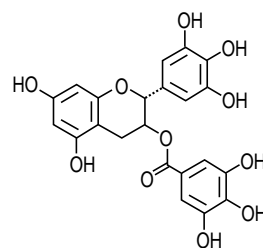

6
galocatechin gallate
GCg(-)

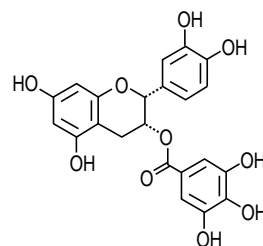

7
epicatechin gallate
ECg(-)

Cadenza CD-C18, 75 x 4.6 mm
0.05 % H₃PO₄ / methanol = 80 / 20
1.0 mL/min, 30 °C
UV at 280 nm
5 uL (0.1%)

Courtesy of Pharma Foods International Co., Ltd.