








Cadenza CD-C18

75 x 4.6 mm

Application

**Food Dyes**

standard

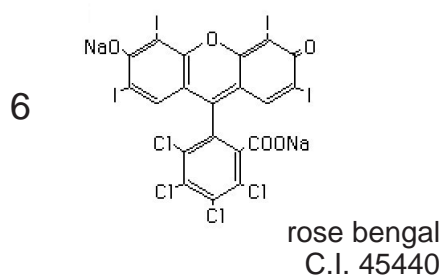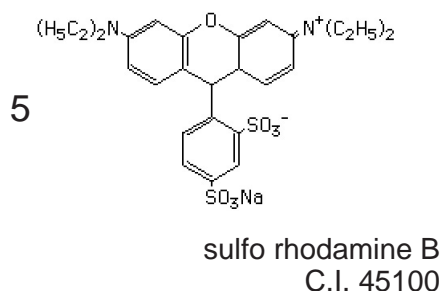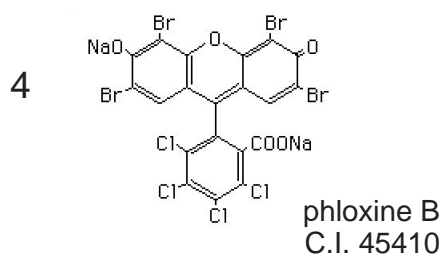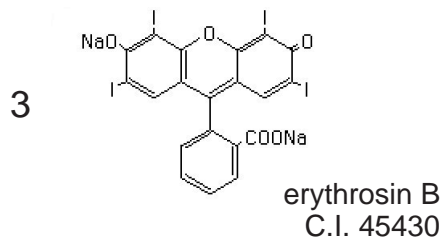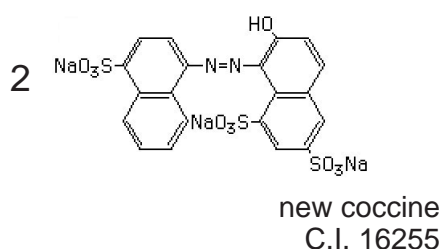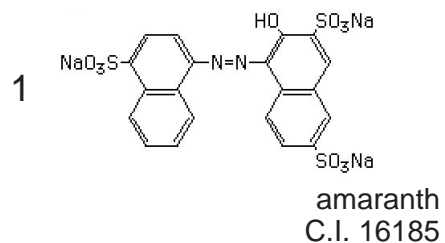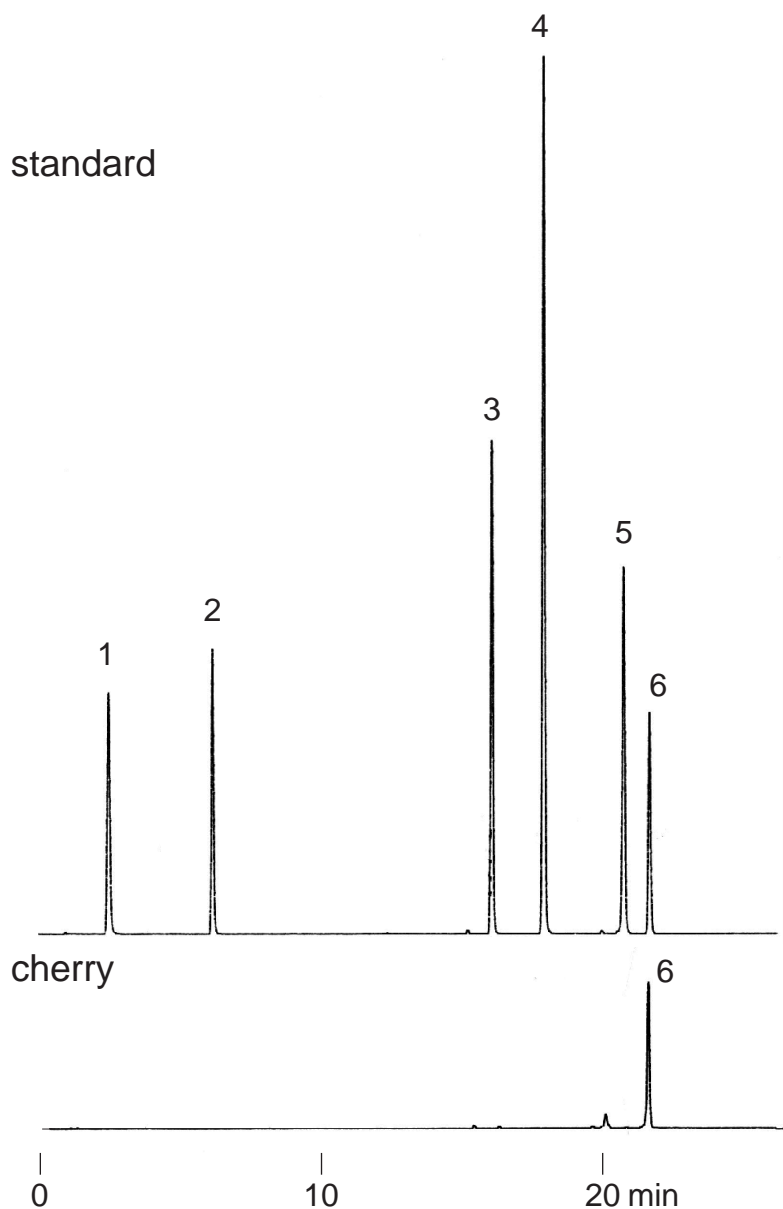

Cadenza CD-C18, 75 x 4.6 mm  
 A : 20mM AcONH4 / methanol = 90 / 10  
 B : 20mM AcONH4 / methanol = 20 / 80  
 0 - 100%B ( 0 - 20 min ), 100%B ( 20 - 23 min )  
 1 mL/min, 30°C, VIS at 530 nm, 0.32 AUFS

Courtesy of S.Kuroda, Bio College Kyoto