








Cadenza CD-C18 Unison UK-Phenyl
 Cadenza CL-C18 Unison UK-C8
 Unison UK-C18 Cadenza HS-C18

75 x 4.6 mm

Application

Polyphenol, Berry

ベリー類のポリフェノール

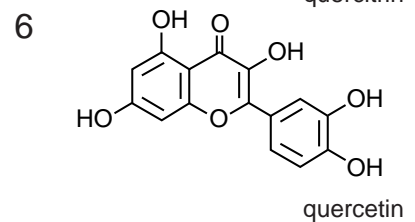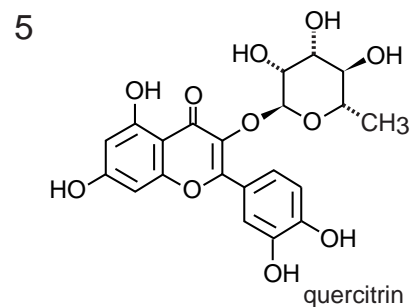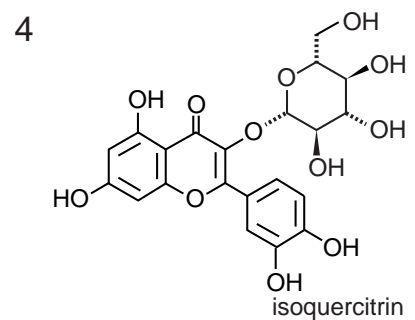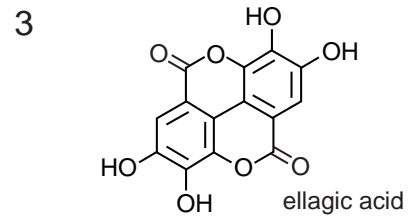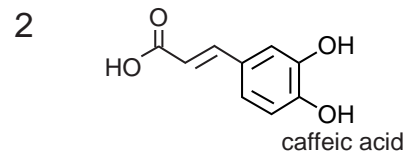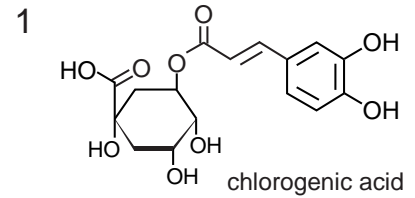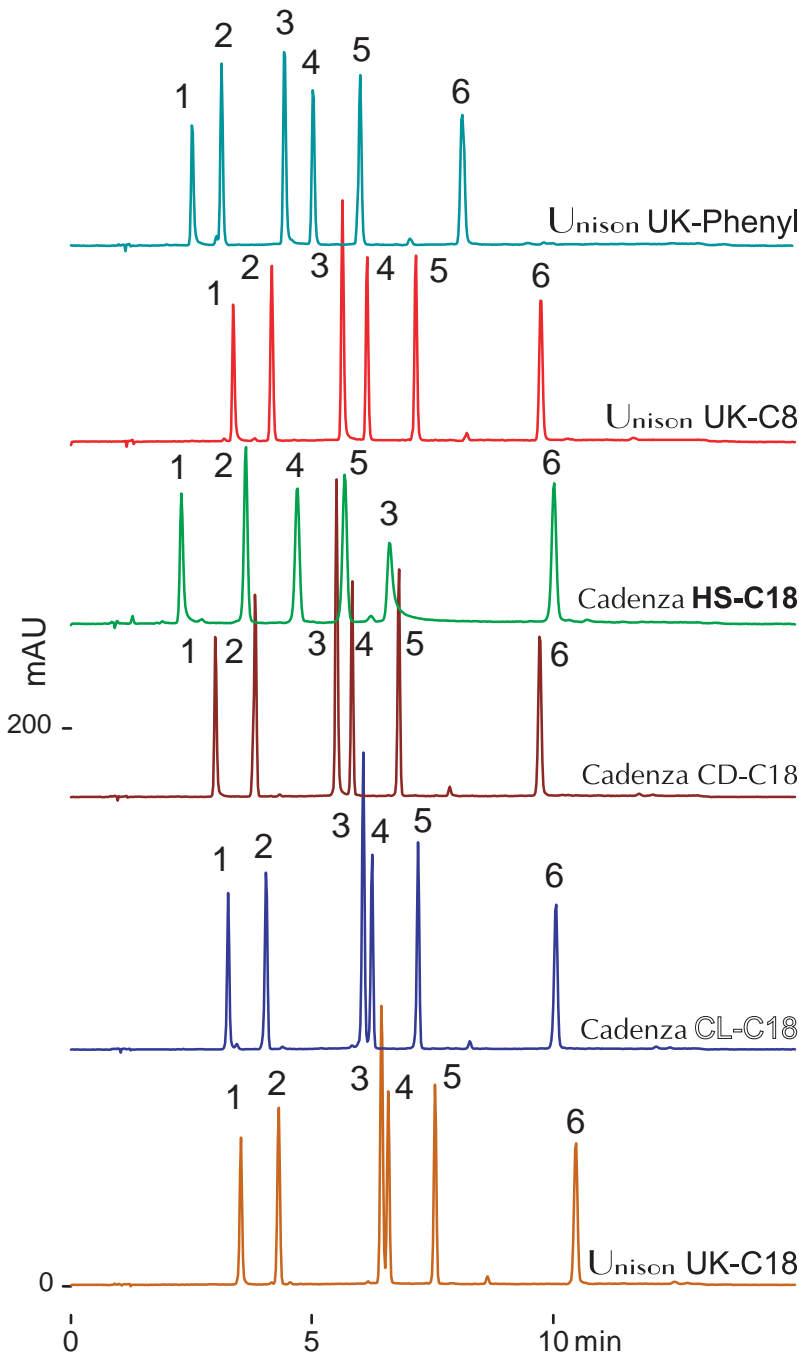

75 x 4.6 mm

A: water /trifluoroacetic acid = 100 /0.1

B: acetonitrile /trifluoroacetic acid = 100 /0.1

10-40%B (0-12min), 1 mL/min

37 deg.C, 260 nm, 2uL (0.1-0.8ug)