


Unison UK-Amino

150 x 2 mm


Application

## Dehydroascorbic Acid and Related Compounds


デヒドロアスコルビン酸と関連化合物

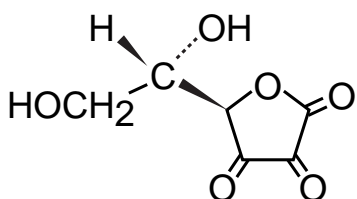


**1** dehydroascorbic acid

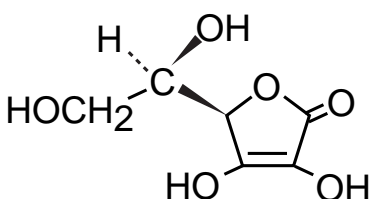

**2** isoascorbic acid  
(erythorbic acid)

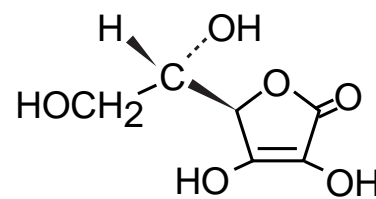


**3** L-ascorbic acid  
(Vitamin C)

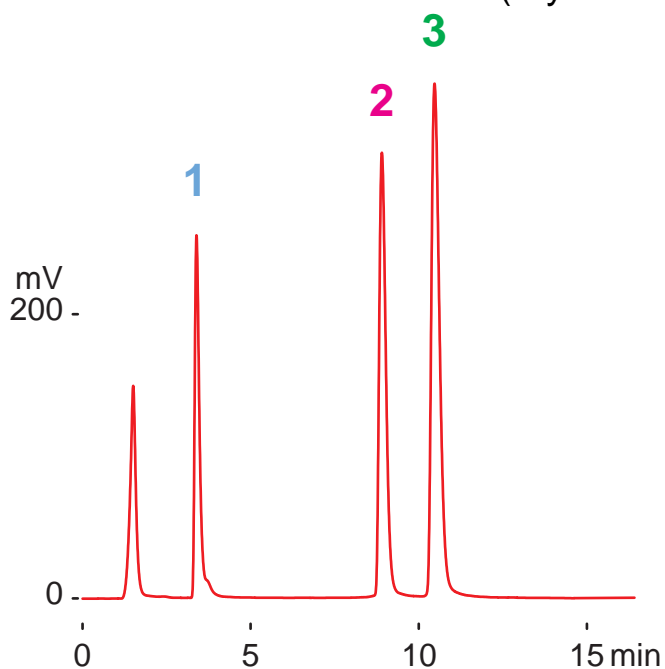

**normal phase**

Unison UK-Amino  
150 x 2 mm  
acetonitrile /water /acetic acid =  
80 / 20 / 5  
0.2mL/min (5MPa),  
37deg.C, ELSD,  
1uL(2-4ug)

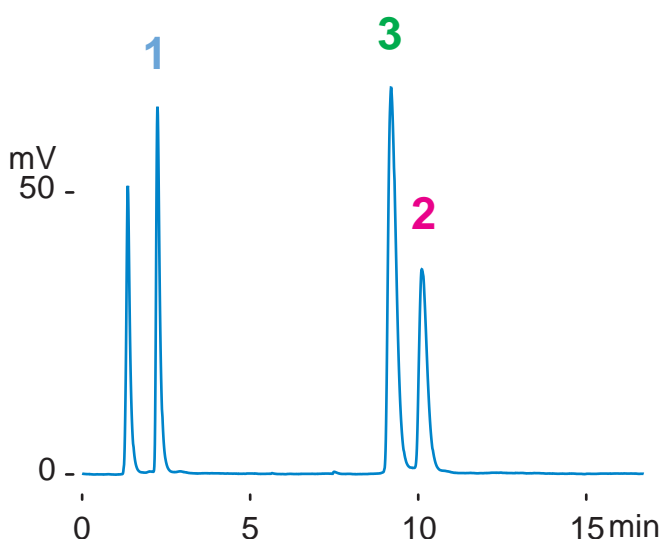

**anion exchange**

Unison UK-Amino  
150 x 2 mm  
water /acetic acid = 100 /0.5  
0.2mL/min (9MPa),  
37deg.C, ELSD,  
1uL(2-4ug)