



Unison UK-Amino  
Cadenza CL-C18

250 x 4.6 mm

Application


### Surfactant (Igepal)

界面活性剤(イゲパール)

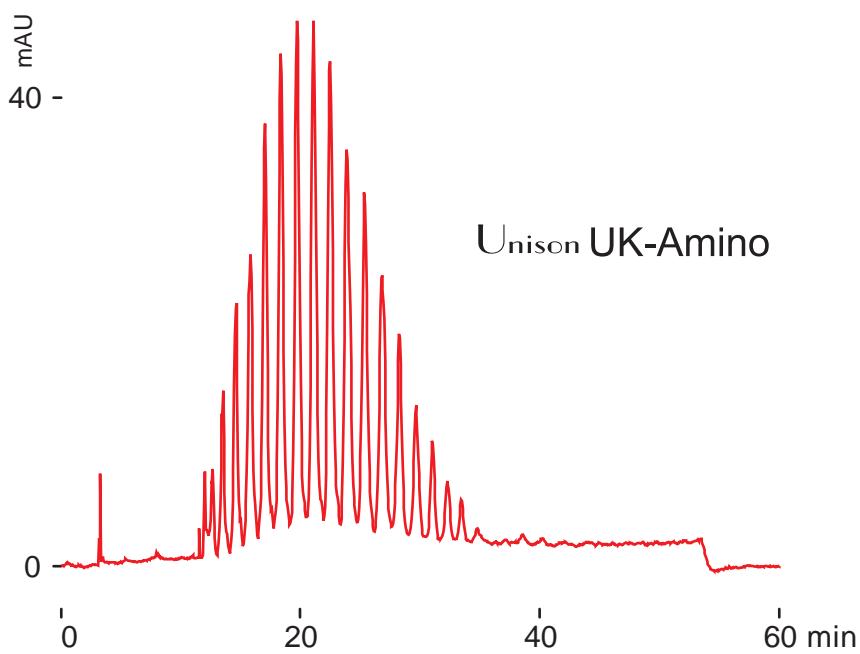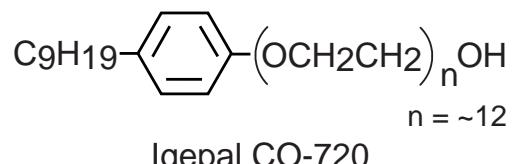

normal phase

Unison UK-Amino  
250 x 4.6 mm,  
A: hexane, B: ethanol  
0.2-10% B(0-50min)  
1 mL/min 37 deg.C  
260 nm, 5 uL (10% solution)

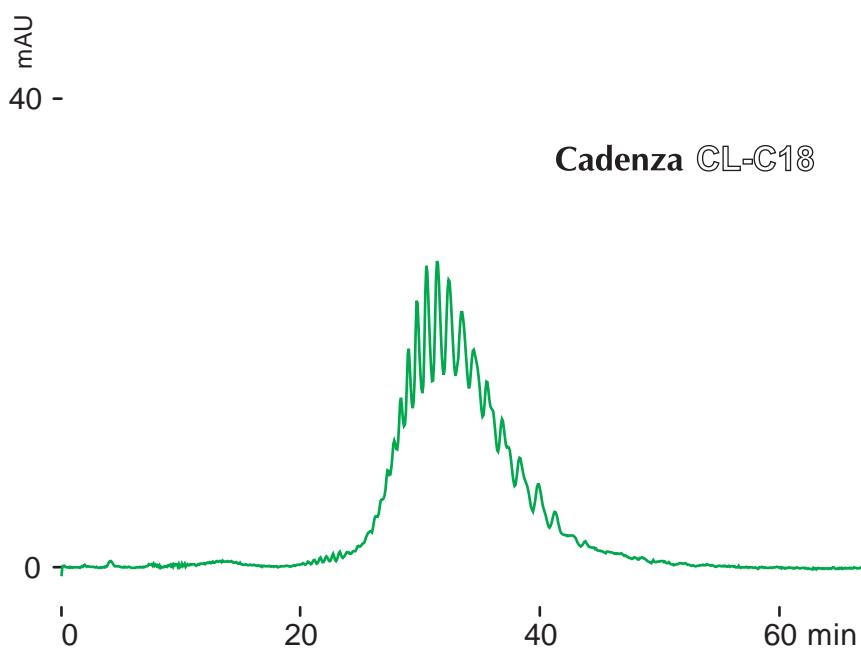

reversed phase

Cadenza CL-C18  
250 x 4.6 mm  
water / acetonitrile = 40 / 60  
0.8 mL/min, 37 deg.C  
260 nm, 5 uL (10% solution)