

Scherzo SM-C18


150 x 3 mm

Application

### Glyphosate

グリホサート


mV  
100 -

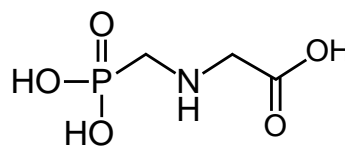

glyphosate

50 -

20 mM ammonium formate / acetonitrile  
= 90 / 10  
9MPa

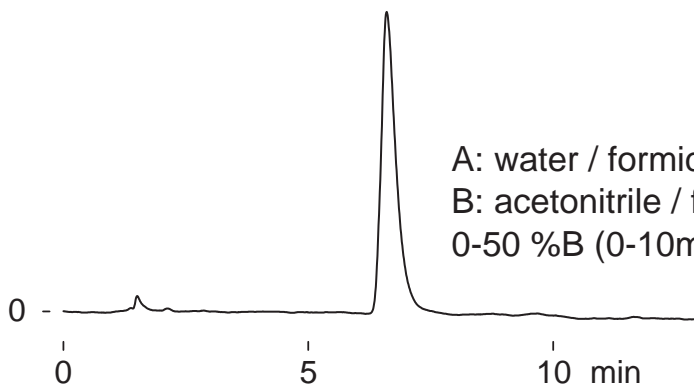

Scherzo SM-C18, 150 x 3 mm

0.4 mL/min, 37 deg.C,ELSD (spray chamber 50 deg.C, drift tube 100 deg.C)

5 uL (5 ug)

Courtesy of Takeo Sakuma, Ph.D, MDS Analytical Technologies (SCIEX), Canada