








Unison UK-Amino

250 x 3 mm

Application

Unsaturated Chondro-Disaccharides

不飽和コンドロニ糖

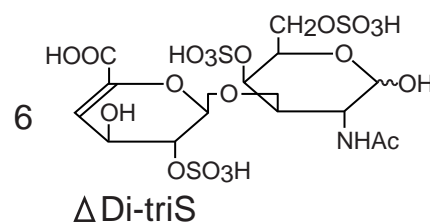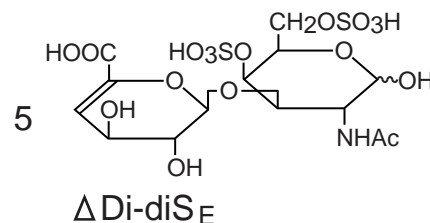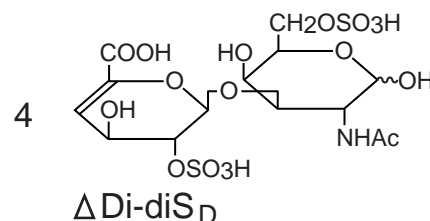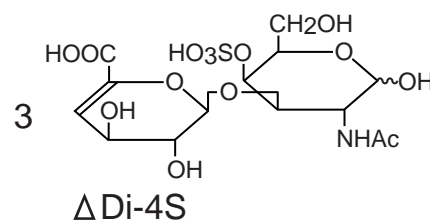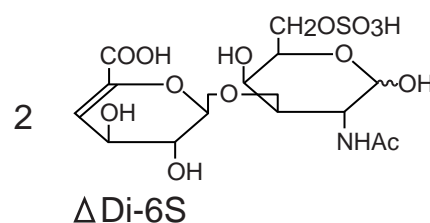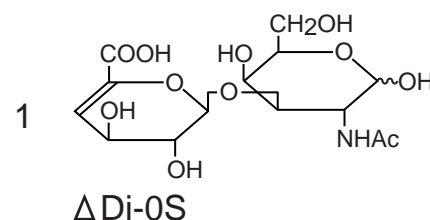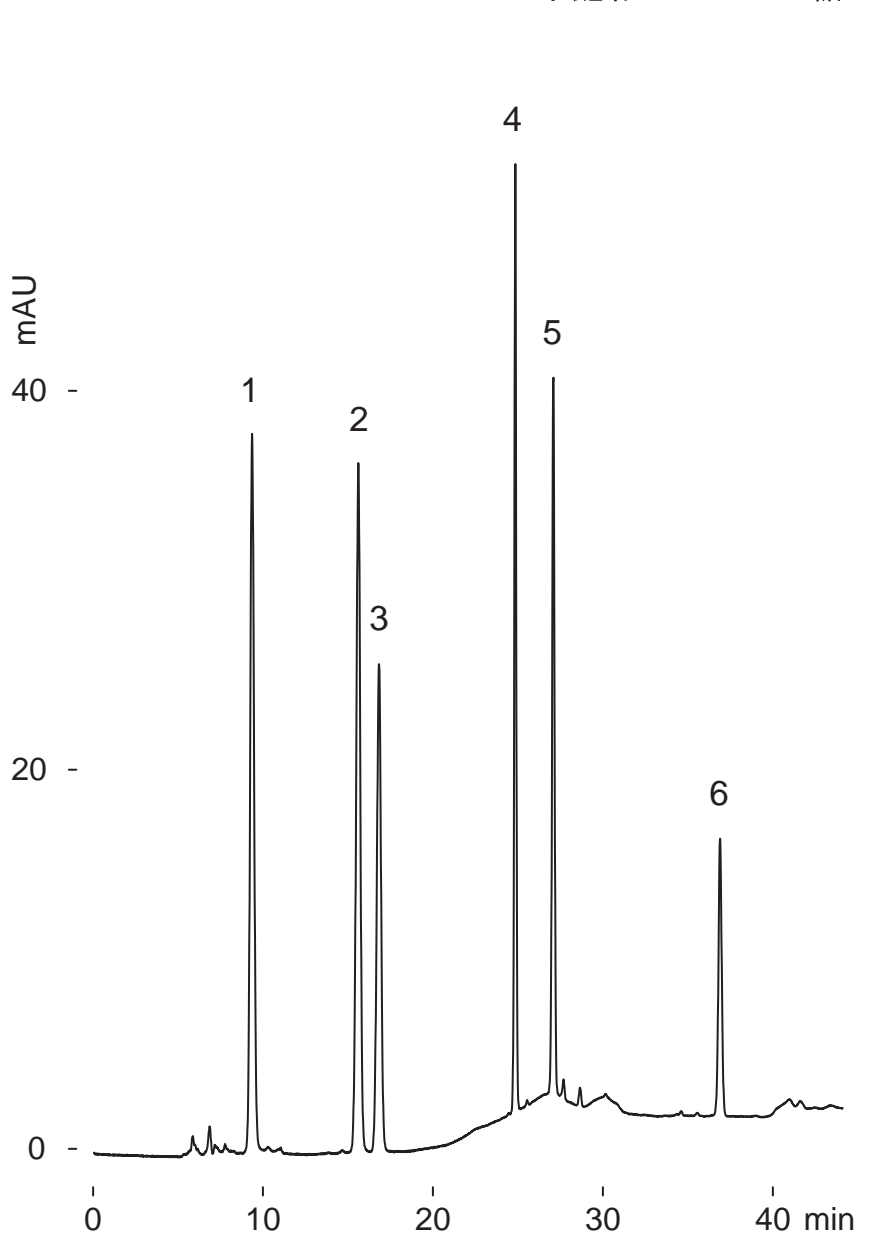

Unison UK-Amino, 250 x 3 mm
 A: 5 mM (NH₄)₂HPO₄ / acetonitrile = 40 / 60
 B: 100 mM (NH₄)₂HPO₄
 2-3 %B (0-10 min), 3-40 %B (10-40 min)
 0.4 mL/min (5 MPa), 50 deg.C, 232 nm
 4 uL (0.26-0.75 ug, 100 mM (NH₄)₂HPO₄)