


Presto FF-C18

250 x 4.6 mm

Application

## Gradient mixer volume affecting HSA separation

ヒト血清アルブミンの分離に与えるグラジエントミキサー容量の影響

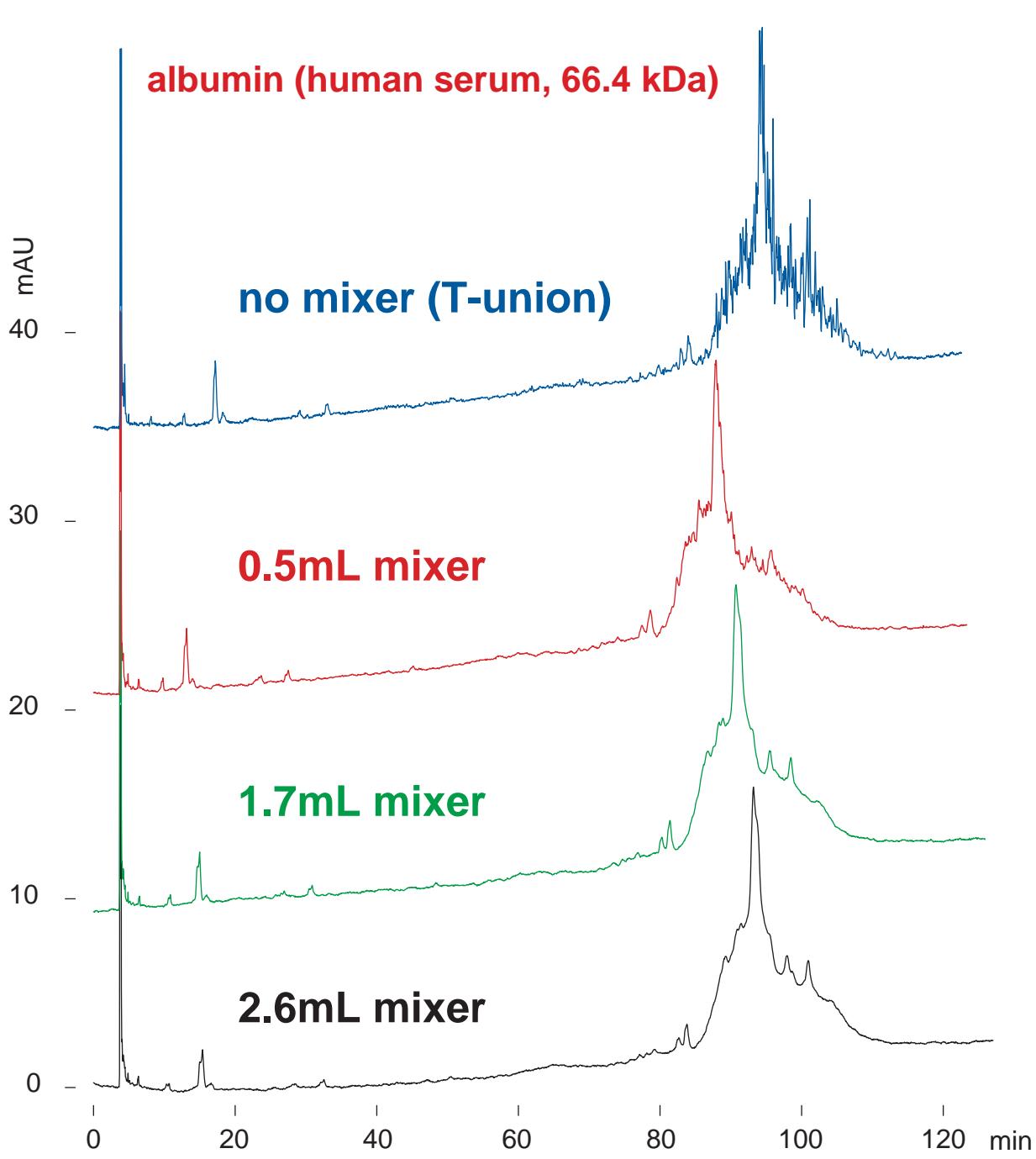

Presto FF-C18, 250 x 4.6 mm

A: water / trifluoroacetic acid = 100 / 0.1

B: acetonitrile / trifluoroacetic acid = 100 / 0.1

30-40 %B (0-120 min)

0.4 mL/min (18 MPa), 80 deg.C, 220 nm, 1.2 uL (12 ug)