


Presto FF-C18

75 x 4.6 mm

Application

Protein separation on non-porous ODS column (80°C, formic acid)

ノンポラスODSカラムによる蛋白質の分離 (80)

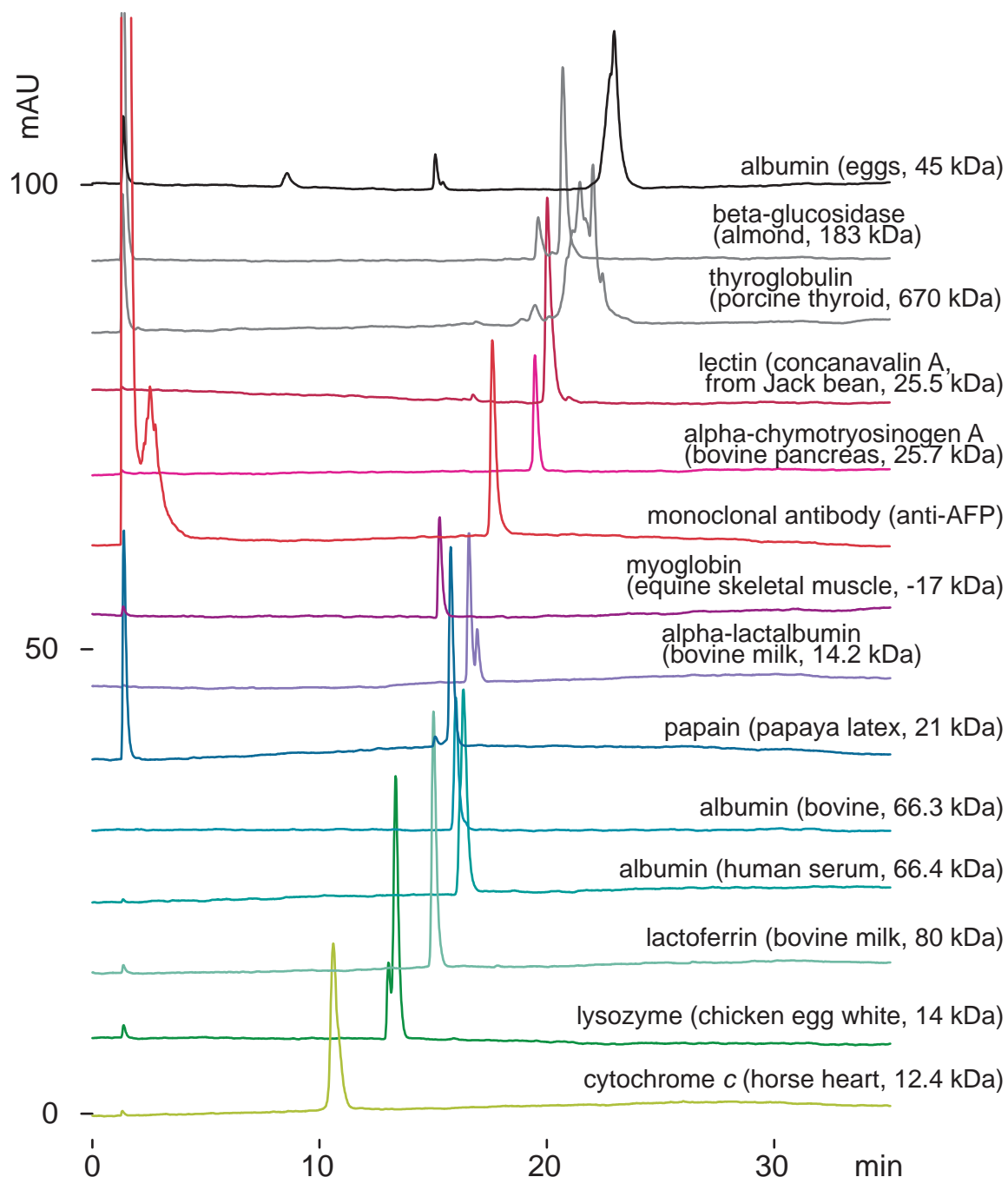

Presto FF-C18, 75 x 4.6 mm

A: water / formic acid = 100 / 0.5, B: acetonitrile / formic acid = 100 / 0.5
10-50 %B (0-30min), 0.4 mL/min(5MPa)

80 deg.C, ELSD(spray chamber 35, drift tube 80 deg.C), 1-14 uL(2-14ug)