



Unison UK-C18 HT

75 x 2 mm

Application

Tadalafil, Sildenafil and Vardenafil

タダラフィル, シルデナフィル, バルデナフィル

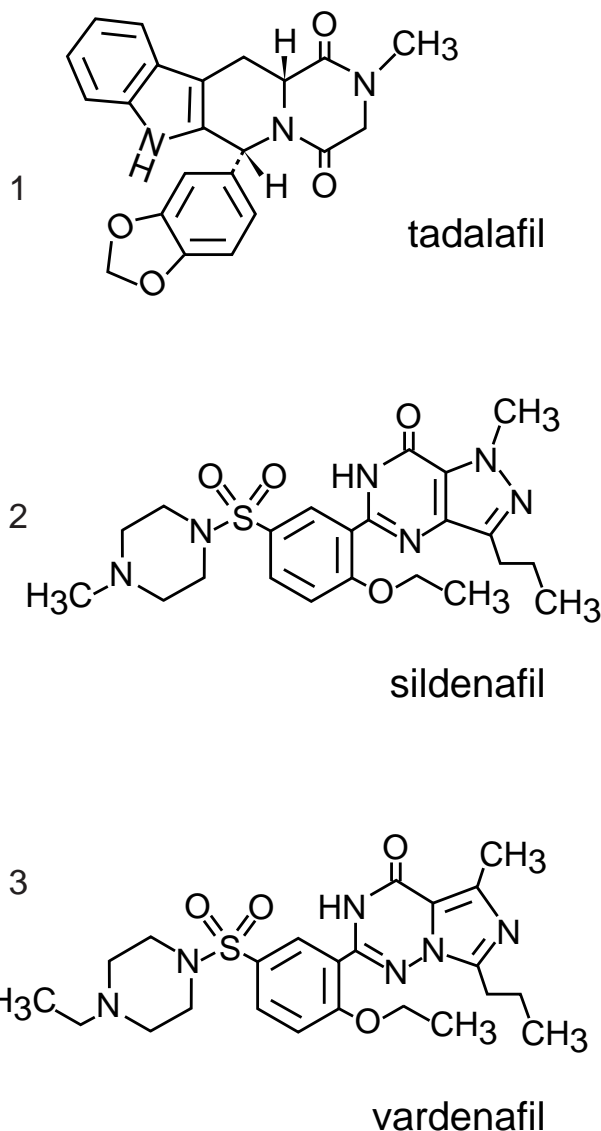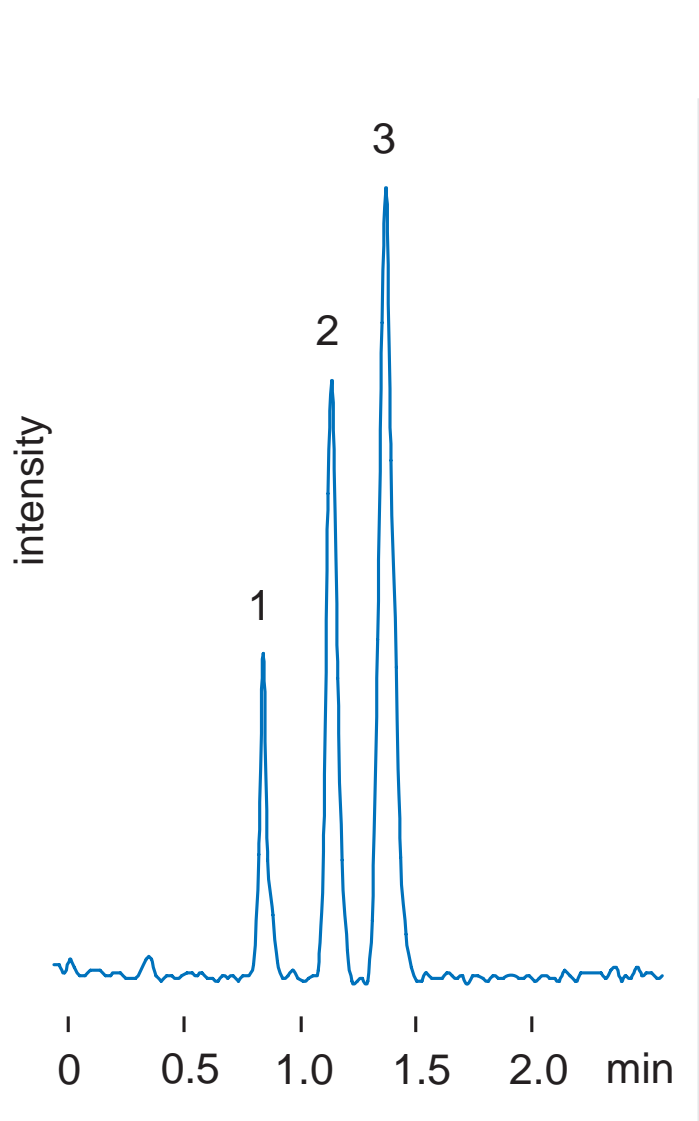

Unison UK-C18 HT, 75 x 2 mm

33mM ammonium acetate / methanol = 30 / 70

0.5mL/min (19-20MPa), 40 deg.C

2uL (2ng-10ng)

LC-MS/MS instrument : SHIMADZU LC-20A + Thermo Fisher SCIENTIFIC LXQ
(ESI, positive)

Courtesy of Kyoto Prefecture Police, Forensic Science Laboratory, JAPAN