

Cadenza CD-C18 HT



100 x 4.6 mm

Application


Comparison between 3um and sub-2um ODS columns

3um と sub-2um ODSカラムの比較

3um Cadenza CD-C18 HT, 100 x 4.6 mm

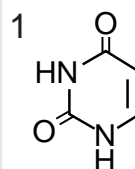
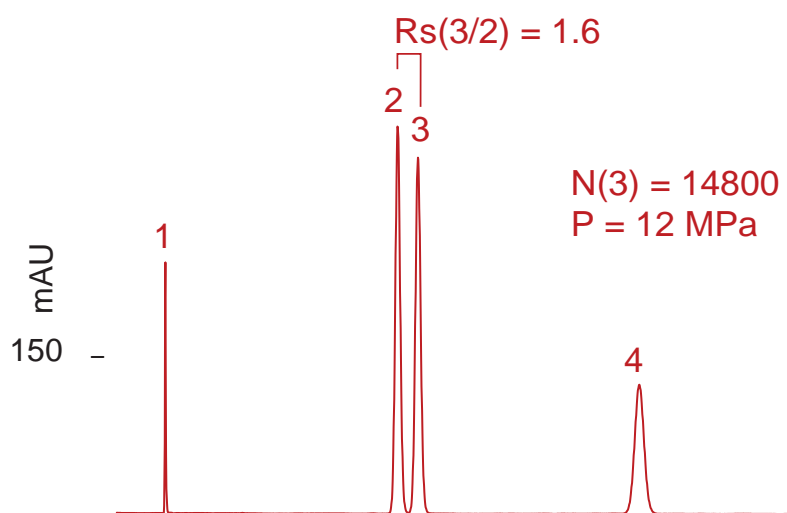


uracil

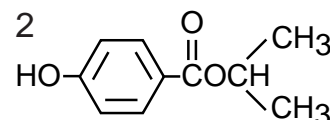


isopropylparaben

1.7um ODS, 100 x 4.6 mm

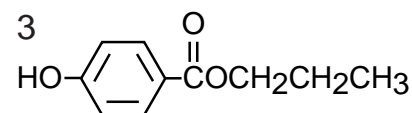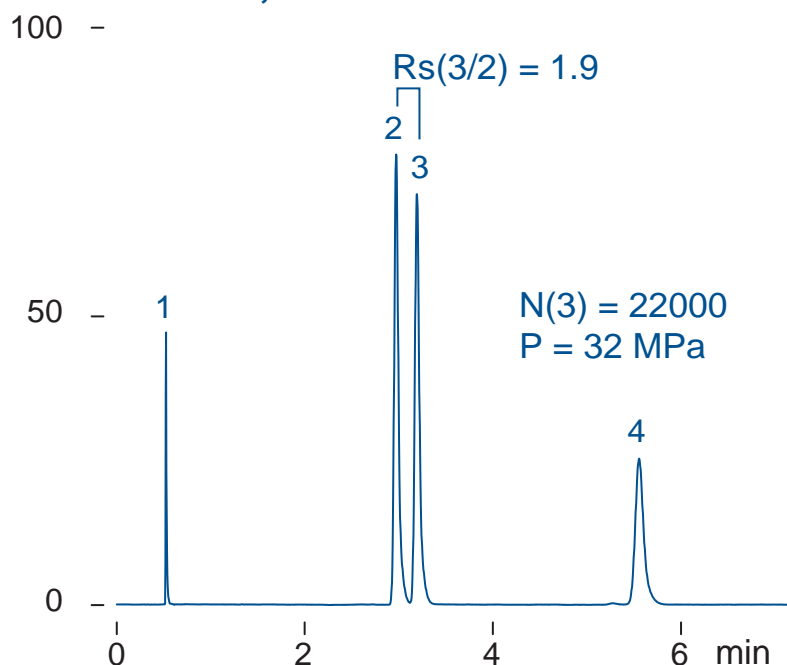

propylparaben

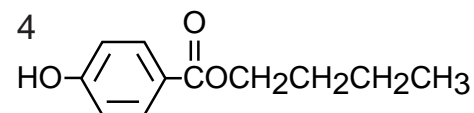

butylparaben

100 x 4.6 mm

water / acetonitrile / acetic acid = 60 / 40 / 0.1

1.8 mL/min, 37 deg.C, 260 nm, 1 uL (0.02- 0.16ug)