

**Scherzo SM-C18**  
**Unison UK-C18**

50 x 3 mm


Application

**Inositol hexaphosphate ( phytic acid )**

イノシトール6リン酸 ( フィチン酸 )

mV  
400 -

200 -

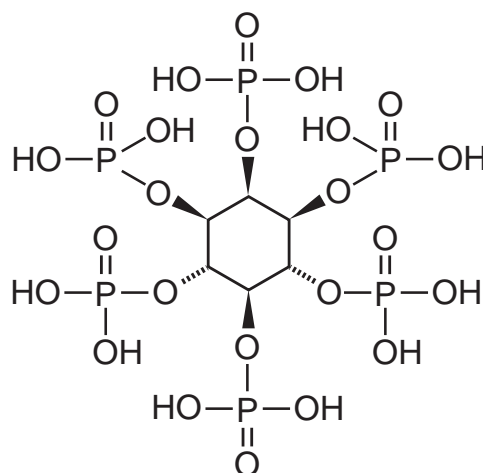

inositol hexaphosphate  
( phytic acid )

**Scherzo SM-C18**



**Unison UK-C18**

50 x 3 mm

A: water / methanol / HCOOH = 80 / 20 / 0.05

B: 50 mM HCOONH<sub>4</sub> / 50 mM NH<sub>4</sub>OH / methanol = 40 / 40 / 20

0-100 %B (0-8 min), 100 %B (8-14 min), 0 %B (14-20 min)

0.4 mL/min (5-6 MPa), 37 deg.C

ELSD (spray chamber 50 deg.C, drift tube 100 deg.C), 1 uL (0.5 %, water)