



Cadenza CW-C18
Unison UK-C18
Unison UK-C8

150 x 3 mm


Application

Coenzyme A (CoA) and acetyl coenzyme A (Acetyl-CoA)

補酵素： コエンザイム A と アセチルコエンザイム A

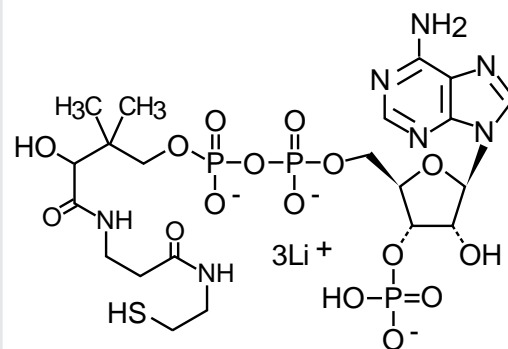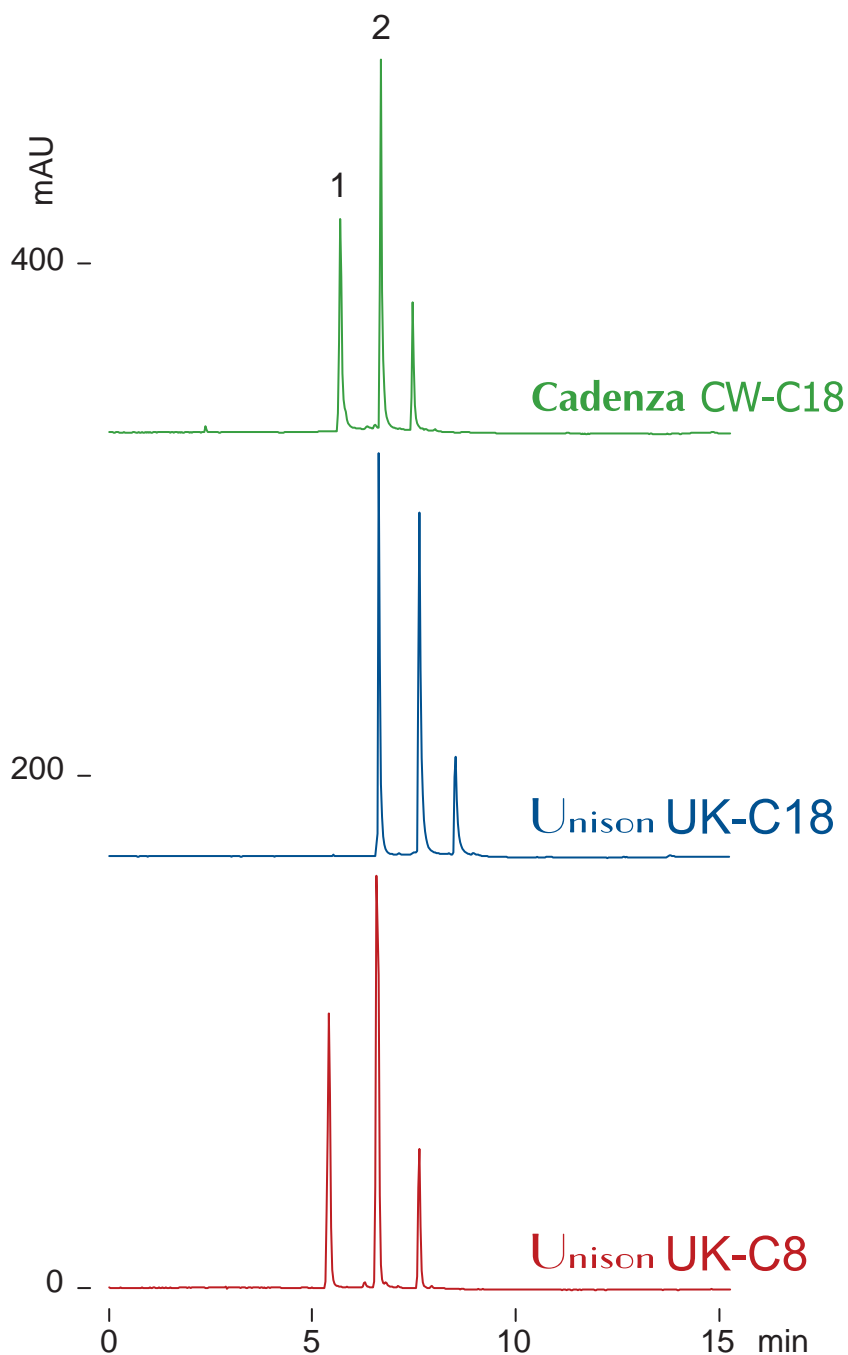

1 coenzyme A (CoA)

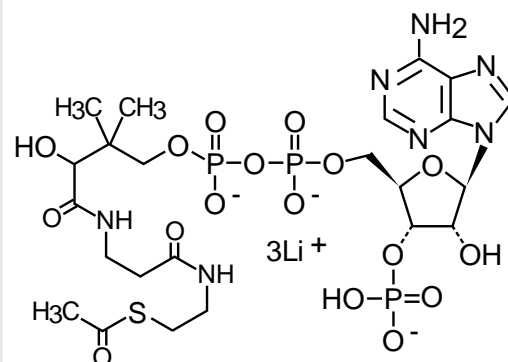

2 acetyl coenzyme A (Acetyl-CoA)

150 x 3 mm

A: 50 mM ammonium formate

B: 50 mM ammonium formate / acetonitrile = 50 / 50

0-60 %B (0-15 min)

0.4 mL/min (8-10 MPa), 37 deg.C, 260 nm, 2 uL (0.3-0.7 ug, 0.5% HCOOH aq.)