

Intrada Amino Acid

100 x 3 mm


Application

LC-MS: 20 standard amino acids (recommended)

20種アミノ酸(標準試料)のLC-MS分析(推奨メソッド)

Recommended method for amino acids analysis

100 x 3 mm

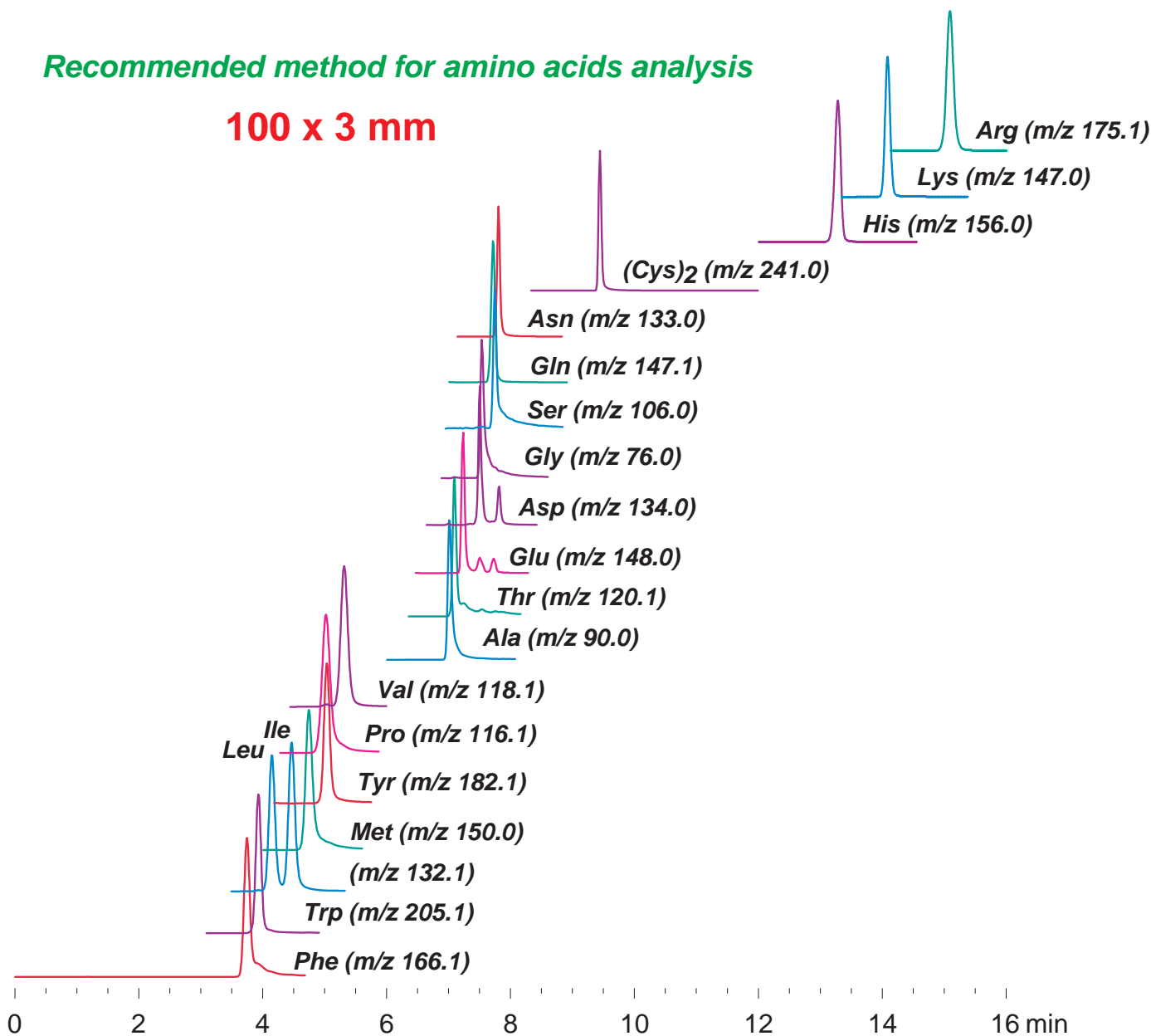

Intrada Amino Acid, 100 x 3 mm

A: acetonitrile / formic acid = 100 / 0.3

B: acetonitrile / 100mM ammonium formate = 20 / 80

20 %B (0-4 min), 20-100 %B (4-14 min), 100 %B (14-16 min)

0.6 mL/min (5MPa), 37°C, 5 uL (1000nmol/mL, 0.1N HCl)

Single Quad. MS: ESI (SIM, positive)