




**Scherzo SS-C18**

100 x 2 mm

Application

**LC-QTOF/MS: Acyl carnitine**

アシルカルニチンのLC-QTOF/MS分析

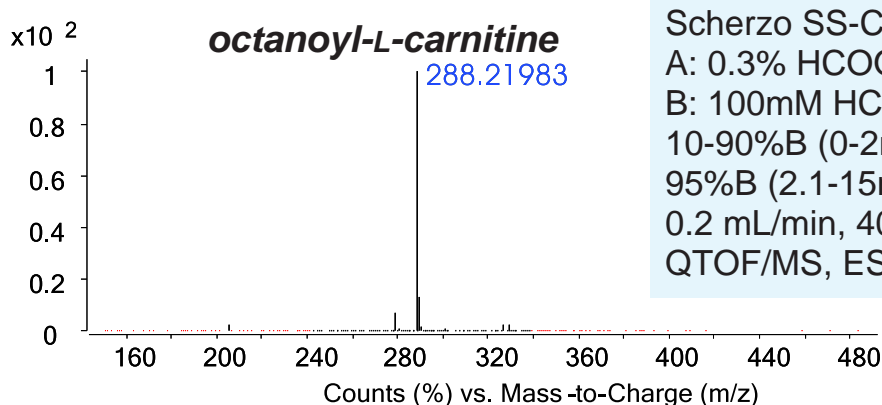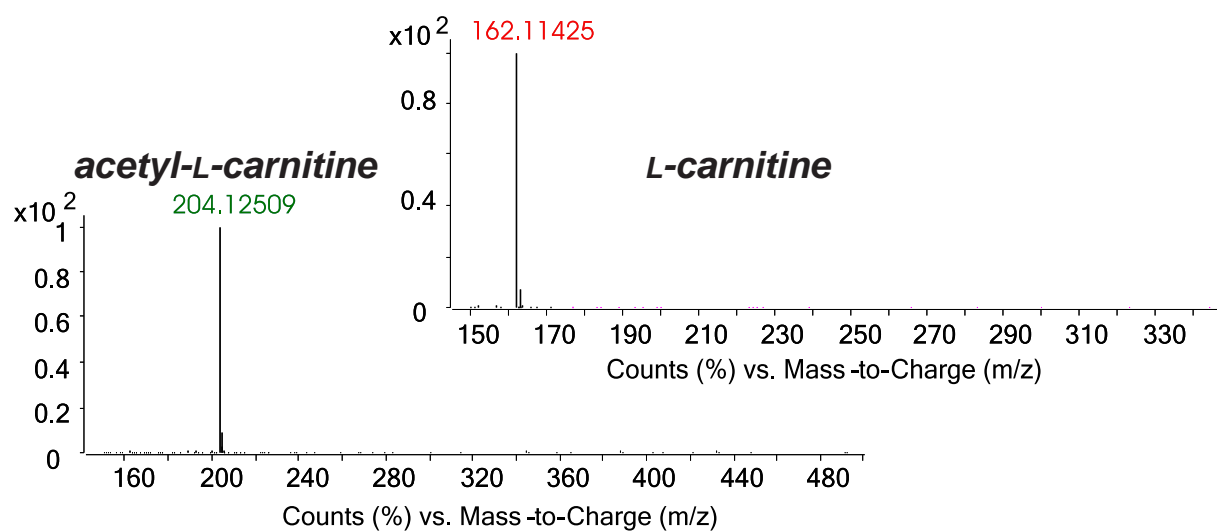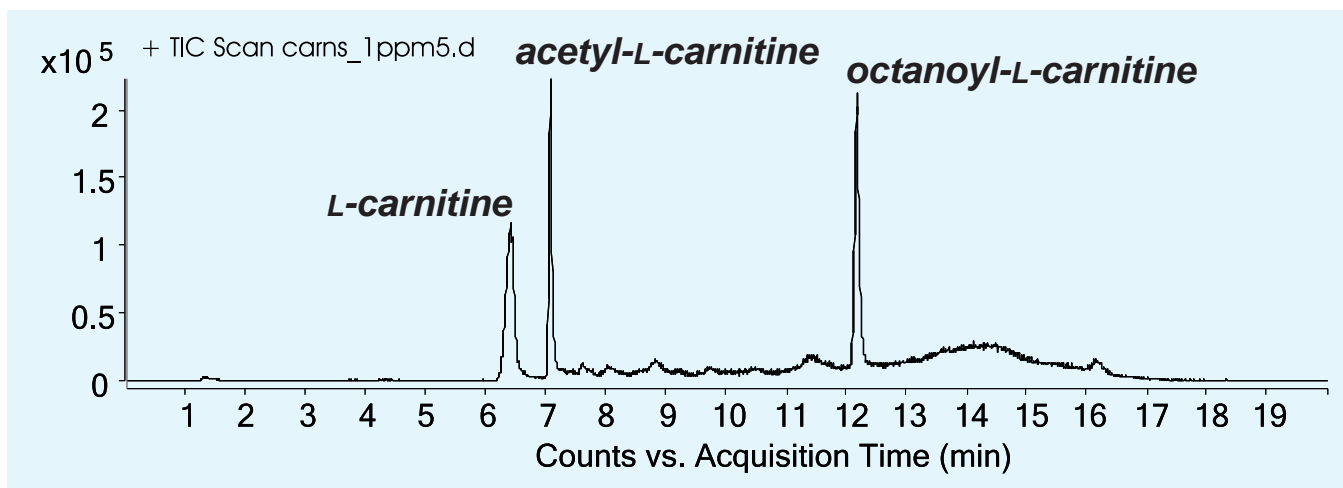

Scherzo SS-C18, 100 x 2 mm  
 A: 0.3% HCOOH  
 B: 100mM HCOONH<sub>4</sub> / ACN (50 / 50)  
 10-90%B (0-2min), 90-95%B (2-2.1min)  
 95%B (2.1-15min), 95-10%B (15-16min)  
 0.2 mL/min, 40°C, 10 μL injection  
 QTOF/MS, ESI, positive

Ref.) I.Nadaoka, E.Hatakeyama, C.Tanada, T.Sakamoto, S.Fukaya, T.Akiba, K.Inoue, Y.Yamano, T.Toyo'oka  
 Analytical Sciences, 30, 839 (2014)

Courtesy of Takashi AKIBA, Wakodo Co., Ltd., JAPAN