


Unison UK-Phenyl  
 Unison UK-C8  
 Unison UK-C18

50 x 3 mm

Application

## LC-MS: fungicides for citrus fruits

防かび剤のLC-MS分析

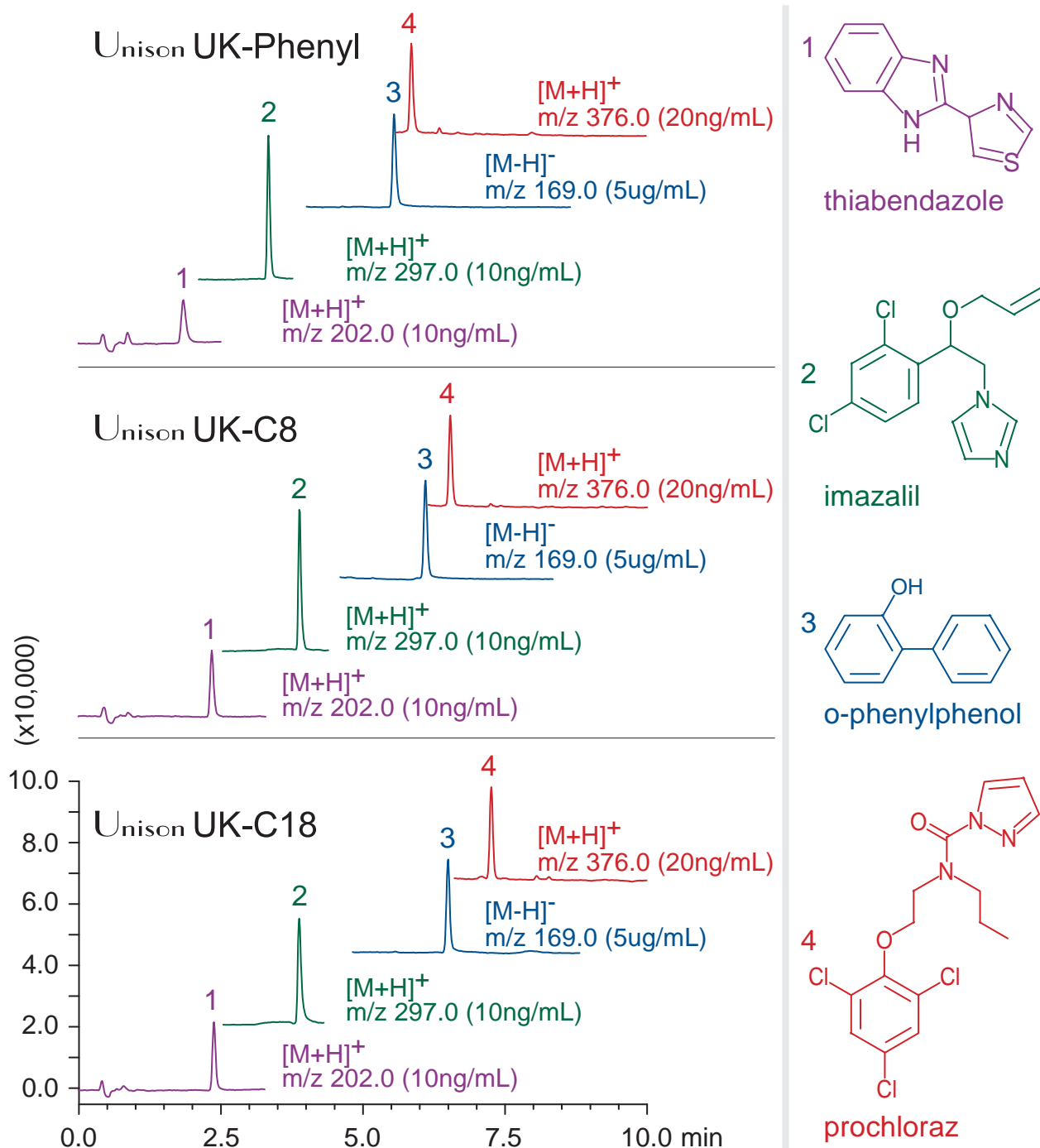

50 x 3 mm

A: water / acetic acid = 100 / 0.1

B: acetonitrile / acetic acid = 100 / 0.1

10-100 %B (0-10 min), 100 %B (10-15 min)

0.4 mL/min (4 MPa), 37°C, 20 uL (10ng - 5ug/mL, 0.1% AcOH+25% ACN)

Single Quad. MS (ESI, positive, negative)