


Unison US-C18

250 x 4.6 mm

Application

Fat-soluble vitamins

脂溶性ビタミン

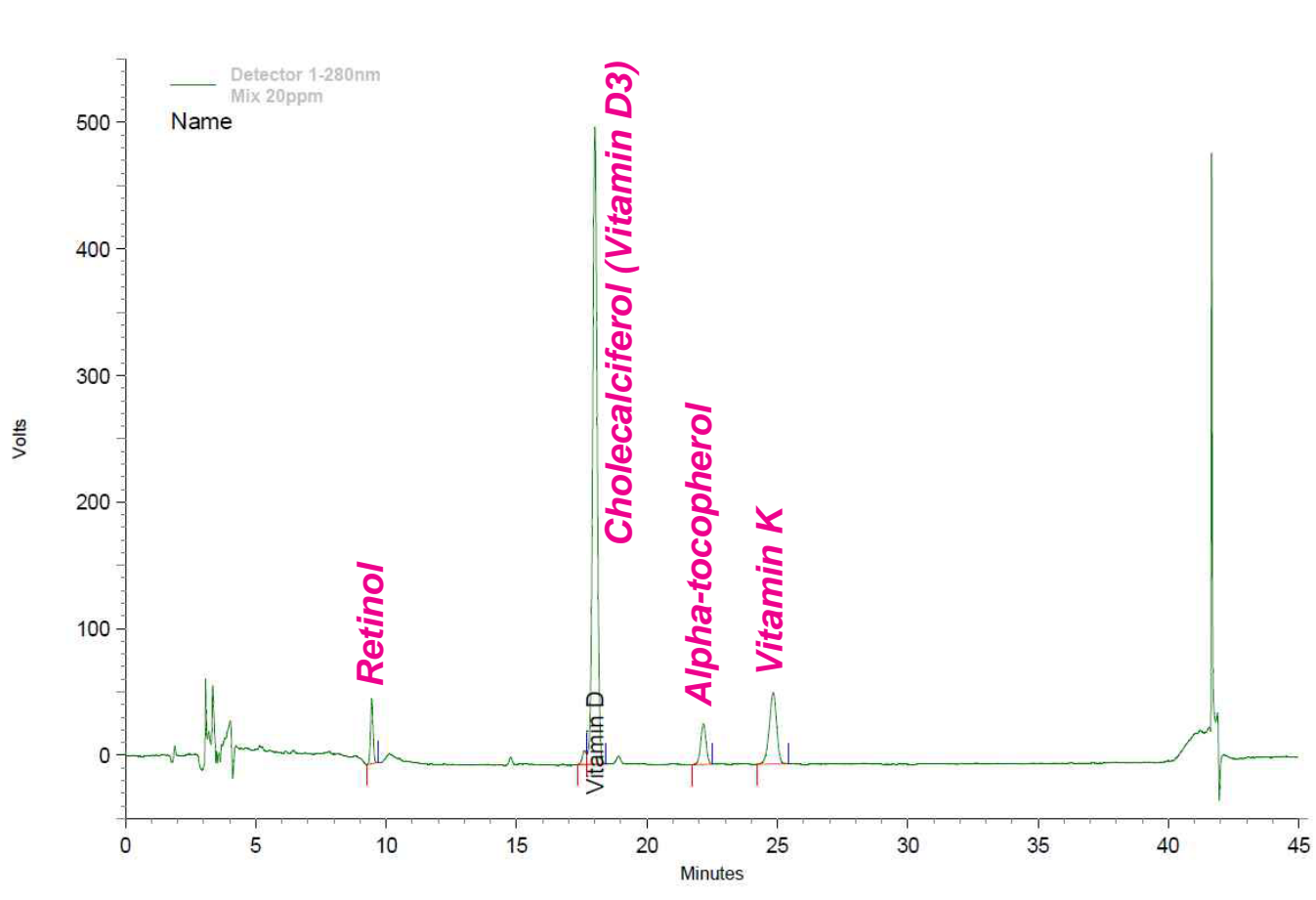

Unison US-C18, 250 x 4.6 mm

A: water

B: acetonitrile / methylene chloride = 90 / 10

85%B (0-5min), 85-100%B (5-6min), 100%B (6-37min)

100-85%B(37-37.1min), 85%B(37.1-45min)

0.8 mL/min, 37°C, 280nm

20 µL (STD 20ppm)

* Alpha-tocopherol, Retinol and Vitamin K dissolved with 2-propanol.
Cholecalciferol dissolved with methanol.

Courtesy of PEAKMAN SP, KOREA