
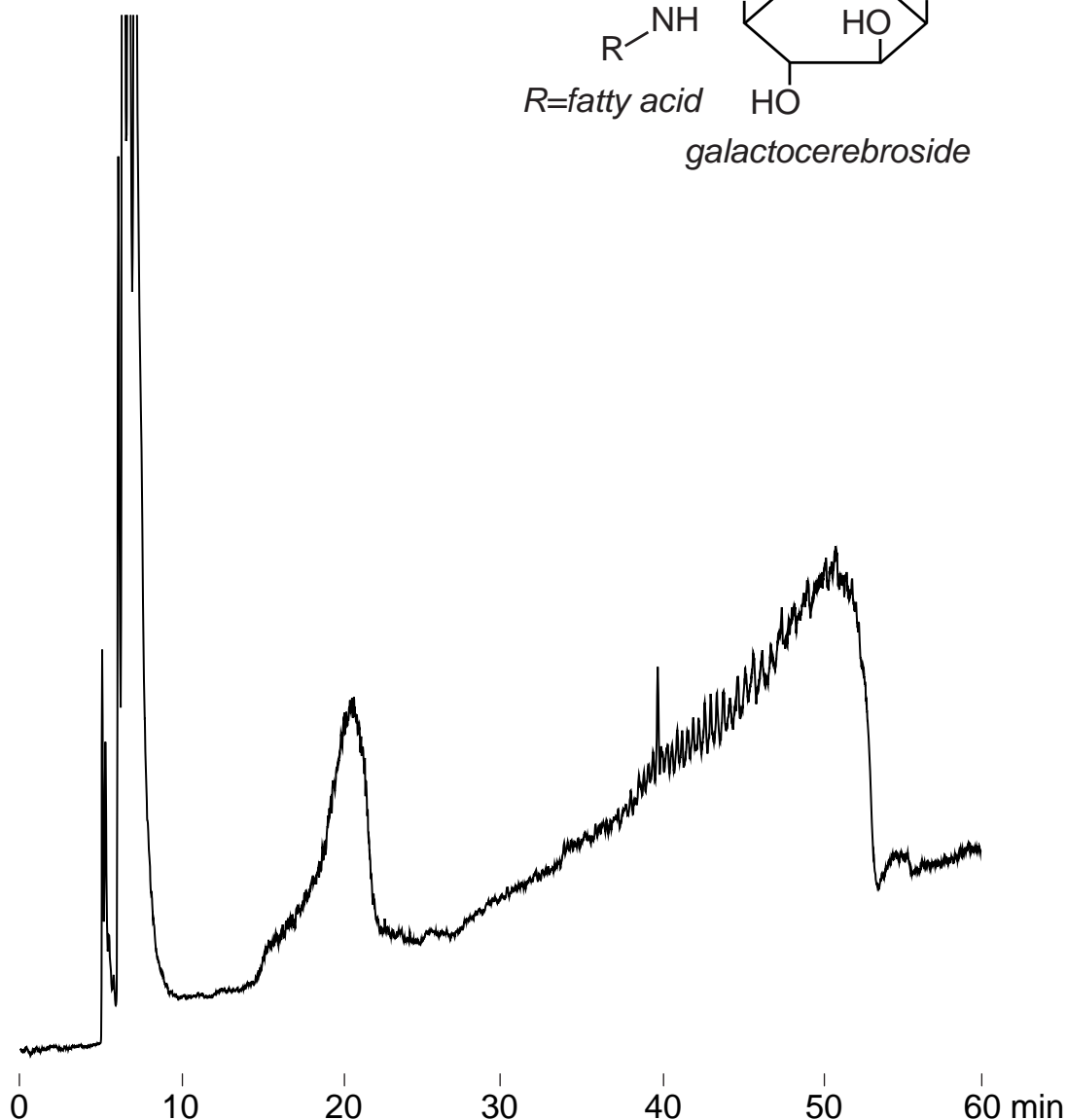


Unison UK-C8

250 x 4.6 mm

Application

Galactosylceramide from bovine brain
 スフィンゴ糖脂質 (ガラクトセレブロシド, 牛脳由来)

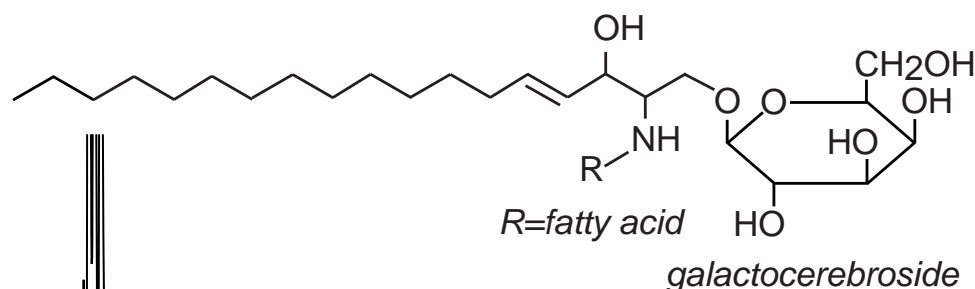

Unison UK-C8, 250 x 4.6 mm

A: acetonitrile / formic acid = 100 / 0.3

B: tetrahydrofuran

0-10%B (0-10min), 10-100%B (10-50min), 100%B (50-60min)

1mL/min (6MPa), 37°C, ELSD (spray chamber 50°C, drift chamber 100°C)

5µL (4mg/mL, chloroform / methanol = 2 / 1)