
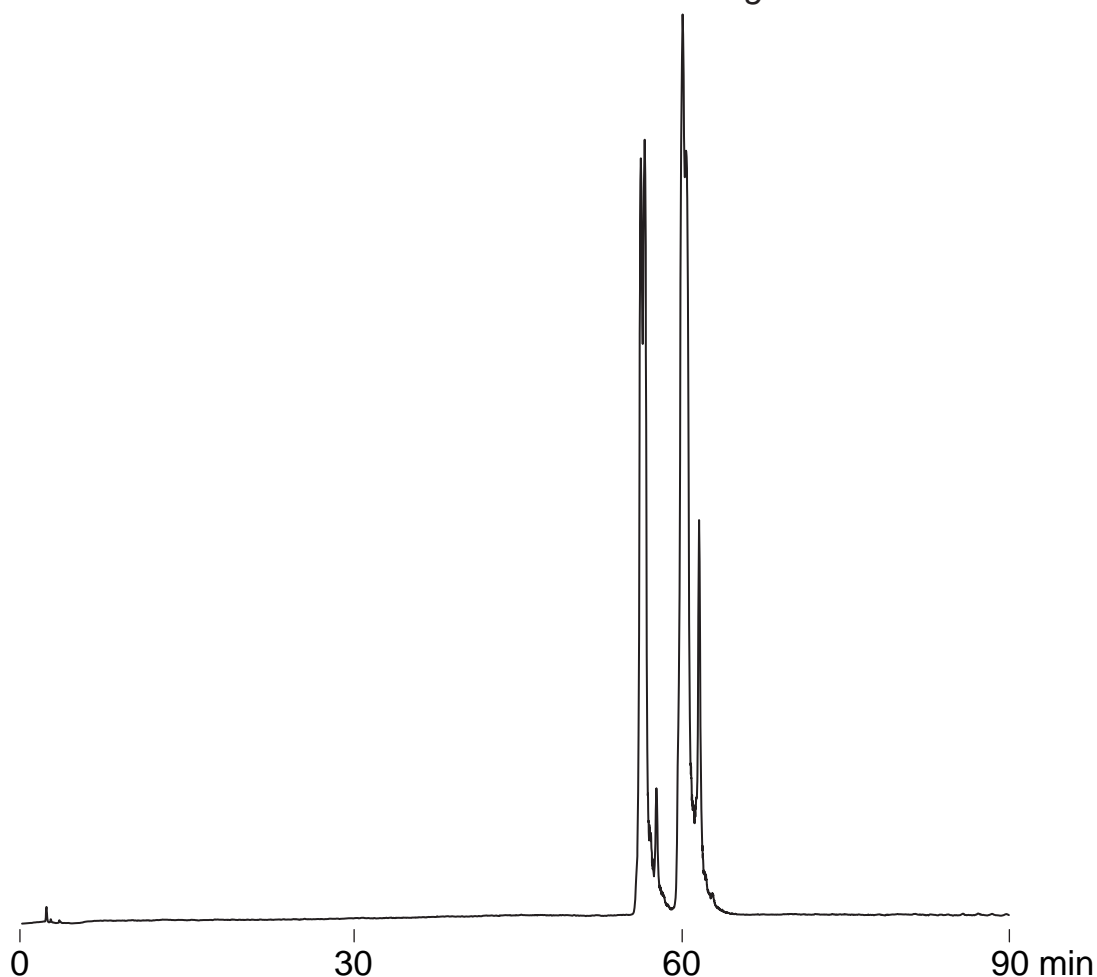


Unison UK-Silica

250 x 4.6 mm

Application

Galactosylceramide from bovine brain
 スフィンゴ糖脂質 (ガラクトセレブロシド, 牛脳由来)

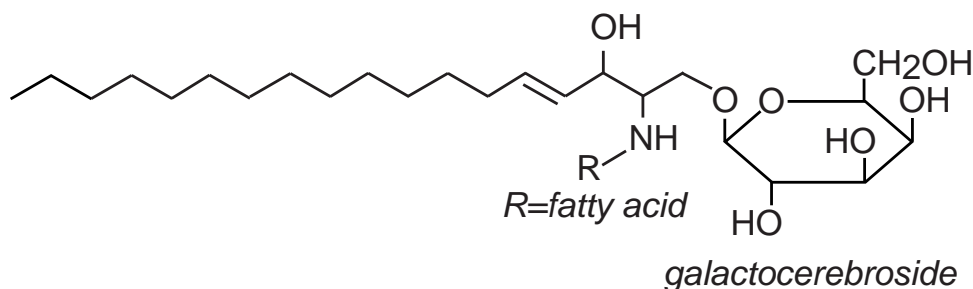

Unison UK-Silica, 250 x 4.6 mm

A: hexane / water / formic acid = 100 / 0.1 / 0.1

B: tetrahydrofuran / methanol / formic acid = 80 / 20 / 0.1

0-60% (0-90min)

1mL/min (4MPa), 37°C, ELSD (spray chamber 50°C, drift chamber 100°C)

5µL (4mg/mL, chloroform / methanol = 2 / 1)