



Unison UK-Phenyl  
Unison UK-C8  
Unison UK-C18

Cadenza CW-C18  
Cadenza CX-C18  
Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

## Acid Green 3 ( Artificial coloring )

ギニアグリーン B (人工着色料)

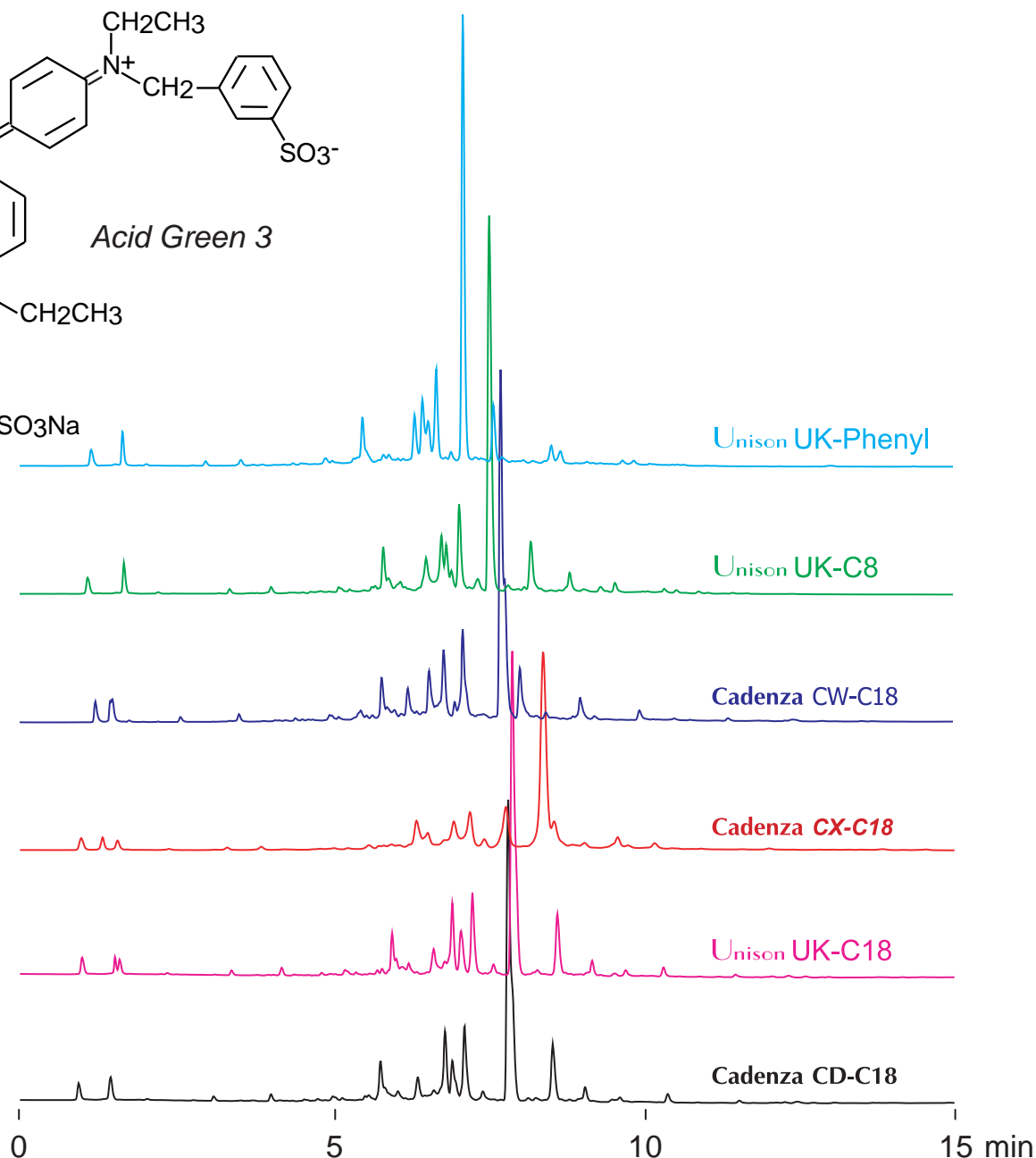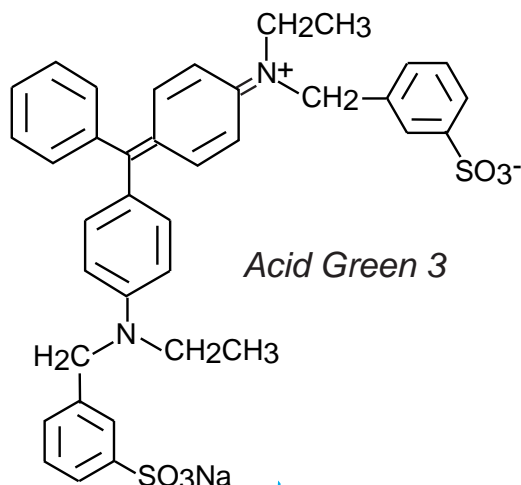

100 x 4.6 mm

A: 100mM ammonium acetate

B: acetonitrile / 100mM ammonium acetate = 90 / 10

20-100%B (0-15min)

1mL/min (9MPa), 37°C, 260nm

10µL (1mg/mL, 100mM ammonium formate)