




Cadenza CW-C18
Cadenza CD-C18

250 x 4.6 mm

HPLC COLUMN APPLICATION

Polysorbate 20 (Nonionic surfactant)

ポリソルベート20 (非イオン性界面活性剤)

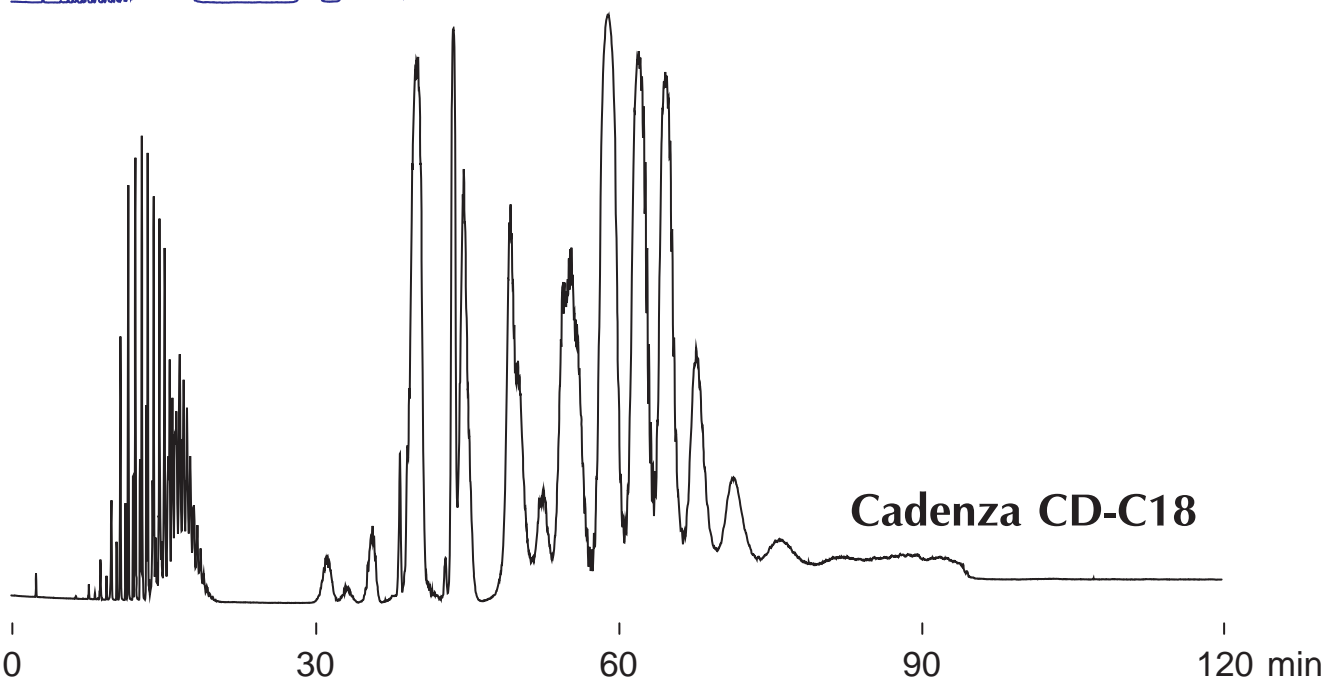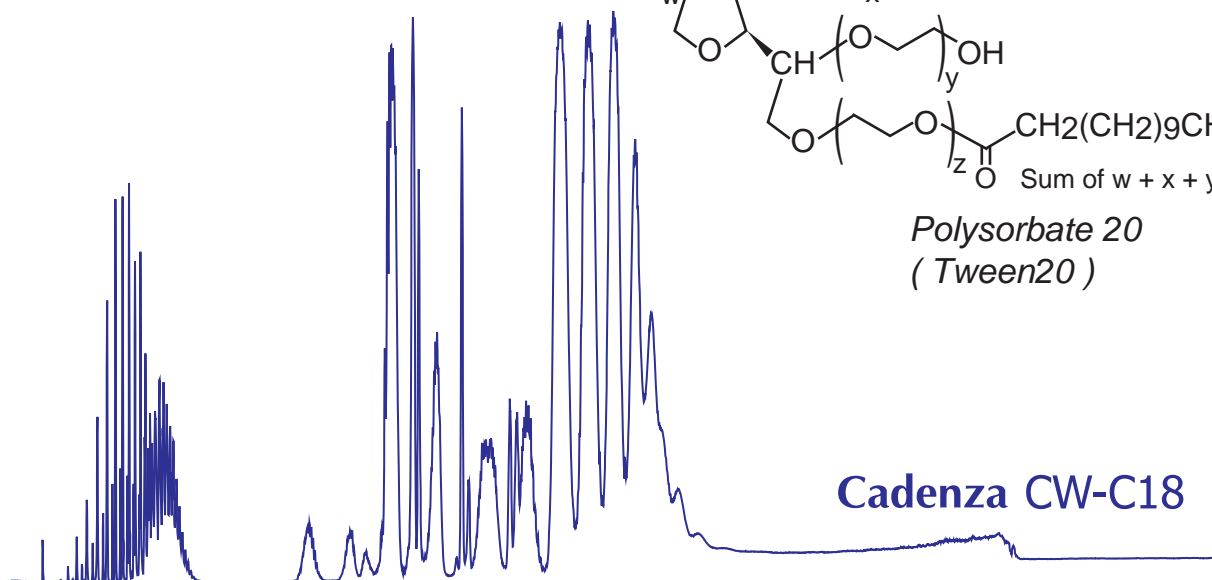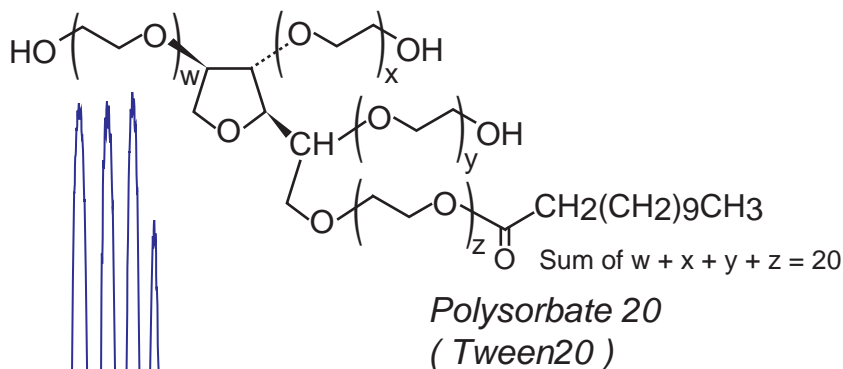

250 x 4.6 mm
A: water / acetic acid = 100 / 0.1
B: acetonitrile / acetic acid = 100 / 0.1
13-100%B (0-60min), 100%B (60-120min)
1.0mL/min (14MPa), 60°C
ELSD (spray chamber 70°C, drift chamber 120°C)
20µL (50mg/mL, 100mM ammonium formate)