

Presto FF-C18


Cadenza CD-C18

250 x 4.6 mm


HPLC COLUMN APPLICATION

Polysorbate 20 (Nonionic surfactant)

ポリソルベート20 (非イオン性界面活性剤)

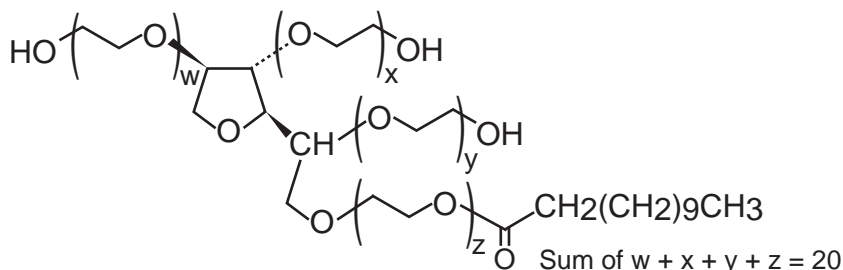


Polysorbate 20
(Tween20)

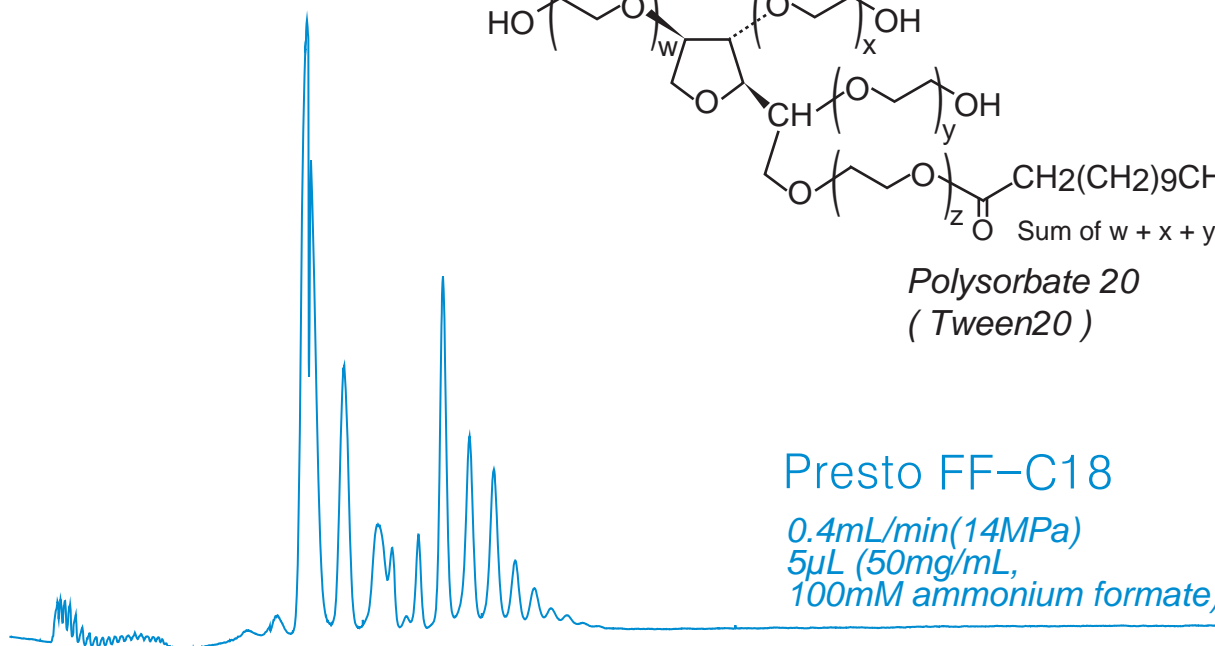

Presto FF-C18

0.4mL/min(14MPa)
5μL (50mg/mL,
100mM ammonium formate)

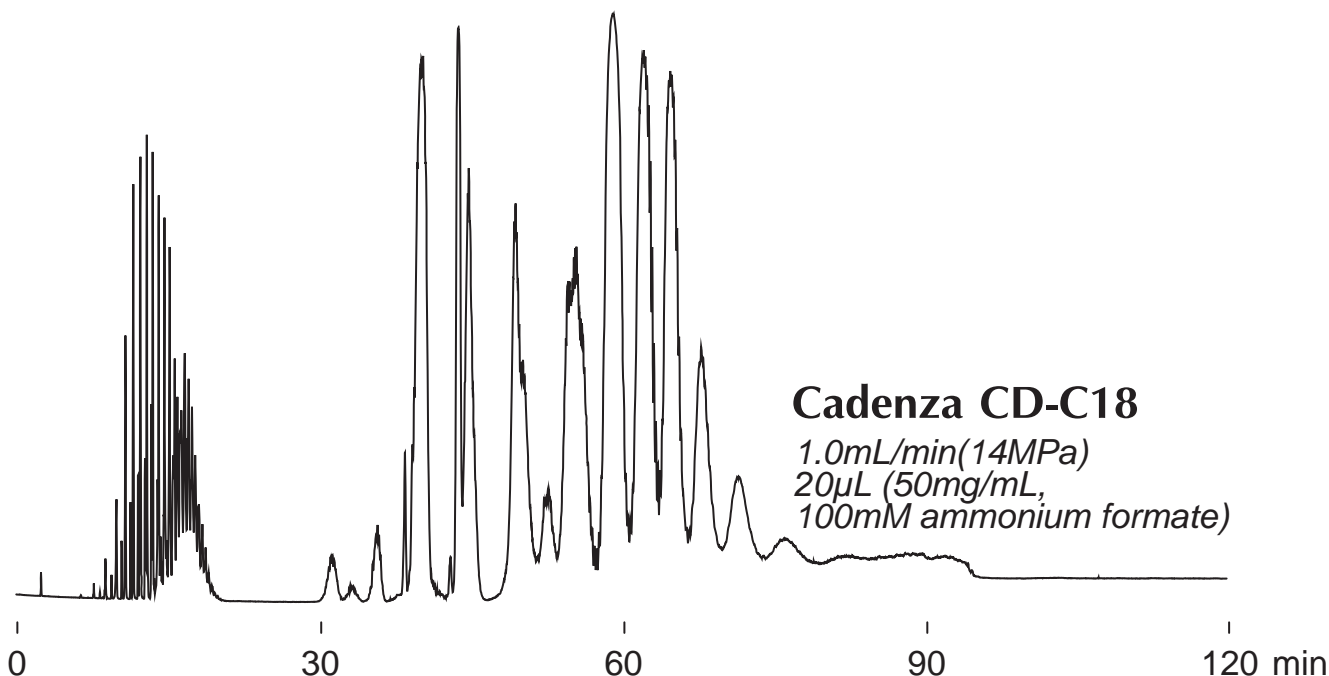

Cadenza CD-C18

1.0mL/min(14MPa)
20μL (50mg/mL,
100mM ammonium formate)

250 x 4.6 mm

A: water / acetic acid = 100 / 0.1

B: acetonitrile / acetic acid = 100 / 0.1

13-100%B (0-60min), 100%B (60-120min)

60°C, ELSD (spray chamber 70°C, drift chamber 120°C)