


Unison UK-C18
Cadenza CD-C18

150 x 4.6 mm


HPLC COLUMN APPLICATION

Chloromandelic acid isomers


クロロマンデル酸異性体

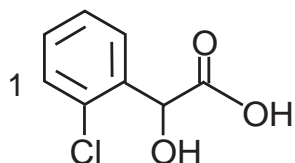


2-chloromandelic acid

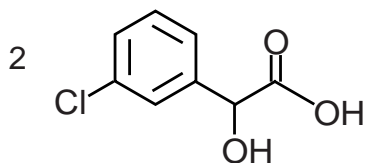

3-chloromandelic acid

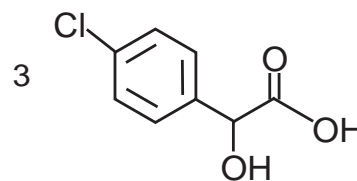

4-chloromandelic acid

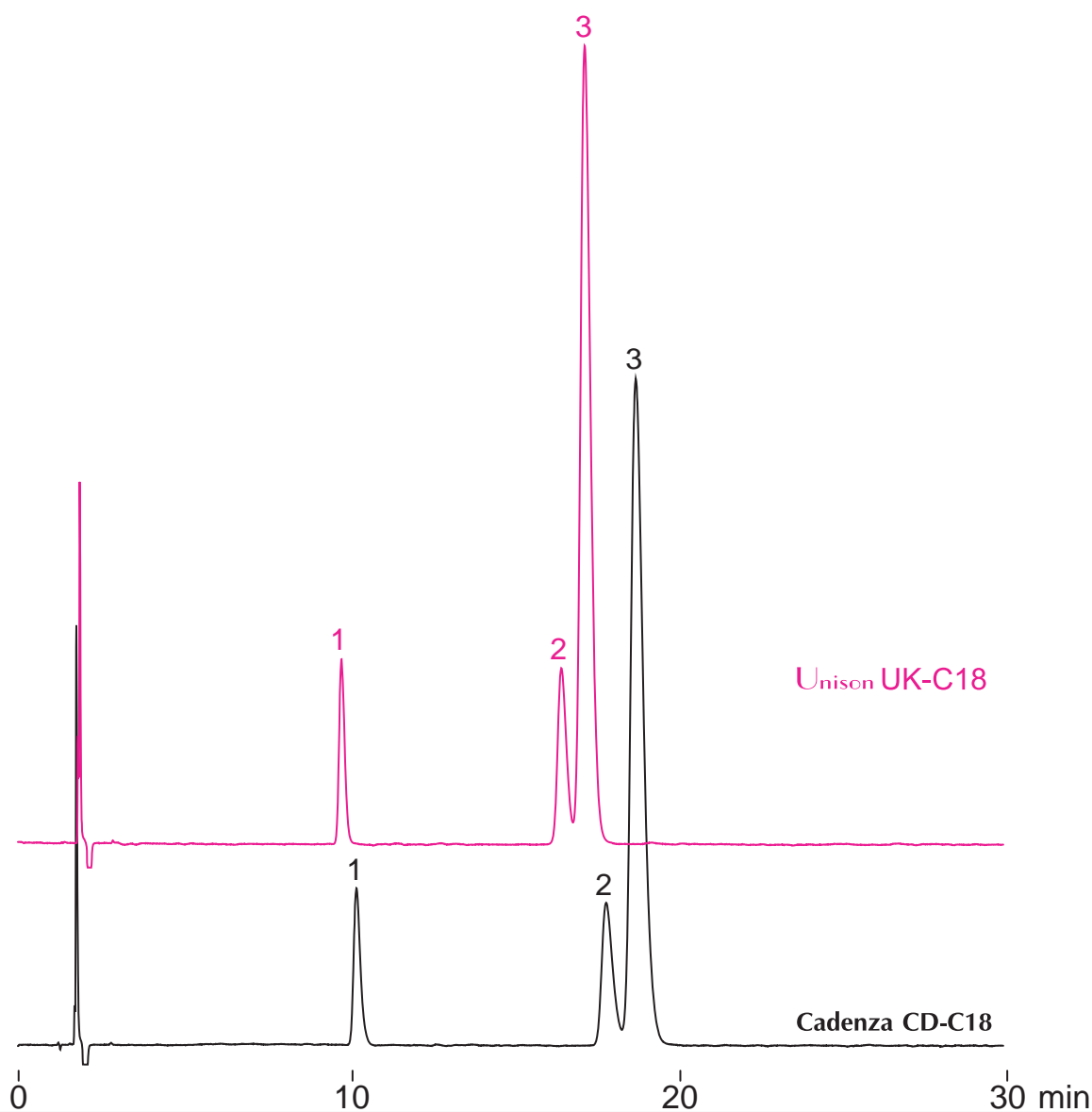

150 x 4.6 mm

acetonitrile / water / formic acid = 15 / 85 / 0.5

1.0mL/min (13MPa), 37°C, 235nm

10µL (1mg/mL, acetonitrile / water / formic acid = 50 / 50 / 1)