



Unison UK-Phenyl

150 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Organophosphorus pesticides (3)

有機リン系殺虫剤 (3) の LC-MS 分析

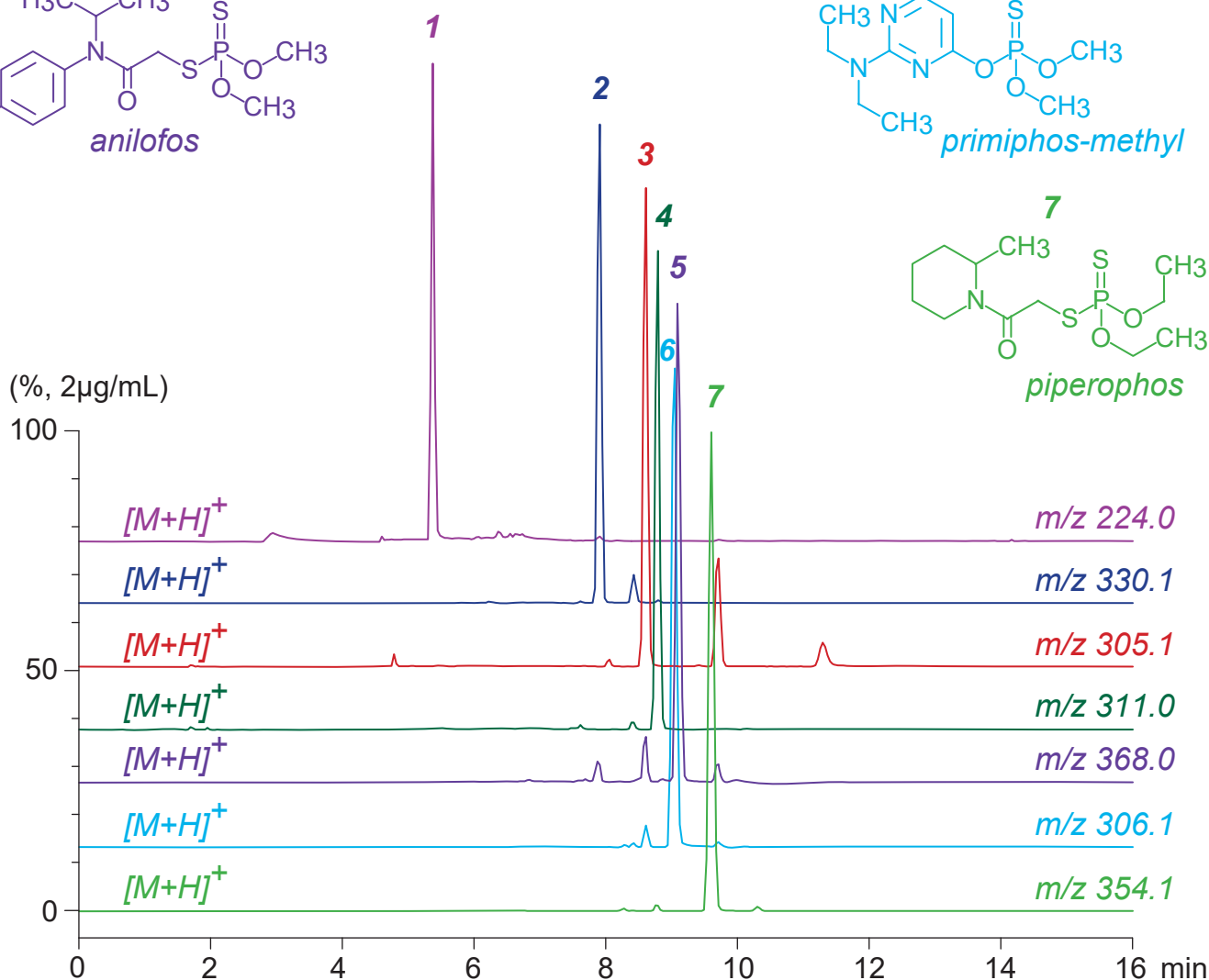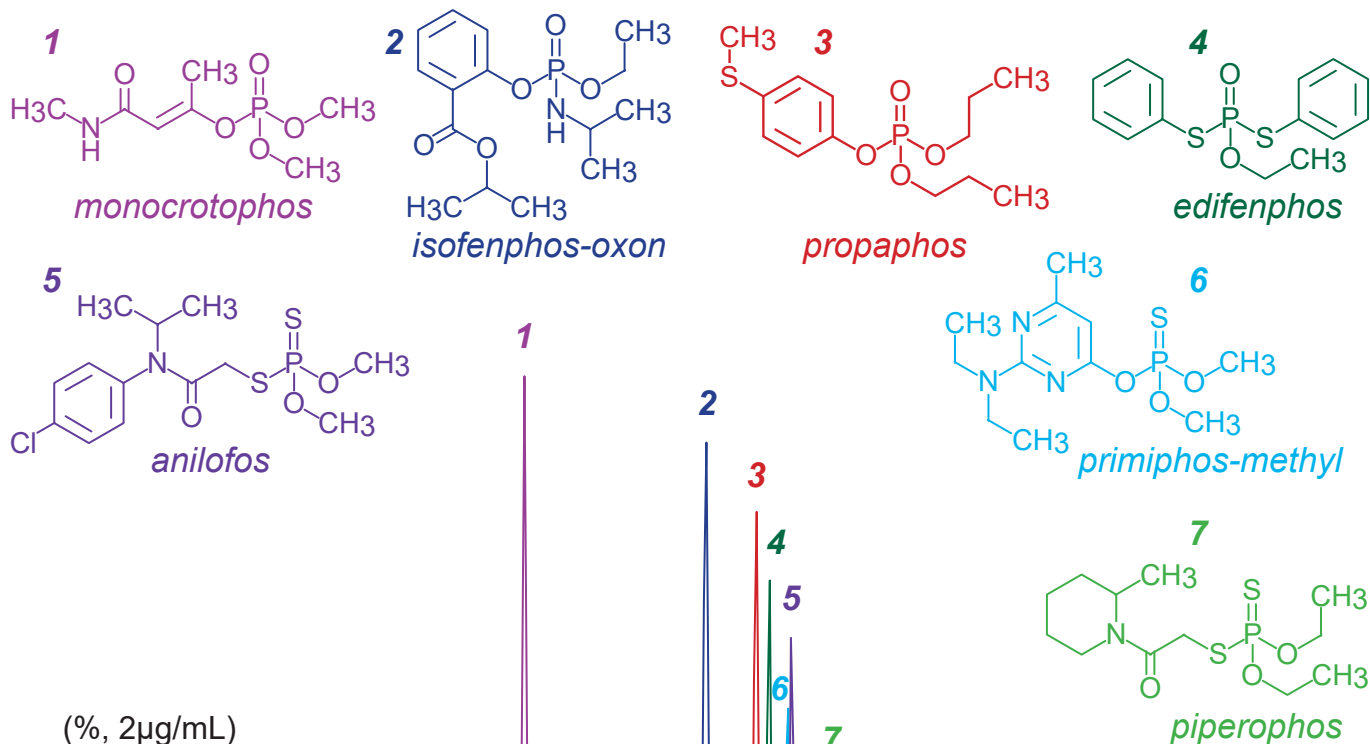

Unison UK-Phenyl, 150 x 3 mm

A: water / formic acid = 100 / 0.1

B: acetonitrile / formic acid = 100 / 0.1

0-65%B (0-5min), 65%B (5-16min)

0.4mL/min (8MPa), 37°C, Single Quad. MS (ESI)

5 μ L (2 μ g/mL, ACN / water / HCOOH = 50 / 50 / 0.1)