

Unison UK-Phenyl Cadenza CW-C18


Unison UK-C8 Cadenza CX-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm


HPLC COLUMN APPLICATION

Coumatetralyl (Coumarin anticoagulant)

クマテトラリル (クマリン系抗血液凝固薬)

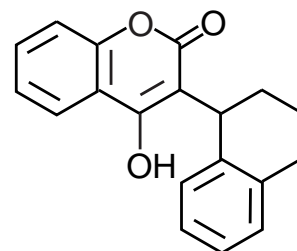

coumatetralyl

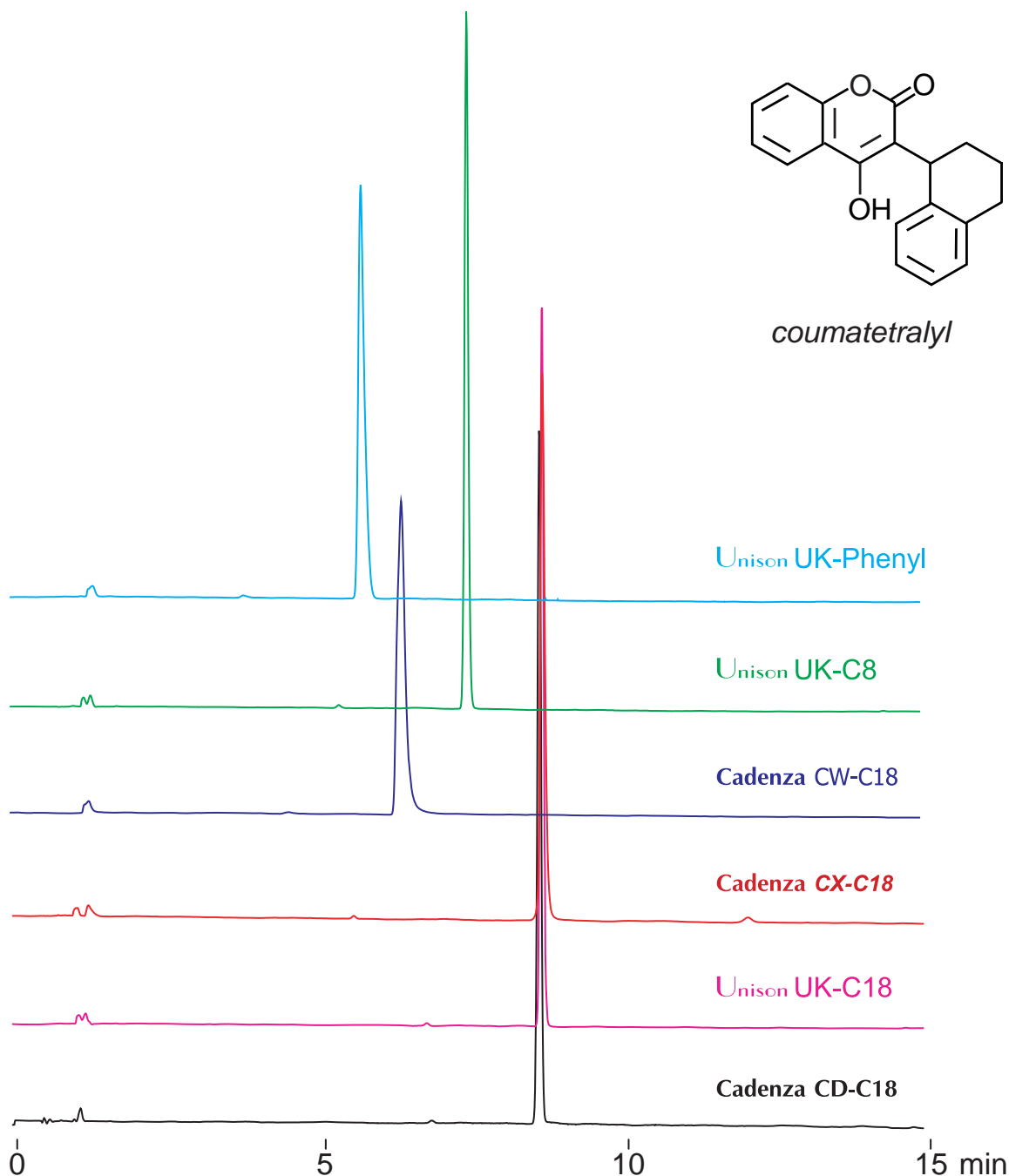

100 x 4.6 mm

A: water / formic acid = 100 / 0.1

B: acetonitrile / formic acid = 100 / 0.1

35-100%B (0-15min)

1.0mL/min (9MPa), 37°C, 280nm

10μL (0.2mg/mL, acetonitrile / water / formic acid = 50 / 50 / 1)