

Unison UK-Phenyl Cadenza CW-C18


Unison UK-C8 Cadenza CX-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

Pitavastatin (Hypolipidemic agent)

ピタバスタチン (脂質異常症治療薬)

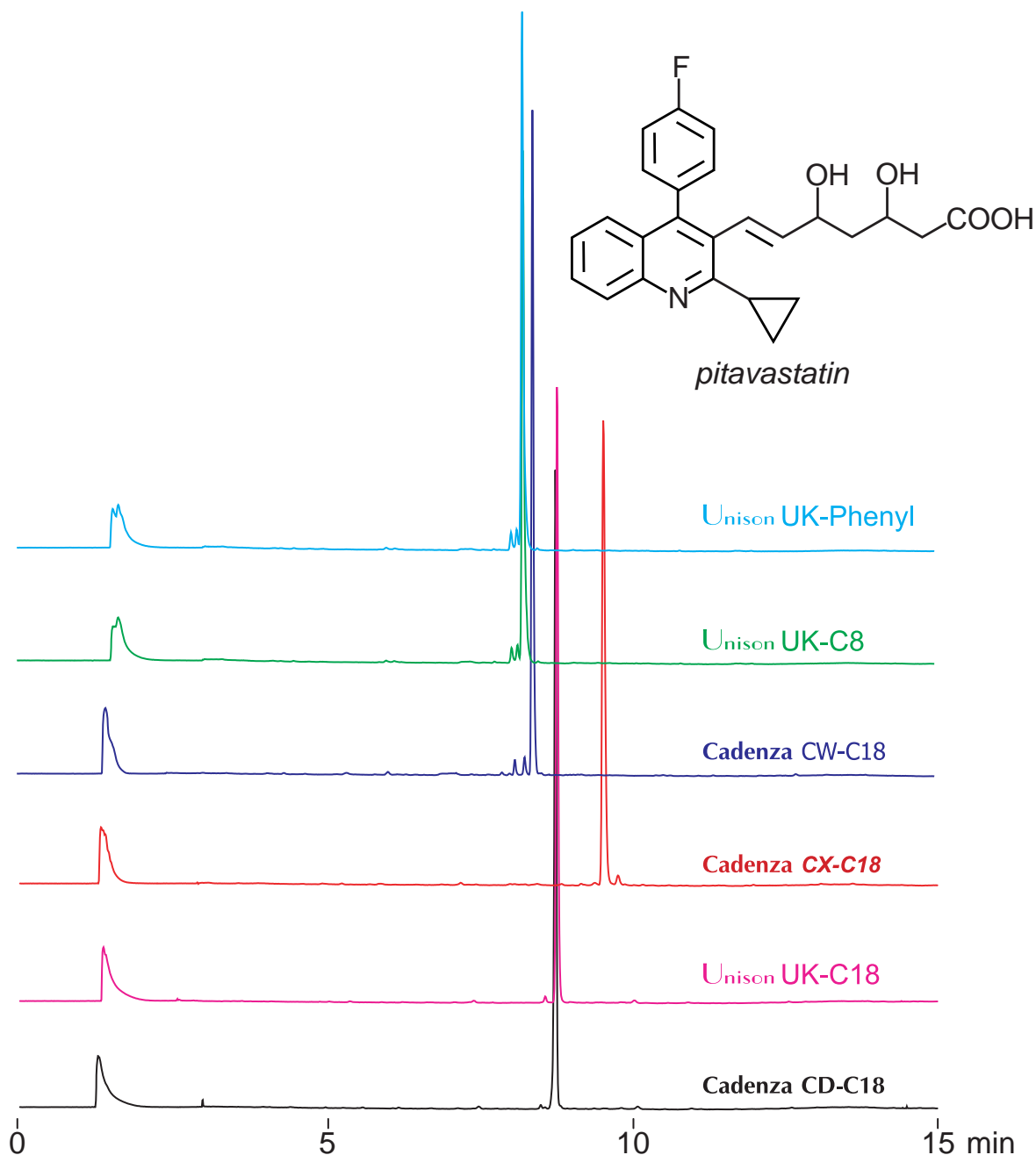

100 x 4.6 mm

A: 100mM ammonium formate

B: 100mM ammonium formate / acetonitrile = 10 / 90

0-100%B (0-15min)

1.0mL/min (10MPa), 37°C, 250nm

10µL (0.05mg/mL, DMSO / 100mM ammonium formate = 75 / 25)