

Unison UK-Phenyl Cadenza CX-C18



Unison UK-C8 Cadenza CW-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

Rhododendrol (Tyrosinase inhibitor)

ロドデンドロール (チロシナーゼ阻害剤)

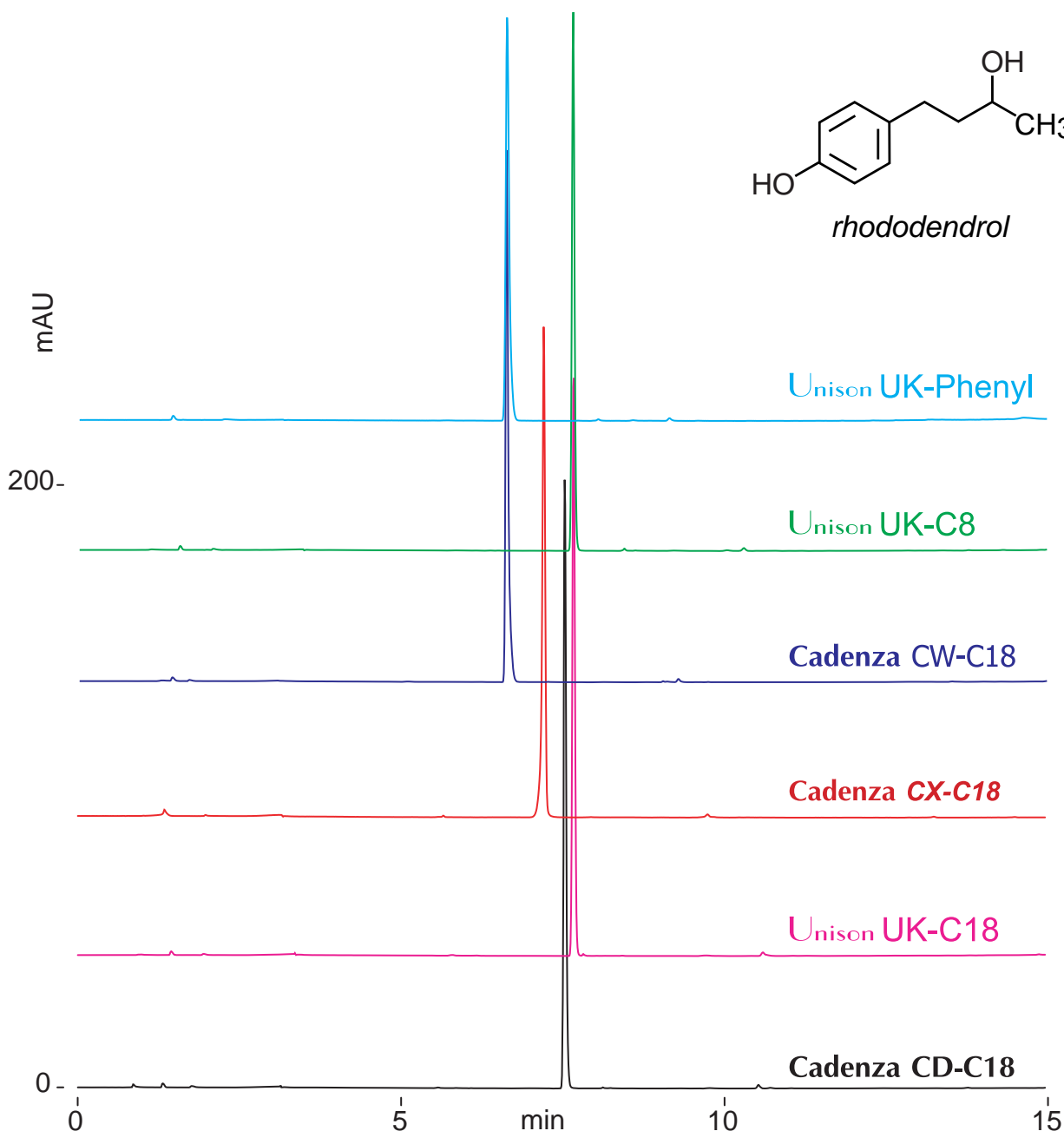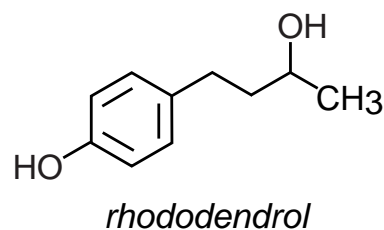

100 x 4.6 mm

A: water / formic acid = 100 / 0.1

B: water / acetonitrile / formic acid = 10 / 90 / 0.1

0-90%B (0-15min)

1.0mL/min (9MPa), 37°C, 280nm

10µL (0.3mg/mL, water / formic acid = 100 / 1)