


Cadenza CD-C18 HT

150 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Antibiotics (5)

抗生物質のLC-MS/MS分析 (5)

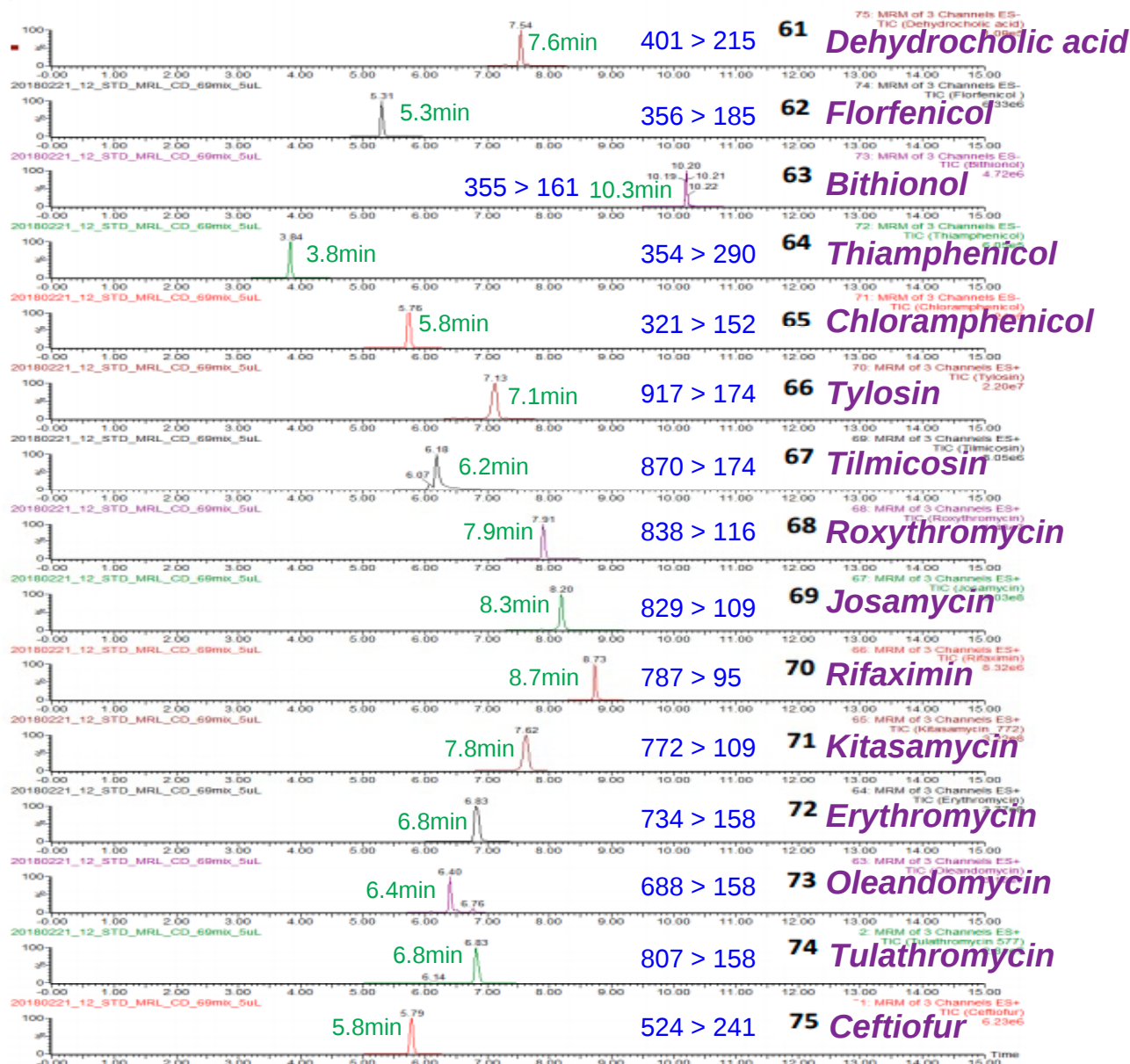

Cadenza CD-C18 HT, 150 x 2 mm

A: 0.1% formic acid in water

B: 0.1% formic acid + 2mM ammonium formate in acetonitrile

10%B (0-1min), 10-40%B (1-6min), 40-95%B (6-10min),

95%B (10-15min), 95-10%B (15-15.1min), 10%B (15.1-20min)

0.4mL/min, 40°C, Triple Quad. MS (ESI, Positive/Negative)

5μL

Courtesy of PEAKMAN SP, KOREA