


Unison UK-C18
 Unison UK-C8
 Unison UK-C1
 Unison UK-Phenyl

150 x 4.6 mm


HPLC COLUMN APPLICATION

Sclareol, (+)-sclareolide (Sesquiterpenes)


スクラレオール, (+)スクラレオリド (テルペン類)

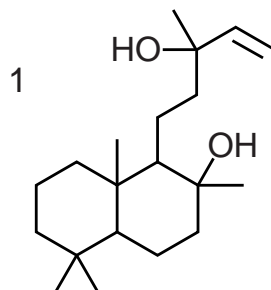

sclareol
(3mg/mL)

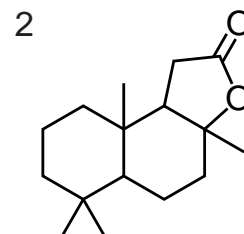

(+)-sclareolide
(17mg/mL)

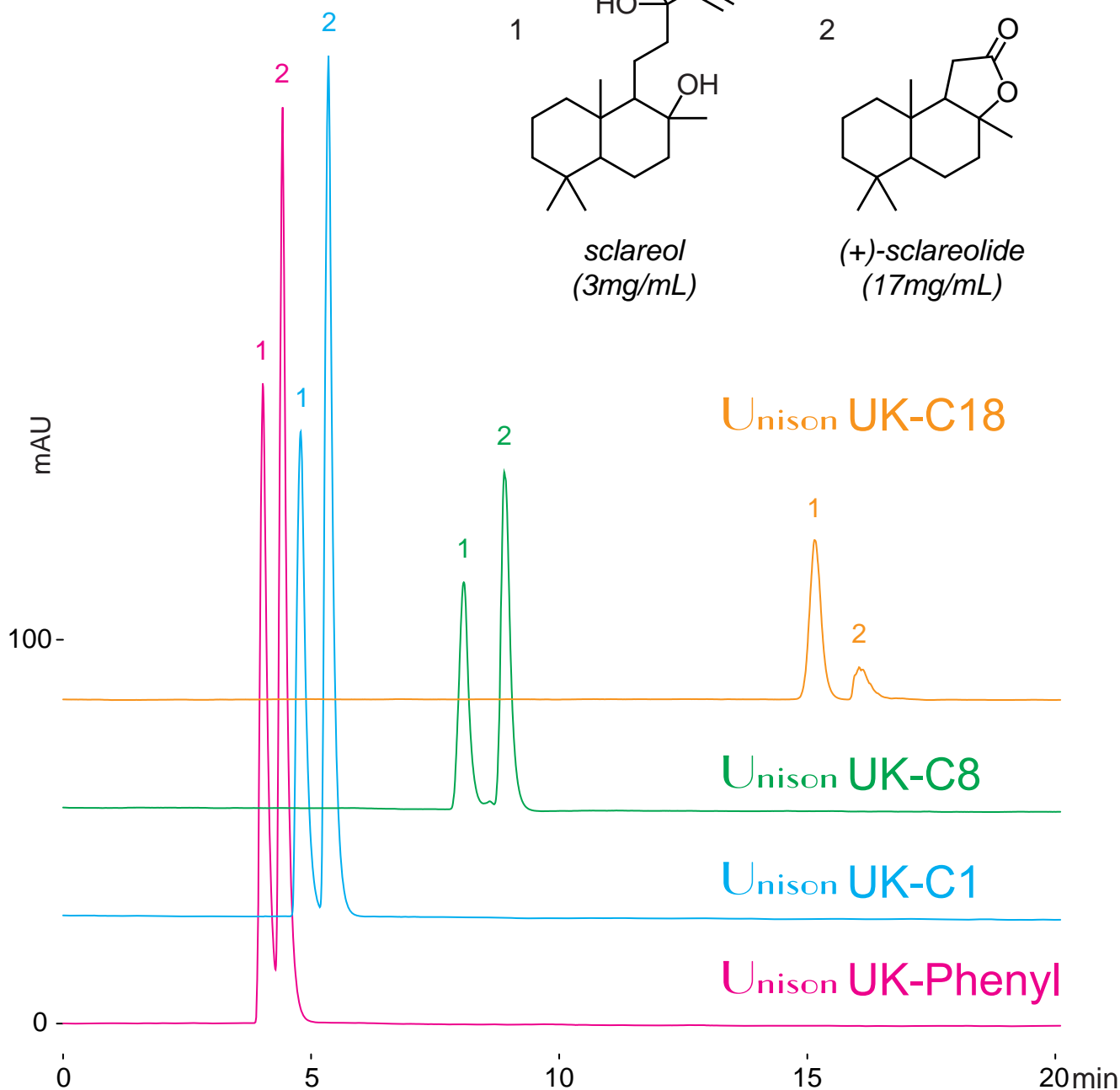

150 x 4.6 mm
 acetonitrile / water / acetic acid = 55 / 45 / 0.1
 1.0mL/min (10-12MPa), 37°C
 ELSD (spray chamber 70°C, drift chamber 70°C)
 10µL (3-17mg/mL, tetrahydrofuran)