

Unison UK-C1

Unison UK-Phenyl

Unison UK-C8

Unison UK-C18

Cadenza CX-C18

Cadenza CW-C18


Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

5,6,11,12-Tetraphenylnaphthacene (Organic light-emitting diode material)

5,6,11,12-テトラフェニルナフタセン
(有機エレクトロルミネッセンス材料)

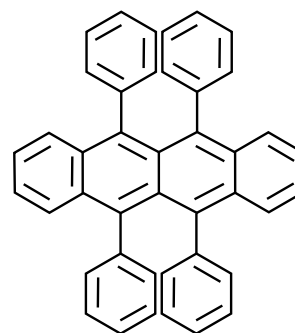

5,6,11,12-tetraphenylnaphthacene

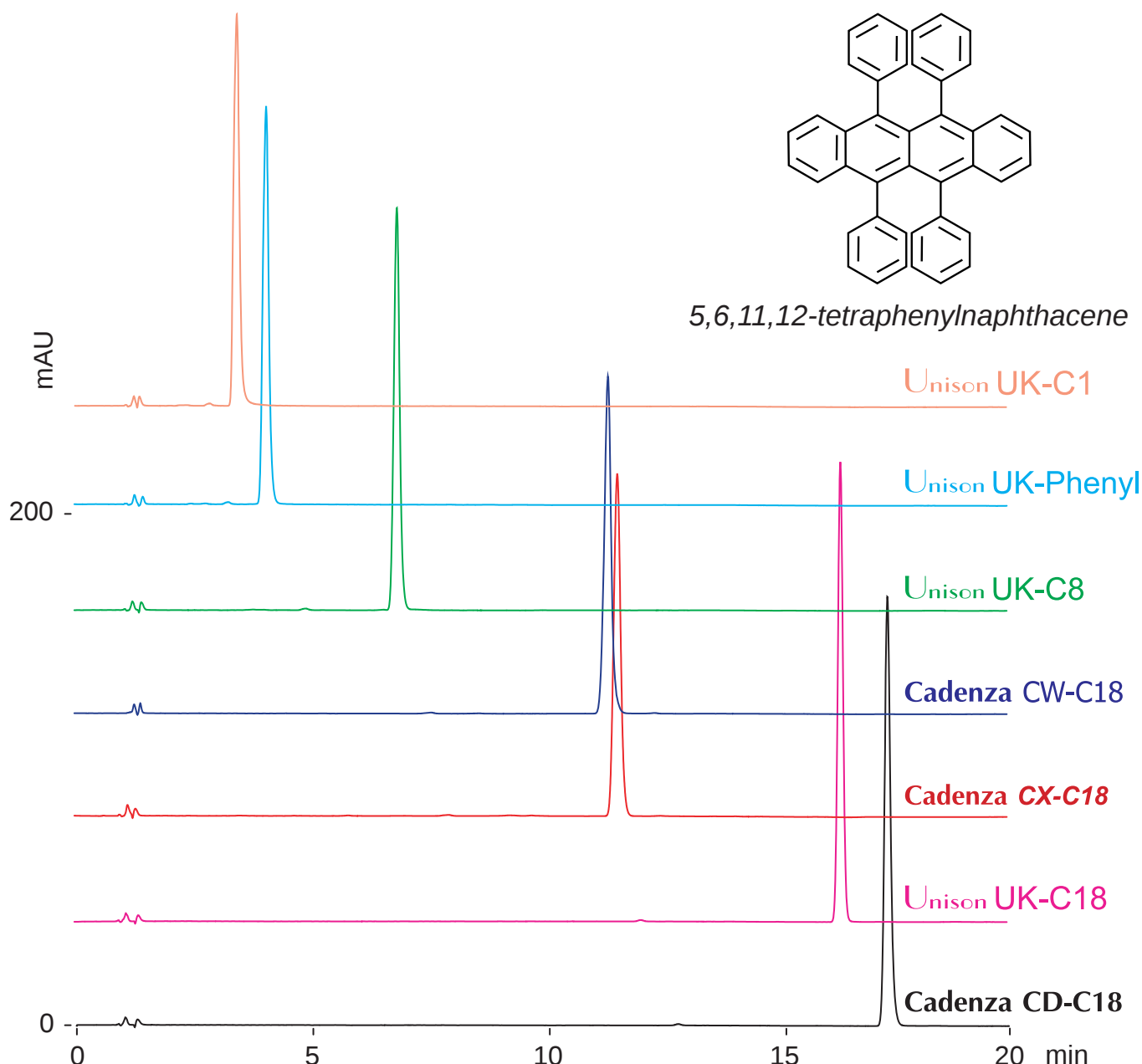

100 x 4.6 mm

A: water / acetonitrile / acetic acid = 50 / 50 / 0.1

B: acetonitrile / acetic acid = 100 / 0.1

60-100%B (0-15min), 100% (15-20min)

1.0mL/min (5MPa), 300nm, 37°C

10μL (0.025mg/mL, acetonitrile / THF /acetic acid= 50 / 50 / 1)