


Unison UK-Phenyl Unison UK-C1
Unison UK-C8 Cadenza CX-C18
Unison UK-C18 Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

N,N'-Bis [4- (diphenylamino) phenyl] -N,N'-diphenylbenzidine (Organic light-emitting diode material)

N,N'-ビス [4- (ジフェニルアミノ) フェニル] -N,N'-ジフェニルベンジジン
(有機EL材料)

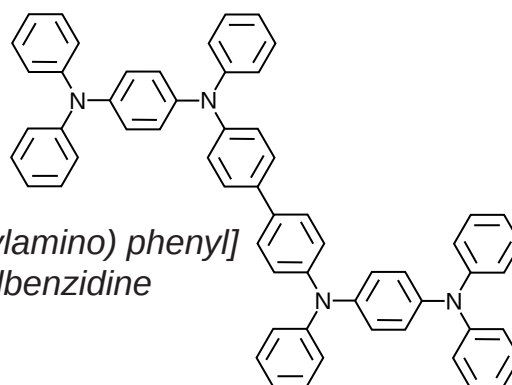

N,N'-bis [4- (diphenylamino) phenyl]
-N,N'-diphenylbenzidine

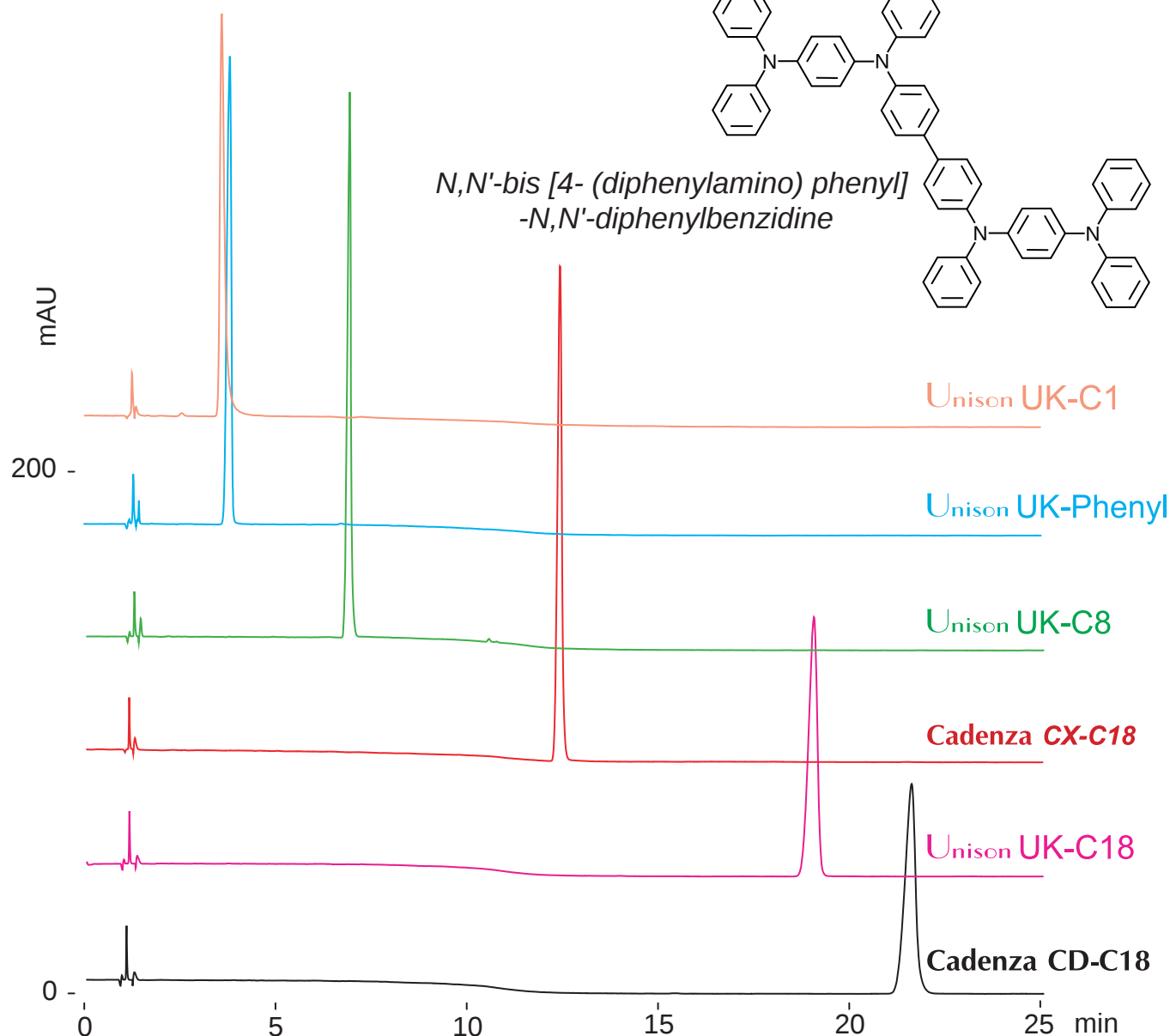

100 x 4.6 mm

A: 100mM ammonium formate / acetonitrile = 50 / 50

B: acetonitrile

70-100%B (0-10min), 100%B (10-25min)

1.0mL/min (6MPa), 37°C, 260nm

10μL (0.2mg/mL, THF / ACN = 50 / 50)