

Unison UK-Phenyl

Nardis ND-RX


Unison UK-C18

100 x 3 mm


HPLC COLUMN APPLICATION

LC-MS: Tributyltin (IV) chloride (Antifouling biocides)


塩化トリブチルスズ (M)(防汚殺生物剤) の LC-MS 分析

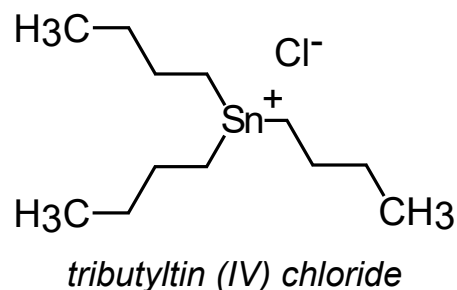


Unison UK-Phenyl

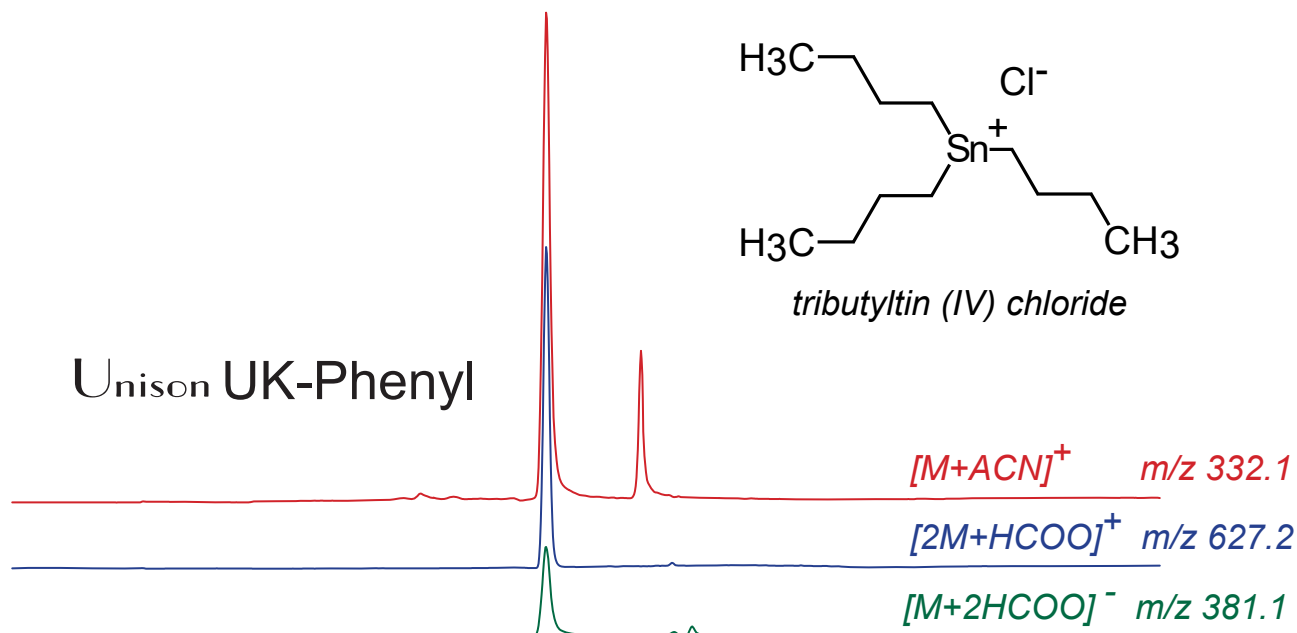

Nardis ND-RX

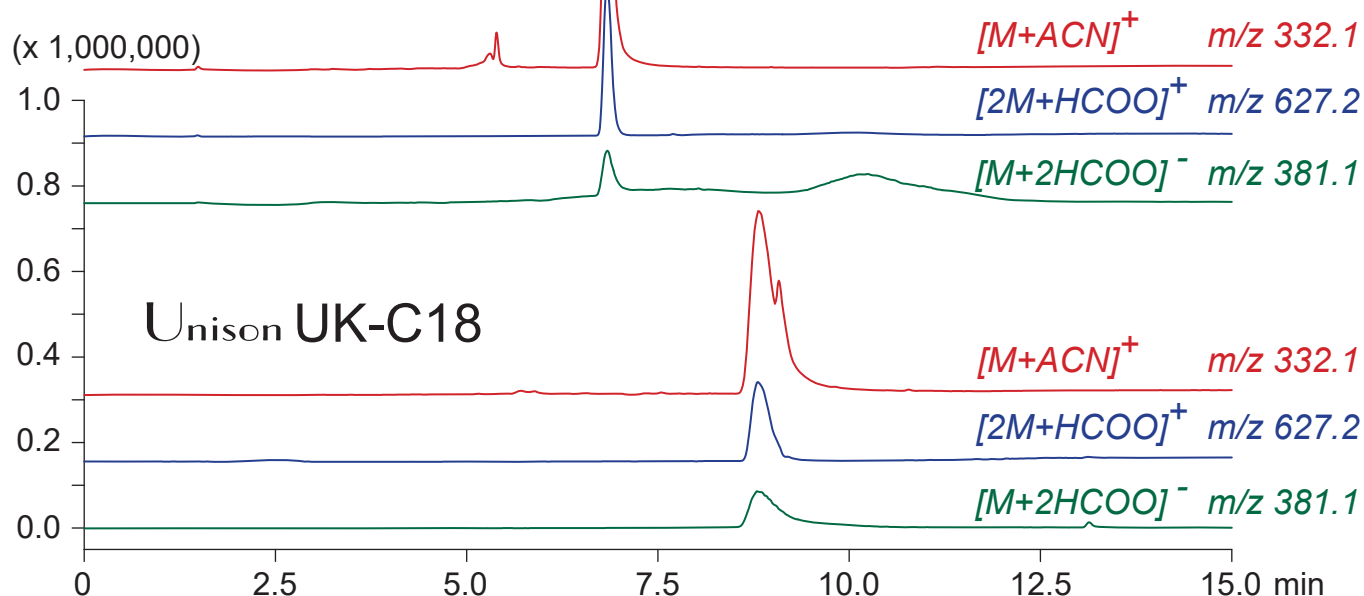

Unison UK-C18

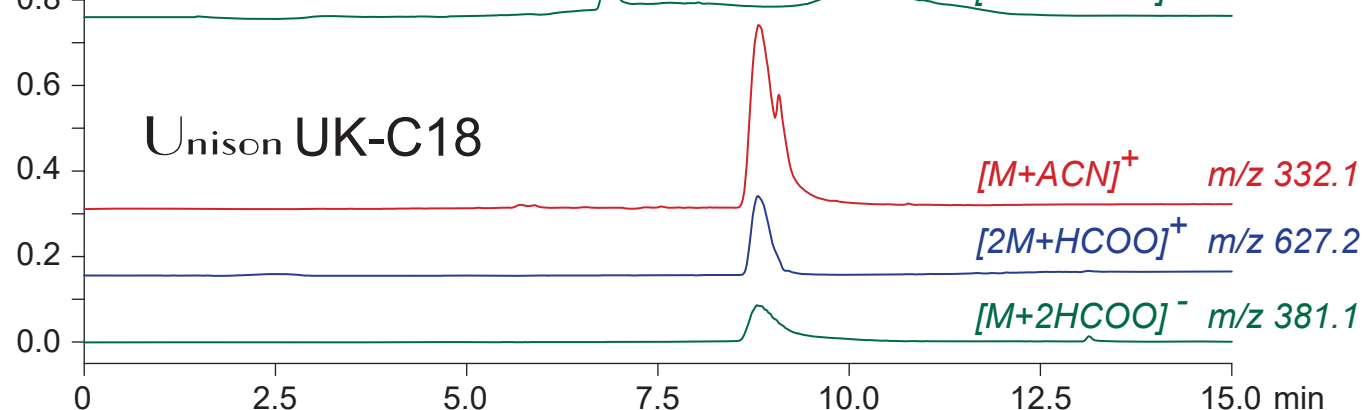

100 x 3 mm

A: water / formic acid = 100 / 0.5

B: acetonitrile / formic acid = 100 / 0.5

0-100%B (0-10min), 100%B (10-15min)

0.4mL/min (6-7MPa), 37°C, Single Quad. MS (ESI)

5μL (10μg/mL, acetonitrile / water / formic acid = 50 / 50 / 0.1)