


Nardis ND-NX

250 x 4.6 mm

HPLC COLUMN APPLICATION

n-octyl-beta-D-thioglucopyranoside,
 n-octyl beta-D-glucopyranoside(Membrane protein solubilizer)

n-オクチル -β -D チオグルコピラノシド、 n-オクチル -β -D -グルコピラノシド
 (膜タンパク質可溶化剤)

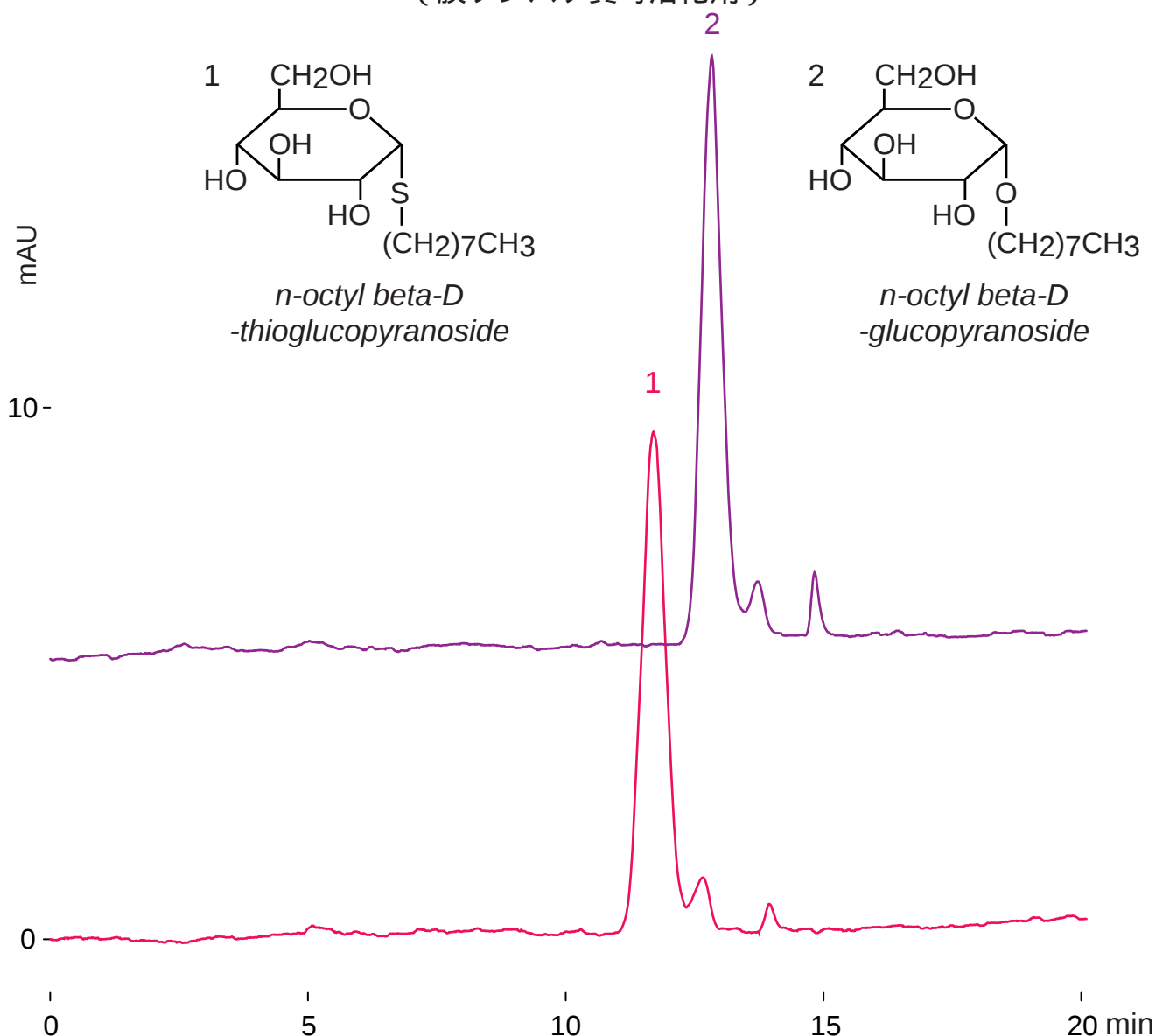

250 x 4.6 mm

A: acetonitrile

B: acetonitrile / 100mM ammonium acetate = 50 / 50

10-15%B (0-20min)

1.0mL/min (7MPa), 60°C

ELSD (spray chamber 70°C, drift chamber 100°C)

10μL (1mg/mL, water)