


Scherzo SM-C18

100 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Trialkylamine

トリアルキルアミン類のLC-MS分析

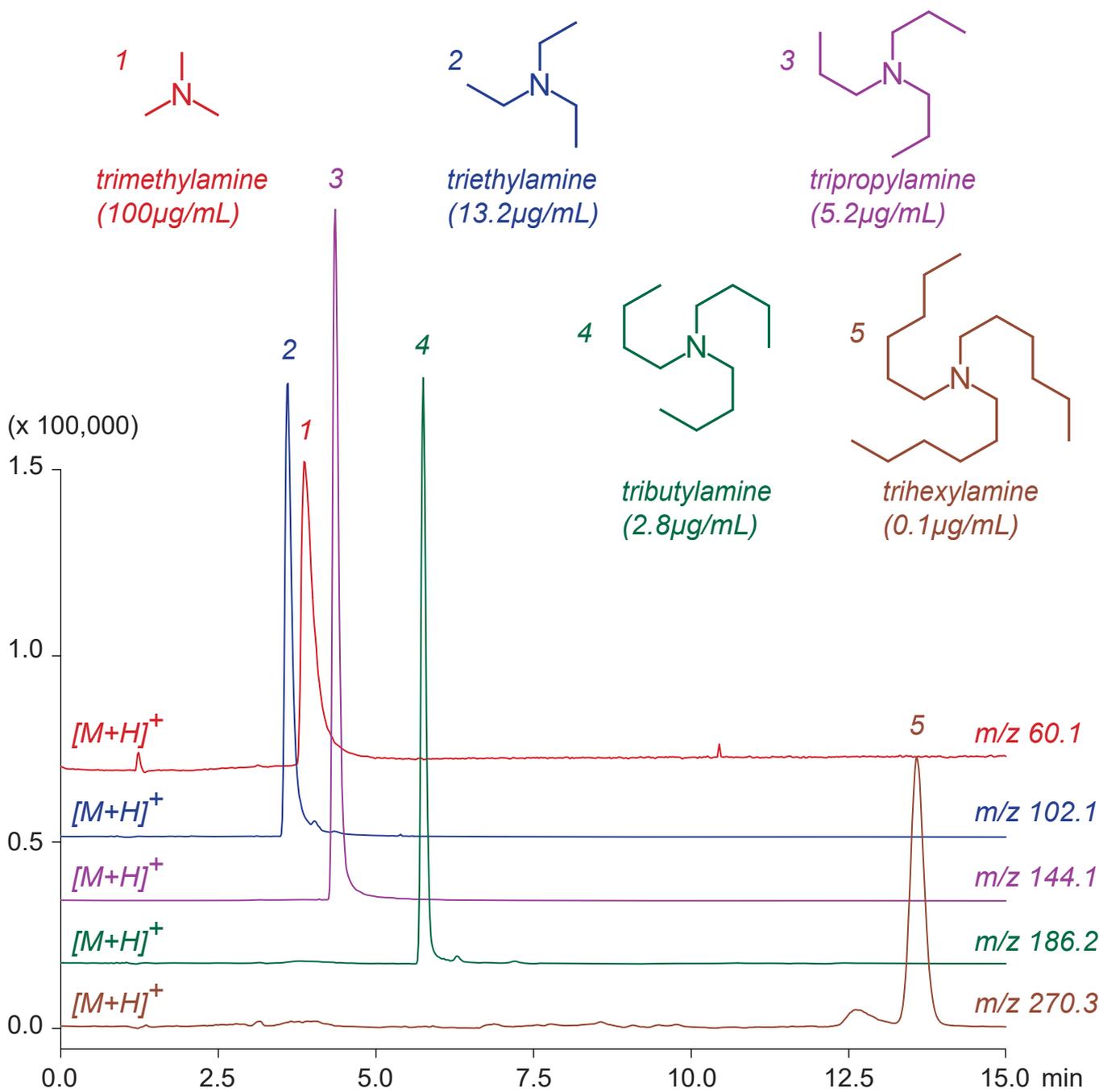

Scherzo SM-C18, 100 x 3 mm

A: 5mM ammonium acetate / methanol = 50 / 50

B: 100mM ammonium acetate / methanol = 20 / 80

0-100%B (0-4min), 100%B (4-15min)

0.4mL/min (13MPa), 37°C, Single Quad. MS (ESI)

5µL (0.1-100µg/mL, EtOH / 100mM AcONH₄ = 50 / 50)