

Unison UK-Amino


Nardis ND-NX

150 x 4.6 mm

HPLC COLUMN APPLICATION

Ribavirin (Antiviral)

リバビリン (抗ウイルス薬)

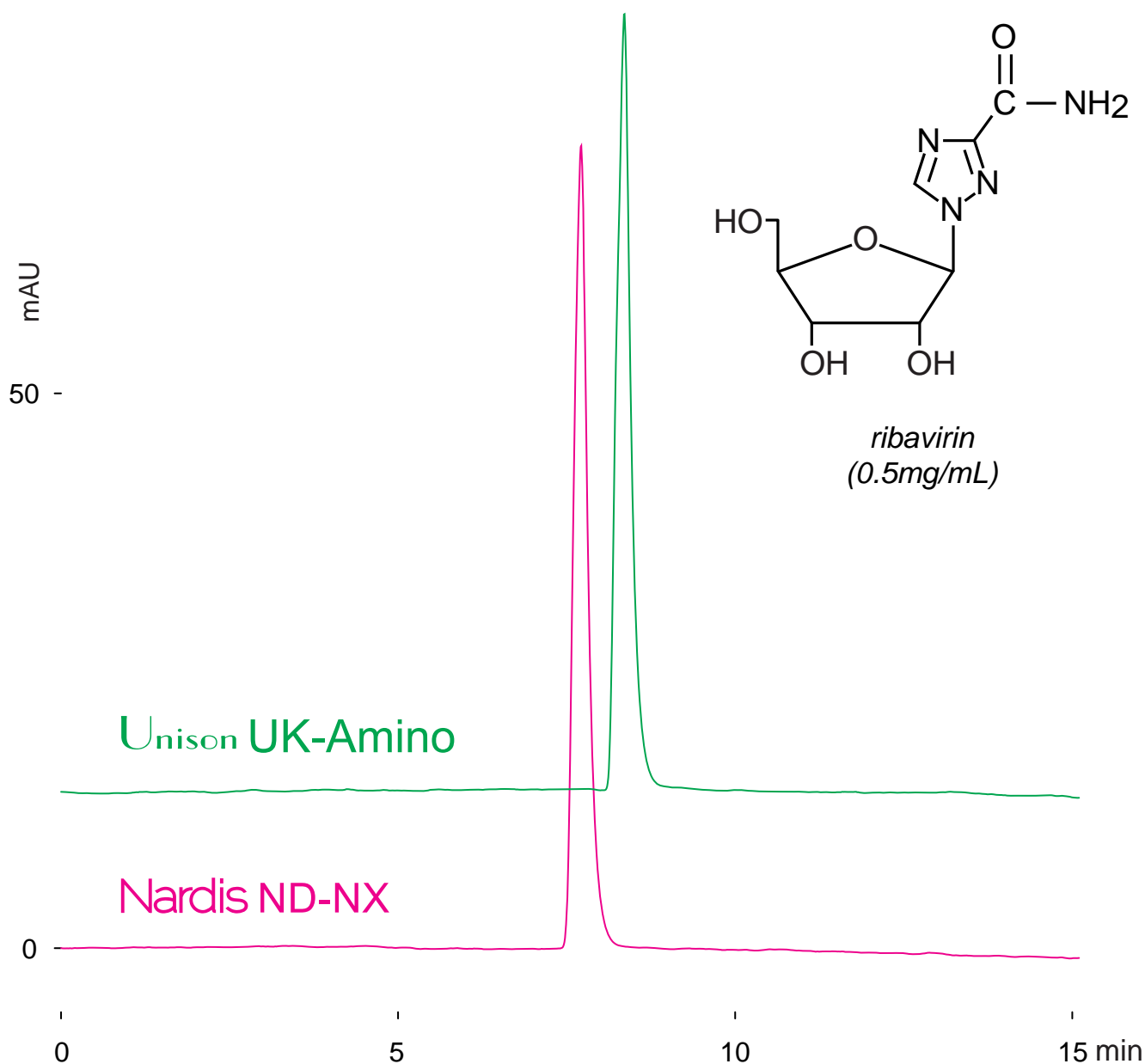

150 x 4.6 mm

A: acetonitrile / formic acid = 100 / 0.5

B: water / formic acid = 100 / 0.5

0-30%B (0-15min)

1.0mL/min (6MPa), 37°C

ELSD (spray chamber 70°C, drift chamber 100°C)

10µL (0.5mg/mL)