


Unison UK-C1
 Unison UK-Phenyl Cadenza CX-C18
 Unison UK-C8 Cadenza CW-C18
 Unison UK-C18 Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

1- (2-Thiazolylazo) -2-naphthol (TAN, Metal Indicator)

1-(2-チアゾリルアゾ) -2-ナフトール (金属指示薬)

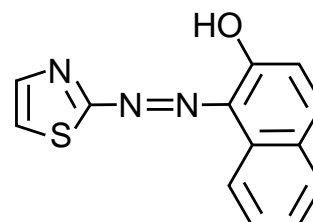

1- (2-thiazolylazo) -2-naphthol (TAN)
 (solution color : orange)

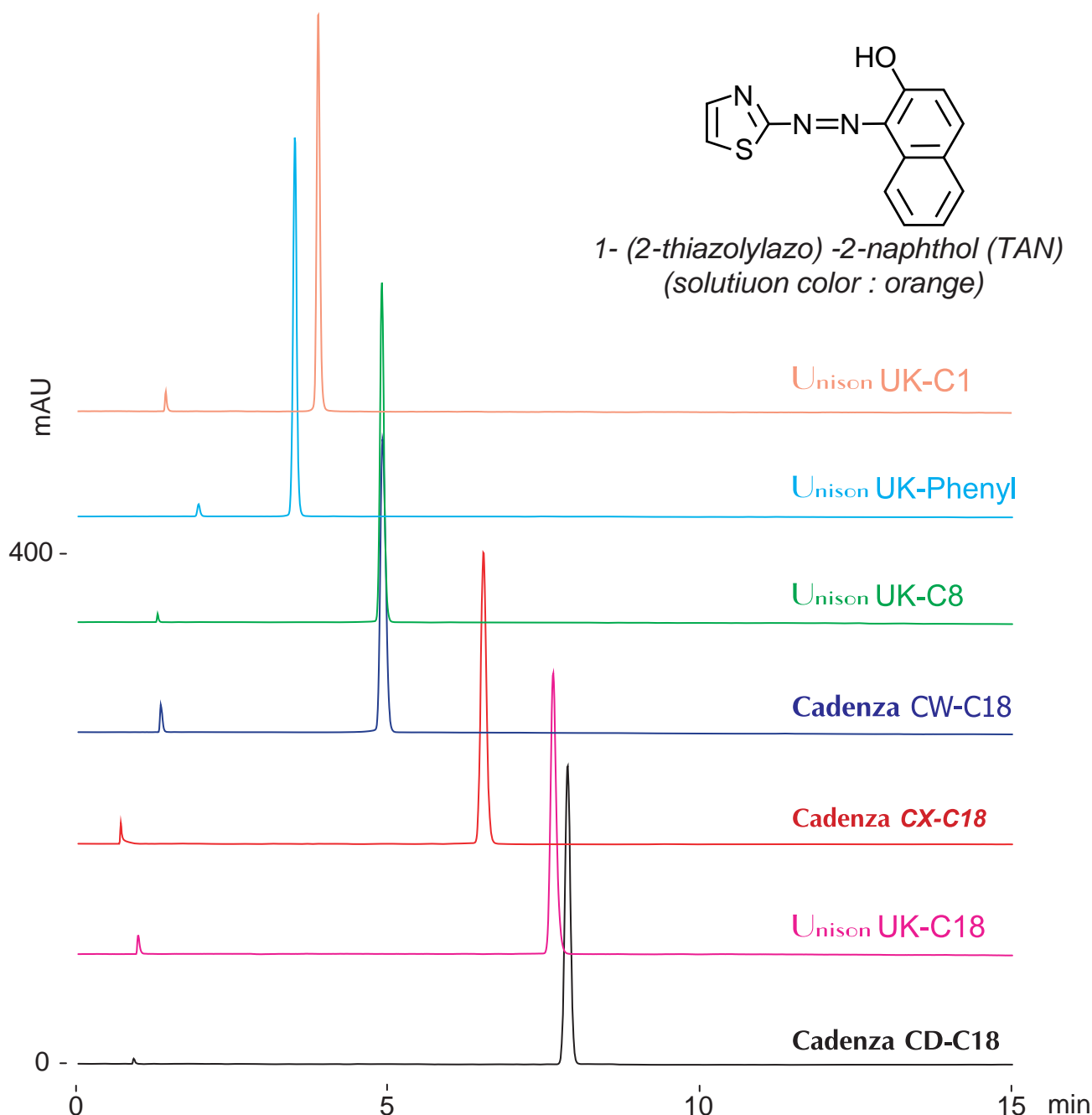

100 x 4.6 mm

A: water / acetonitrile / formic acid = 50 / 50 / 0.5

B: water / acetonitrile / formic acid = 10 / 90 / 0.5

0-100%B (0-15min)

1.0mL/min, 37°C, 450nm

10µL (0.05mg/mL, water / acetonitrile / formic acid = 50 / 50 / 1)