


Unison UK-Silica

150 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Guanidine hydrochloride salt

グアニジン塩酸塩のLC-MS分析

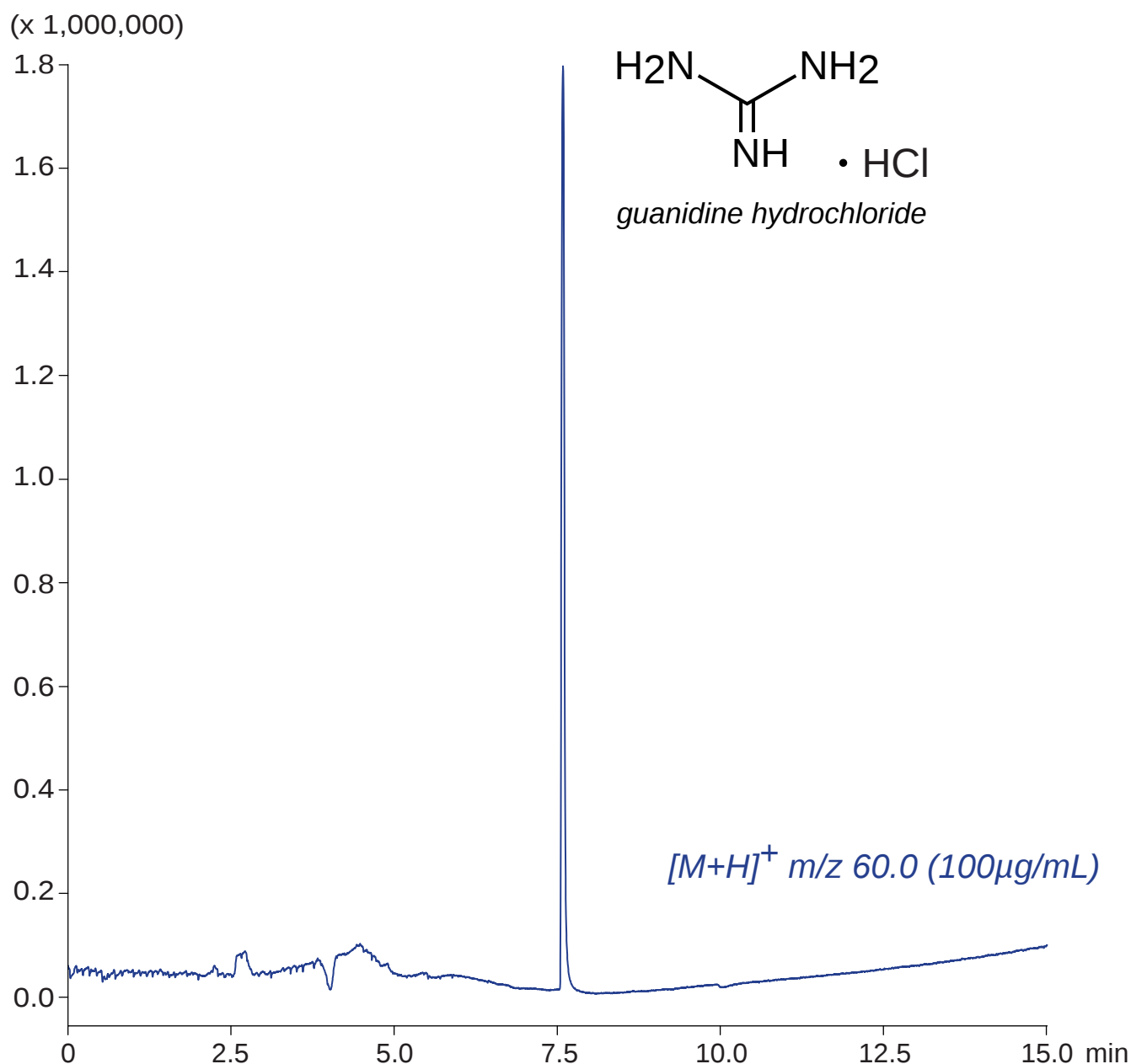

Unison UK-Silica, 150 x 3 mm

A: acetonitrile / water / formic acid = 95 / 5 / 0.1

B: acetonitrile / water / formic acid = 10 / 90 / 0.1

0-100%B (0-15min)

0.4mL/min (3MPa), 37°C, Single Quad. MS (ESI)

2µL (100µg/mL, water / formic acid = 100 / 1)